**INSPECTION SYSTEMS Series**

# INSTRUCTION MANUAL

Model
IS-01

Version: 1

Issued: 1.0.0

---

**IS-01**

Compact Inspection Video Microscope

**INSTRUCTIONS**

This instruction manual is for the operation guide, troubleshooting and maintenance to IS-01 compact Inspection Video Microscope. Please study this manual thoroughly before operating, and keep it with the instrument. The manufacturer reserves the rights to the modifications by technology development. On the basis of operation ensured, technical specifications may be subject to changes without notice

---

# Contents

## Hardware

1. Getting Ready	4
1.1 Operation	4
1.2 Maintenance	4
1.3 Components	5
2. Assembling	7
2.1 Assembling Scheme	7
2.2 Assembling Steps	8
3. How to Use	11
3.1 Adjust the Focusing Tension	11
3.2 Set Illumination	11
3.3 Place the Specimen	12
3.4 Adjust the Focus	12
3.5 Lock the Magnification	12
3.6 Use the Button and Interface	13
4. Troubleshooting	14
5. Optical Data	15

## Software

6. Overview	16
6.1 Software Start and Quit	16
6.1.1 Software Start	16
6.1.2 Software Quit	16
6.2 Main Commands	17
6.2.1 Command introduction	17
6.2.2 Command's icon introduction	17
7. Operation Instructions	19
7.1 Local browse	19
7.2 Colour and exposure adjustment	19
7.2.1 Exposure Adjustment	20
7.2.2 Colour Adjustment	20
7.3 Calibration	20
7.4 Drawing mark tool	25
7.5 AWB (Automatic White Balance)	26
7.6 WDR Function	26
7.7 Setting Button	27
7.8 Image real-time transmission	28
7.8.1 WiFi transmission	28
7.8.2 USB Transmission	30
7.9 Image / Video Switch	31
7.9.1 Image function	31
7.9.2 Video Function	31
7.10 Zoom Picture	32
8. Troubleshooting	33

---

# 1 Getting Ready

## 1.1 Operation

- 1) Do not expose the microscope in the sun directly. The microscope ought to be placed where is dry and clean. Avoid high temperature and violent vibration.
- 2) As the microscope is a high precision instrument, always operate it with care, and avoid physical shock or crash during the transportation.
- 3) To keep the image clear, please do not leave fingerprints or stains on the surfaces of lens.
- 4) Never rotate the left and right focusing knobs in reverse direction; otherwise the microscope will be damaged.

## 1.2 Maintenance

- 1) Keep all the surface of the lens clean. Wipe the lens gently with a soft lens tissue or blow away the dust. Carefully wipe off oil or fingerprints on the lens surfaces with tissue moistened with a little of 3:7 mixture of alcohol and ether.
- 2) Do not use organic solution to wipe the surface of other components, especially the surface of plastic components. Please use the neutral detergent if necessary.
- 3) Never disassemble the microscope yourself, otherwise it will influence its function or damage it.
- 4) After using, cover the microscope with a dust cover to prevent it from the dust, and store it in a place free from moisture to avoid rusting.

To keep the performance of the microscope, please check it periodically. For more details, please contact the agents nearby.

### 1.3 Components


---

## 2. Assembling

### 2-1 Assembling Scheme

Below is the Assembling Scheme to describe how to assemble the components, and the numbers denote the assembling order.

- ❖ Before assembling, make sure there is no dust, dirt or other materials which will disturb it. Assemble carefully and do not scrap any part or touch the glass surface.
- ❖ Preserve the hexagon spanner, and it will be used when changing the parts.

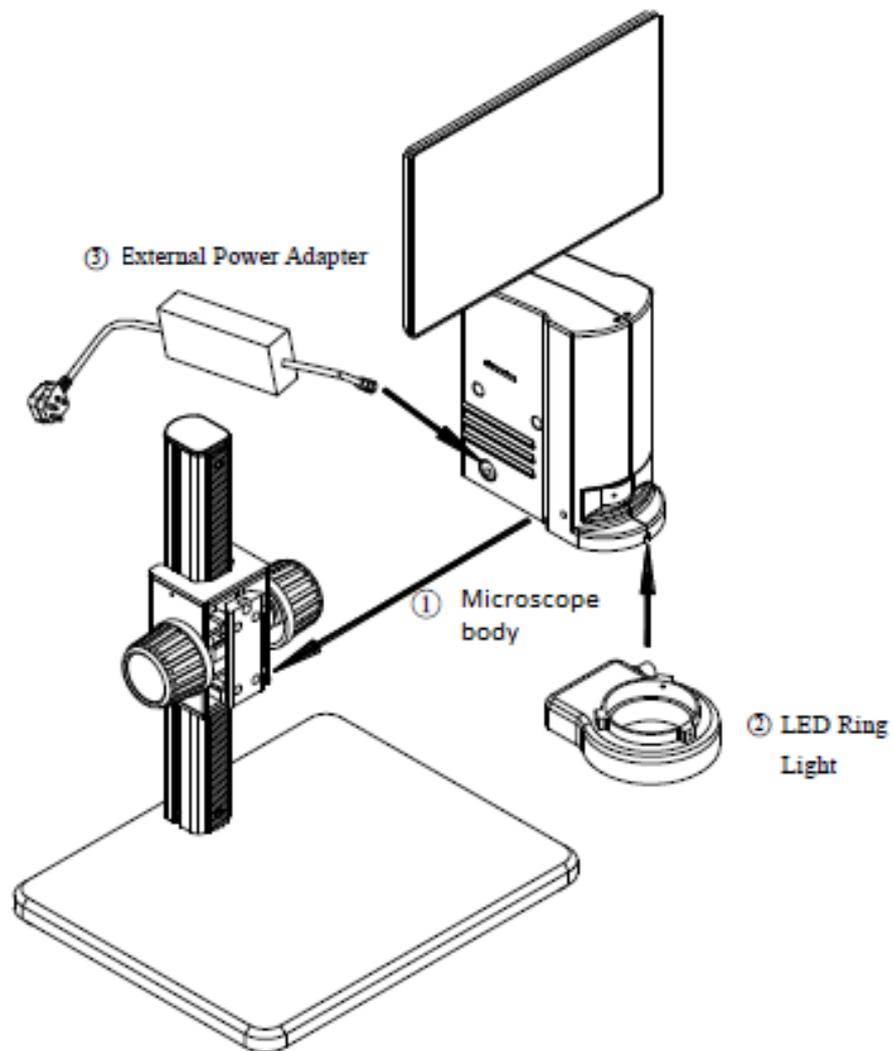## 2.2 Assembling Steps

Fig. 1

### 2-2-1 Assemble the extension ring (Fig. 1-2)

1. Unscrew the three screws on the bottom of the stand in order to remove the microscope arm.
2. Install the extension ring on the stand, then reinstall the microscope arm.
  - The extension ring is used to increase the length of the microscope arm, thus increasing the working distance.

Fig. 2

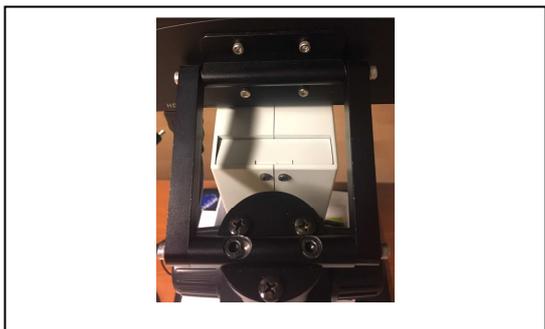
Fig. 3

### 2-2-2 Assemble the HDMI monitor (fig.3-4)

1. Unscrew the three screws on the top of the microscope arm, then remove the cap.
2. Install the provided cap by screwing the three screws.

3. Install the rectangular HDMI monitor holder using the two provided screws.
4. Assemble the HDMI monitor on the rectangular bracket using the provided screws.

Fig. 4

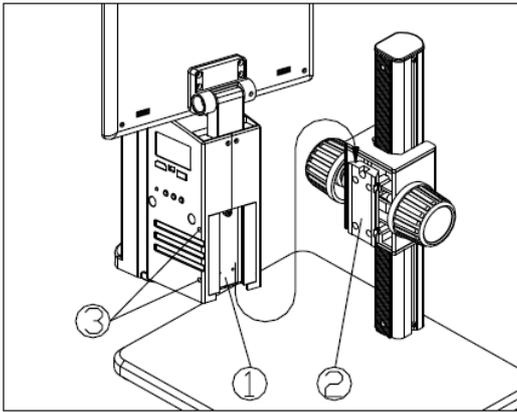
Fig. 5

### 2-2-3 Assemble the Microscope Body (Fig. 5)

1. Match the dovetail interface ① of the microscope body with the dovetail groove ② of the focus bracket group, and insert it from top to bottom as the direction shown in the figure.
2. Tighten the M4 inner hexagon screw on the body group ③ with a M4 inner hexagon spanner (side length 2mm).

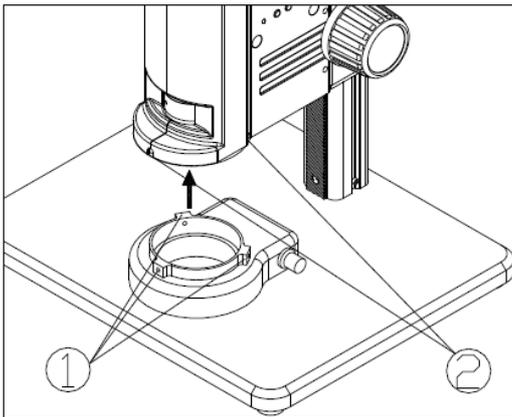
Fig. 6

### 2-2-4 Assemble the LED Ring Light (Fig. 6)

1. Move the LED ring light close to the microscope body according to the arrowhead direction, align the three screw holes ① with the thread groove ②, and lock the screw with a M4 inner hexagon spanner (side length 2mm).

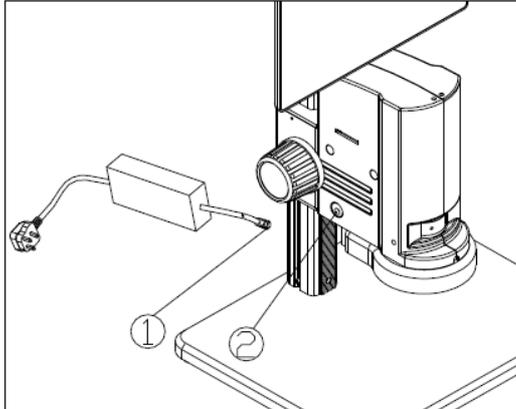
Fig. 7

### 2-2-5 Connect the Power Cord (Fig. 7)

1. Insert the external adapter plug ① into the power supply adapter socket ② of the microscope body to the bottom, and the power will be switched on.
- ❖ Don't use excessive force when the power cord is bended or twisted, otherwise it will be damaged.
  - ❖ Use the special wire supplied by our company. If lost or damaged, choose an external power adapter with the same specifications (power cord/charger).

Fig. 8

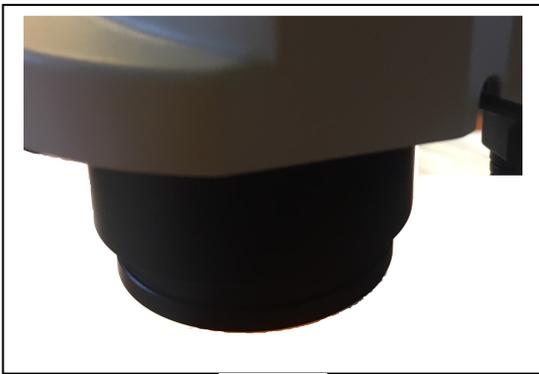
Fig. 9

Fig. 10

### 2-2-6 Installing Additional lenses (Fig. 8-9-10)

It is possible to install any of the available additional lenses (0.5x or 1.4x). To do this, please follow these steps:

1. Remove the LED ring light from the bottom of the microscope body. (Fig. 8)
2. Using the provided Allen wrench, loosen the screws placed on the front, on the right and on the left sides of the microscope body, in order to remove the LED ring light holder.(Fig. 9)
3. Unscrew the existing lens installed on the microscope and replace with the desired additional lens. (Fig. 10)
4. Reverse the disassembly procedure to reinstall the removed parts.

### 3. How to Use

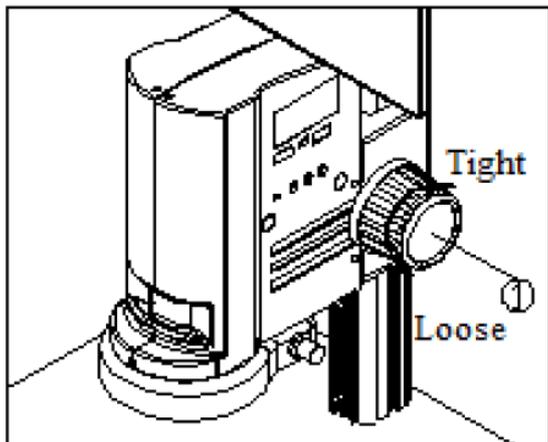
Fig. 11

#### 3-1 Adjust the Focusing Tension (Fig. 11)

1. To adjust the focusing tension, grip the left knob and rotate the right knob ①. It will be tighten when rotated in clockwise, while loosen in counter-clockwise.
2. Adjust the focusing tension to appropriate position, to prevent the microscope body declining with the bracket when observing, and also make focusing more comfortable.

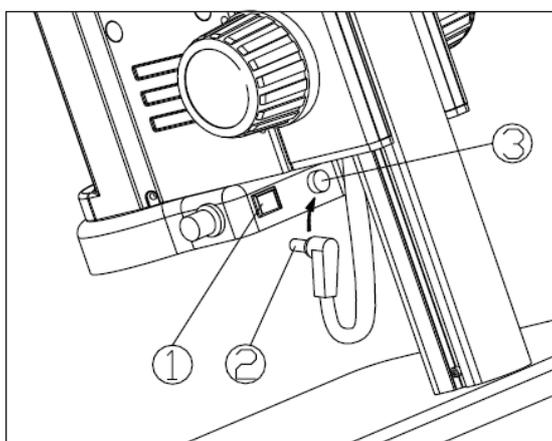
Fig. 12

#### 3-2 Set Illumination (Fig. 12-13)

1. Insert the power cord plug ② of body group into the ring light interface ③, then turn on the switch ① to start the light source.

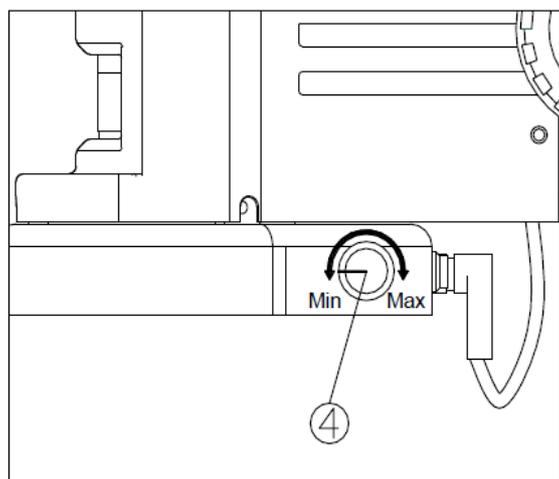
Fig. 13

2. Rotate the light adjustment knob ④ to adjust the illumination brightness.

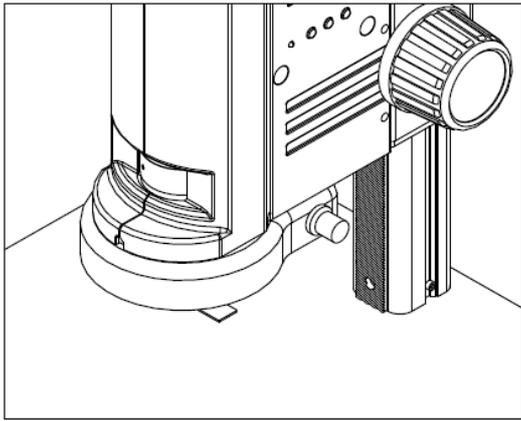
Fig. 14

### 3-3 Place the Specimen (Fig. 14)

1. Put the specimen on the base and let the observation point exactly below the microscope.

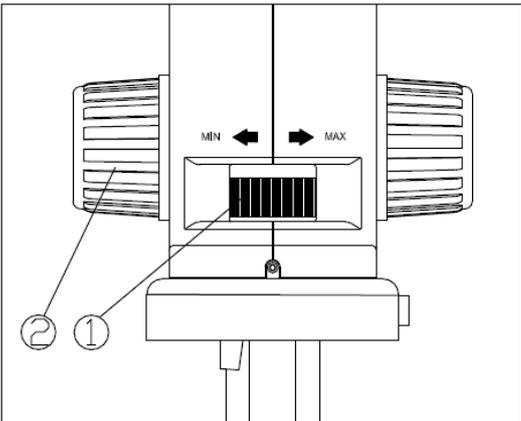
Fig. 15

### 3-4 Adjust the Focus (Fig. 15)

1. Rotate the zoom adjustment ring ① to the maximum magnification.
2. Observe the output image, if it is not clear, rotate the focusing knob ② to make it sharp.
3. Rotate zoom adjustment ring to the minimum magnification. Observe the output image until it is clear.

❖ **High low magnification parfocality is set to best position when this product is produced.**

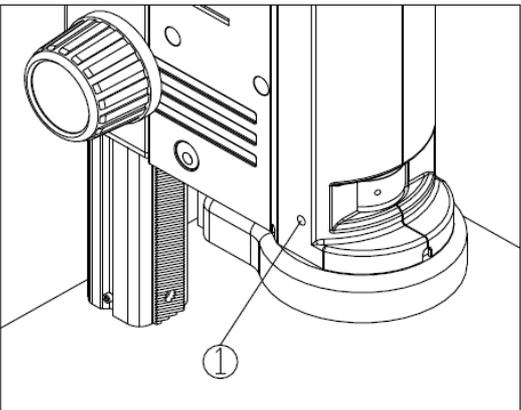
Fig. 16

### 3-5 Lock the Magnification (Fig. 16)

1. If use a fixed magnification for long term or use with shake, tighten and fix the magnification lock-screw ① with a M4 inner hexagon spanner (side length 3mm).

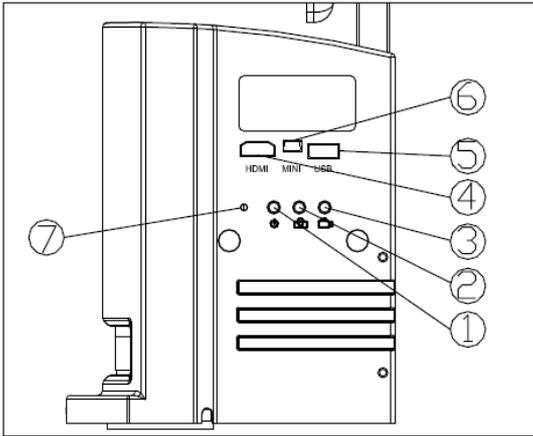
Fig. 17

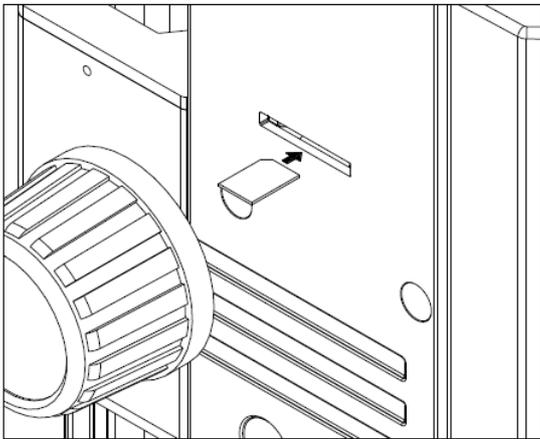
Fig. 18

### 3-6 Use the Button and Interface (Fig. 17-18)

1. Press the ON-OFF key ① for 3-5 seconds, the indicator ⑦ will turn from orange to green, which means the microscope is starting up.

❖ **It means the camera is working when the indicator flashes.**

2. Camera button ②: when the camera is working, press once to capture picture.
3. Video button ③: when the camera is working, press down for the first time to start recording, and press again to end recording.
4. HDMI cable interface ④, connect the HDMI cable to output the high definition image signal.
5. USB interface ⑤: connect the USB cable with the devices as mouse and USB disk.
6. MINI interface ⑥: connect the MINI cable to output the USB function image.
7. SD card is as shown in the figure, face the chamfer to the left and insert it to the bottom. (Fig. 18)

## 4. Troubleshooting

As the performance of microscope cannot play fully due to unfamiliar operations, the table below can provide some solutions.

<b>Problem</b>	<b>Cause</b>	<b>Solution</b>
1) Stain or dust is observed in the video.	Stains have accumulated on the specimen	Clean the specimen.
	Stains have accumulated on the CCD surface.	Clean the CCD surface.
2) The video image is unclear.	Stains have accumulated on the lens surface.	Clean the lens.
	Focusing is not correct.	Adjust the focusing.
3) The video image doesn't display	The power switch is not ON.	Turn on the switch button of microscope body.
	The external power cable is not well connected.	Check cable connections.
4) The image is too bright or too dark.	The LED ring light illumination has a wrong setting.	Adjust the brightness of LED ring light.
	Display properties setting	Open the display settings menu to adjust the brightness.
5) The display is no responding and flashes back.	Too many operations in the same time and some disorder, which results in error feedback of system.	Restart it.
6) The focus knob is not flexible	The focusing knob is too tight.	Loosen the focusing knob
7) The image is not clear due to the dropping of the microscope body.	The focusing knob is too loose.	Tighten it.
8) Microscope body cannot zoom	The zoom adjustment ring is locked.	Loosen the lock-screw
9) LED does not lights up when powered on.	No power.	Check cable connections
	The LED light is burnt out	Replace with a suitable light source.
10) LED light is burnt out suddenly.	Voltage is too high.	Use a suitable external power adapter.
11) The illumination brightness is not enough.	Voltage is too low.	Use a suitable external power adapter.

## 5. Optical Data

Objective	Optical Magnification	On-Screen Magnification	Horizontal Field of View (mm)	Vertical Field of View (mm)	W.D. (mm)
0.5x	0.35x	8.4x	26	16	170
	0.5x	12x	18.2	11.2	
	0.75x	18x	12.12	7.4	
	1x	24x	9.1	5.6	
	1.5x	36x	6.06	3.72	
	2x	48x	4.54	2.8	
	2.5x	60x	3.64	2.24	
1x	0.7x	16.8x	13	8	135
	1x	24x	9.1	5.6	
	1.5x	36x	6.06	3.7	
	2x	48x	4.55	2.8	
	3x	72x	3.03	1.86	
	4x	96x	2.27	1.4	
	5x	120x	1.82	1.12	
1.4x	0.98x	23.52x	9.28	5.71	100
	1.4x	33.6x	6.5	4	
	2.1x	50.4x	4.33	2.64	
	2.8x	67.2x	3.25	2	
	4.2x	100.8x	2.16	1.32	
	5.6x	134.4x	1.62	1	
	7x	168x	1.3	0.8	

Calculation done on 11,5" HDMI screen


---

## 6. Overview

### 6.1 Software start and quit


#### 6.1.1 Software Start

Click the icon application in the main menu to start.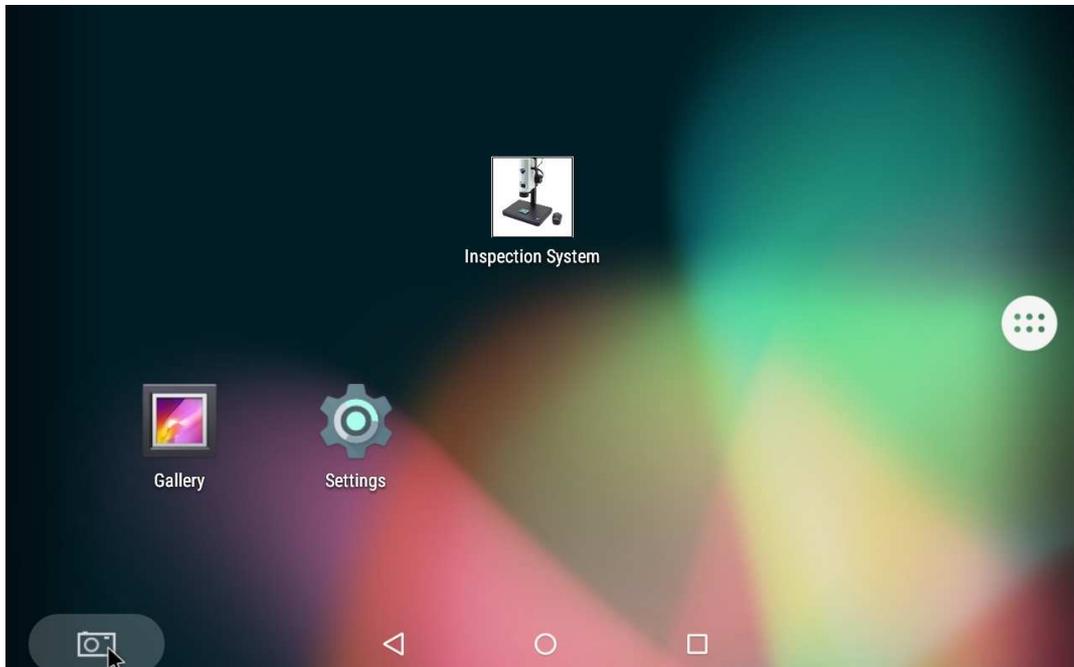

#### 6.1.2 Software Quit

Click the background button, all opened software in android system will be appeared.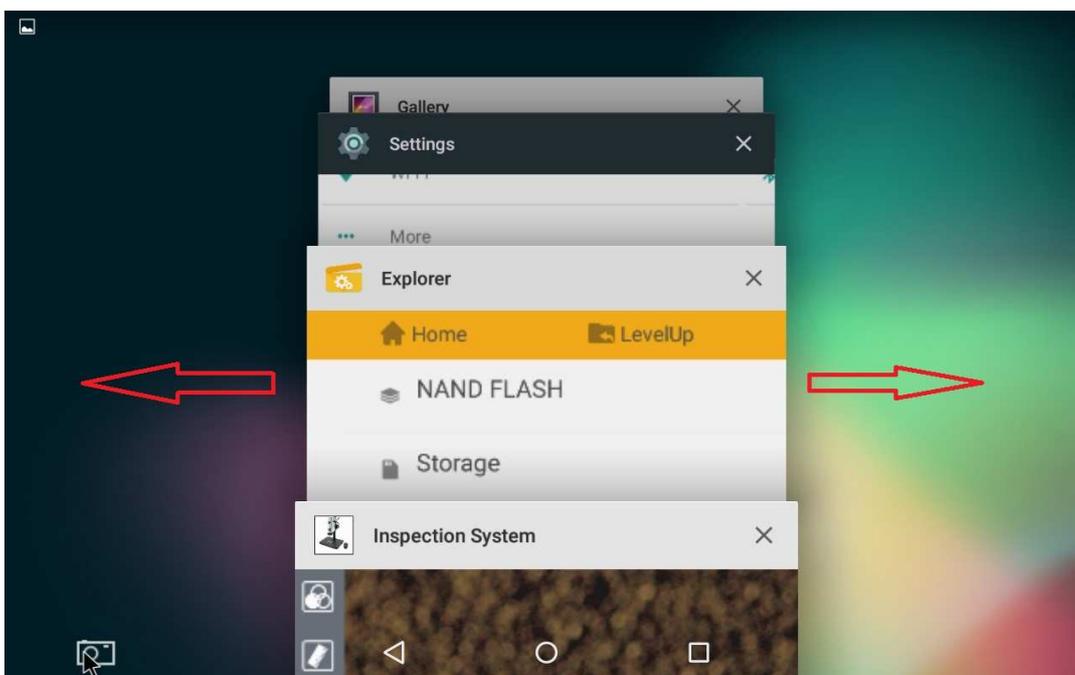


If need to quit, just click the exit button "X" in the top right corner of this software. Or gently swipe to the left or right to exit the program.

---

## 6.2 Main Commands

### 6.2.1 Command introduction

Start the program. Once started the program shows the following screen: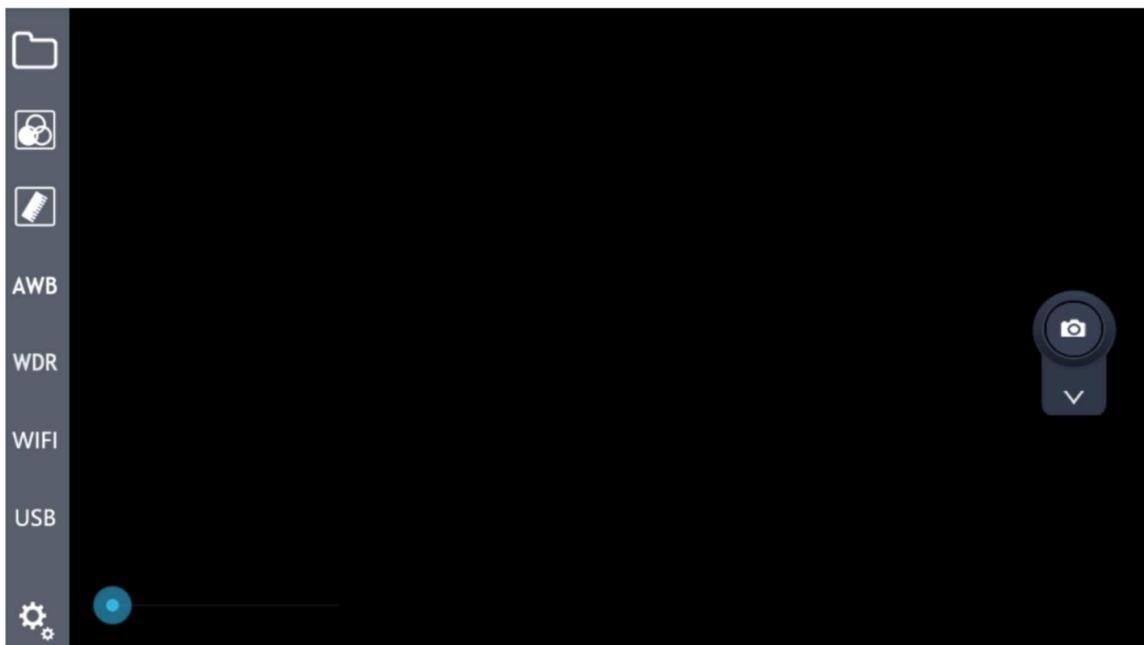

### 6.2.2 Commands icon introduction

A brief introduction to the current default icons:

 : Local browse

 : Colour and exposure adjustment

 : Drawing mark tool

 : White balance lock

 : WDR


---

 : WiFi output hotkeys

 : USB output hotkeys

 : Settings


 : Image / video switch hotkeys

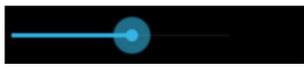 : Image zoom

## 7 Operation Instructions

### 7.1 Local browse

Click on the  button in the left menu bar, will jump to the local gallery. Click on the corresponding folder to view and edit the pictures and videos files: (delete images, colour effect transformation, etc.)

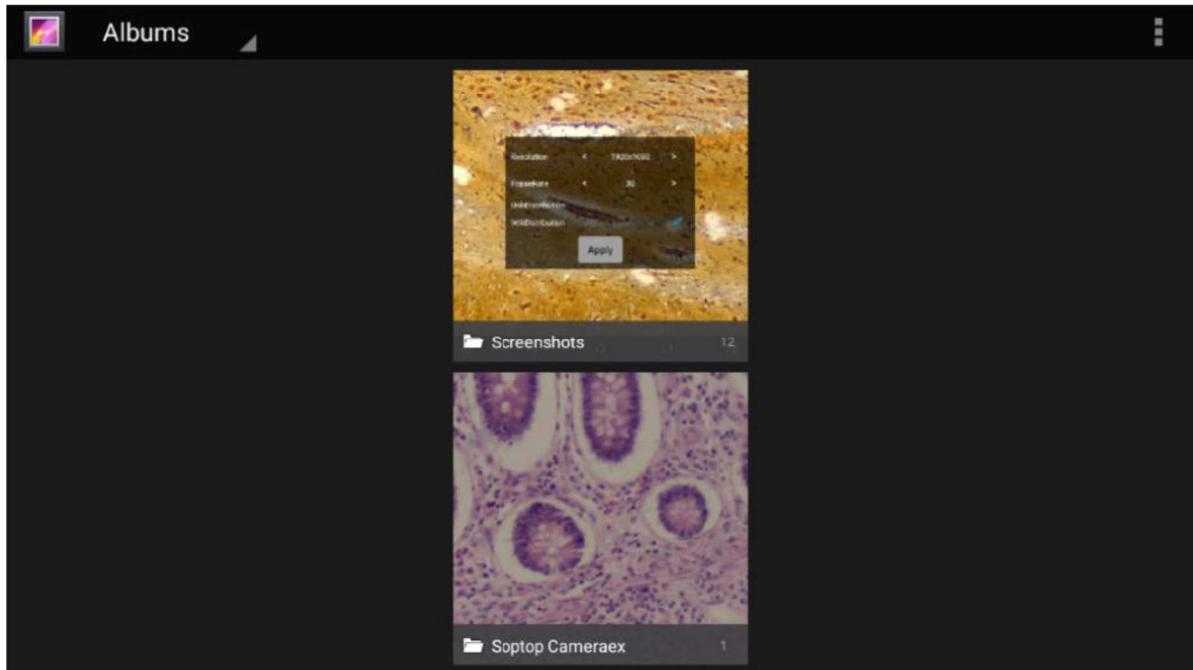### 7.2 Colour and exposure adjustment

Click on the  button in the left menu bar, will pop up a translucent view window in the main screen, including exposure adjustment and colour adjustment, as shown in the figure:

Exposure Adjustment

Colour Adjustment

## 7.2.1 Exposure Adjustment

Exposure adjustment region can select the exposure mode, Manual or Automatic exposure (AutoExposureLock “OFF” or “ON”), Exposure Compensation, Gain, Integration Time.

## 7.2.2 Colour Adjustment

Colour adjustment region affects the visual contrast adjustment under the different content through the microscope.

Colour adjustment includes brightness, contrast, saturation, chrominance, hue.

## 7.3 Calibration

In order to perform precise measurements, a system calibration is needed.


Click on the ruler button . A window will appear:


Click on “Edit” button


A new window will appear.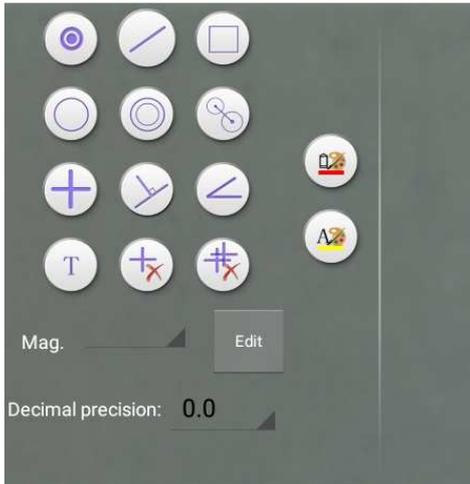


Click on "Edit" and a new window appears: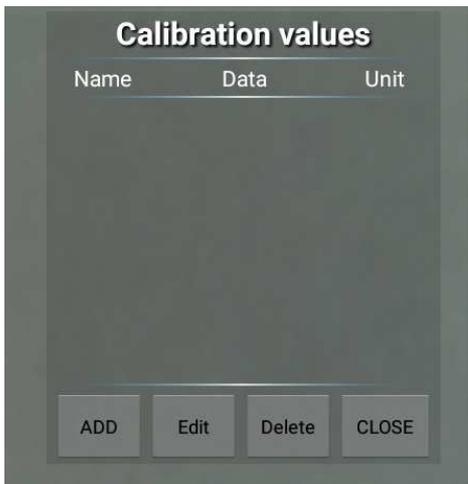


Now put a ruler or a known reference under the microscope and focus.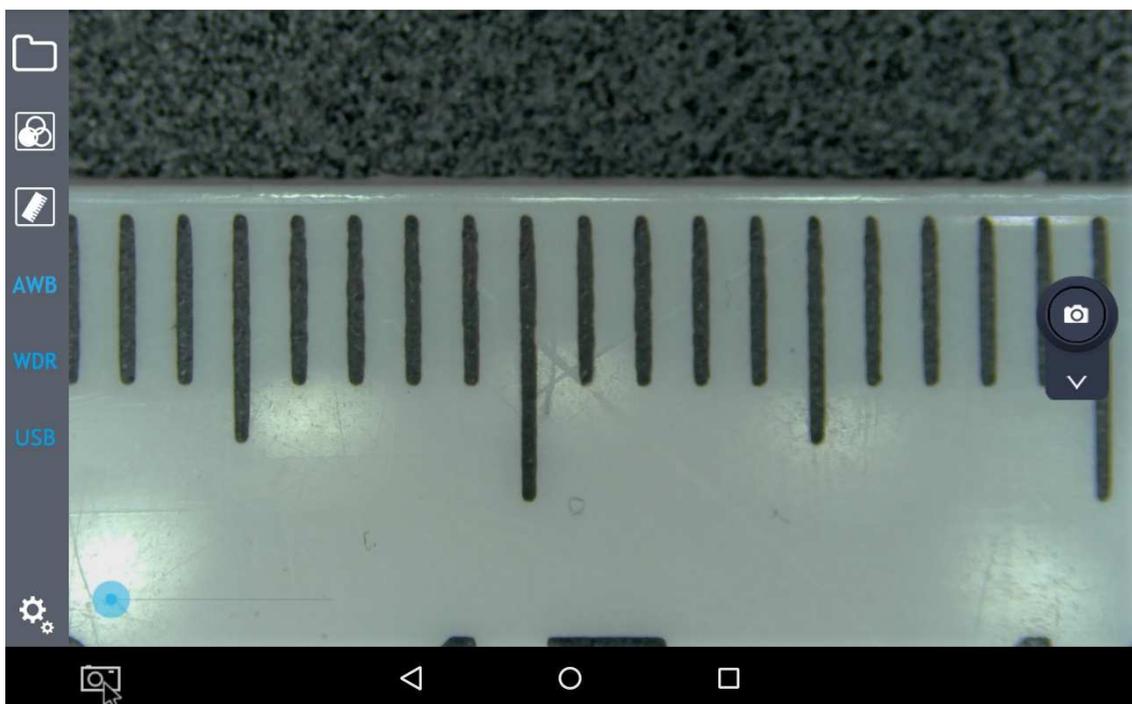


Click on "Add":

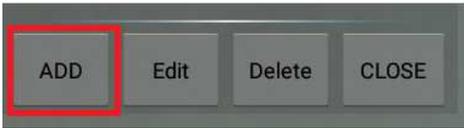
A new window will appear:

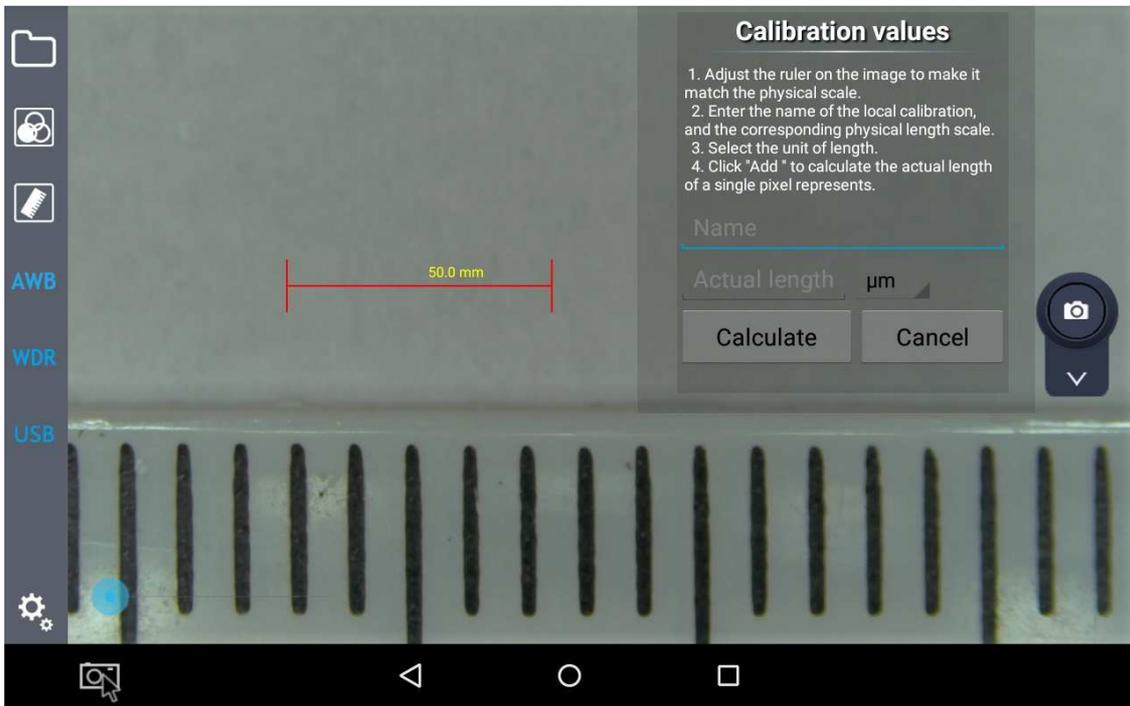
Move the red marker on the ruler

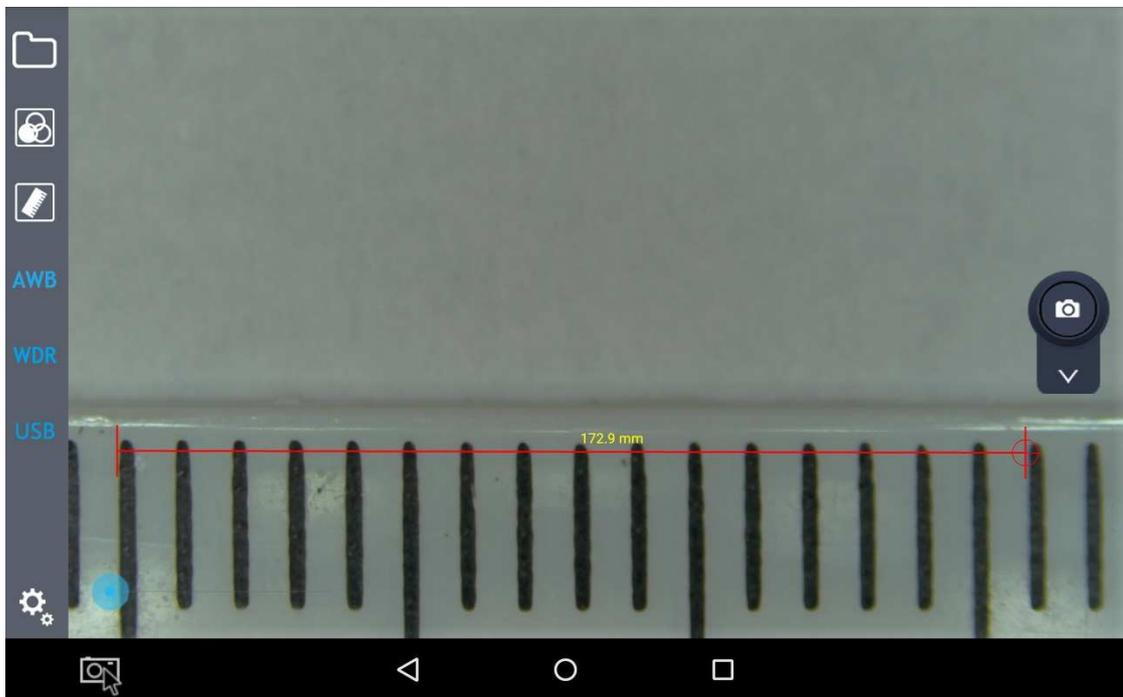and then click on "Name"

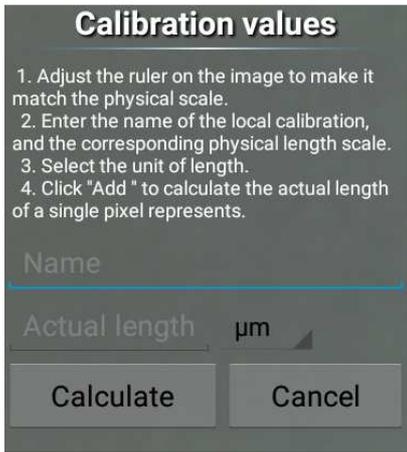An input dialog will appear where the user must input the actual magnification:

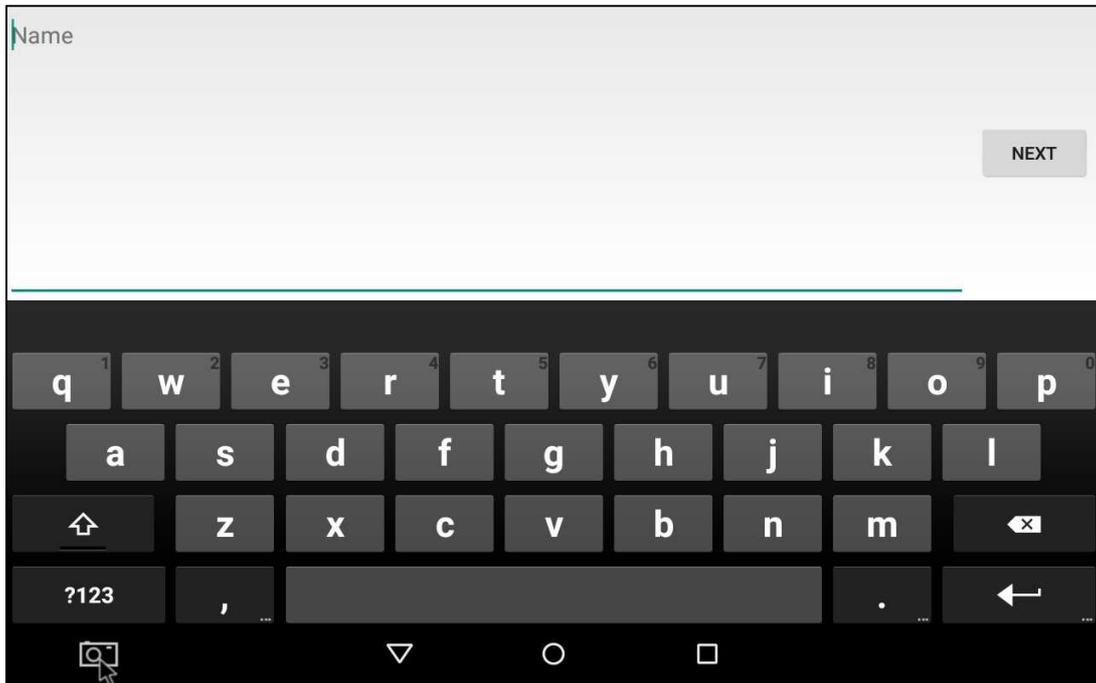
Insert the magnification (e.g. 0,7x) and click on "Next".  
Another input dialog will prompt to insert the length of the red marker (e.g. 12 mm)


Click on "DONE".

You will return to the previous dialog.  
Now select the unit ( $\mu\text{m}$  or mm)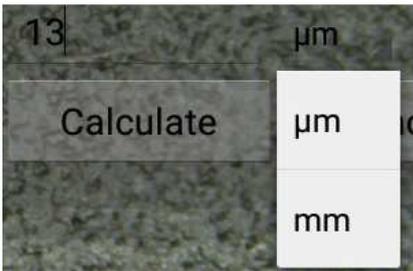


At the end of all the operations click on "Calculate". A dialog will appear: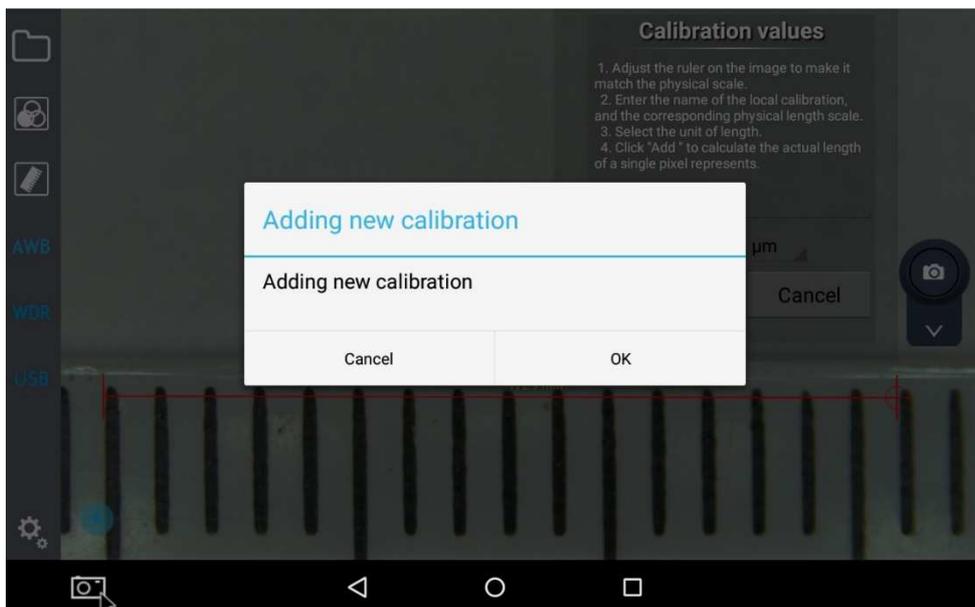

Press “OK” and the first calibration point has been saved.

Now repeat all these steps for each magnification of the microscope.

When all the magnifications will be calibrated, the calibration list will be available:

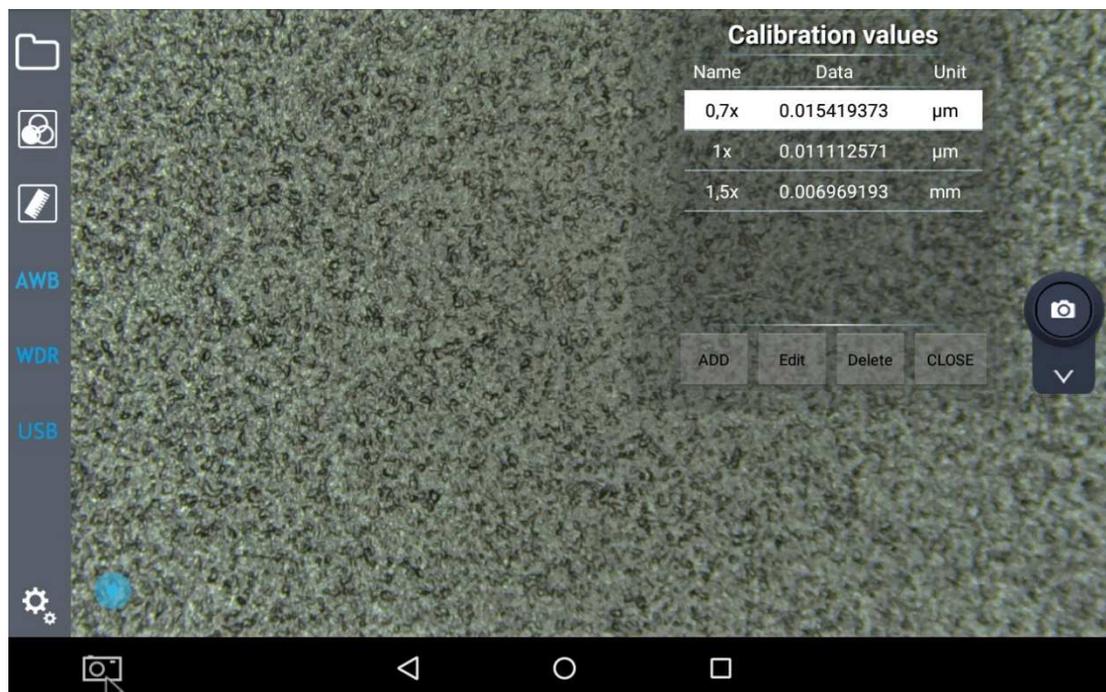Before performing any measurement, it is important to open the related calibration according to the magnification in use.

#### 7.4 Drawing mark tool

On the displayed images the user can have many useful function.


These can be activated by clicking the ruler button .

Click on drawing mark tool button on the left menu bar, and then many kinds of drawing marker will appear in the top right of frame.


Commands are the following: point , straight line , rectangle , circle , concentric circle , center distance , cross line , perpendicular line , angle , text , delete single item (delete last step or delete after mark) , delete all .

By clicking the line with mark , or character colour , another option window will appear showing a colour palette with six colours (yellow, black, white, green, red and blue)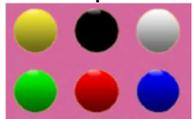

By selecting one of these colours will change the colour of the selected item. When use image function, the user can save the marked picture into the local folder along with the image (the marked video will not be saved when recording).

The effect of drawings on the image can be like this: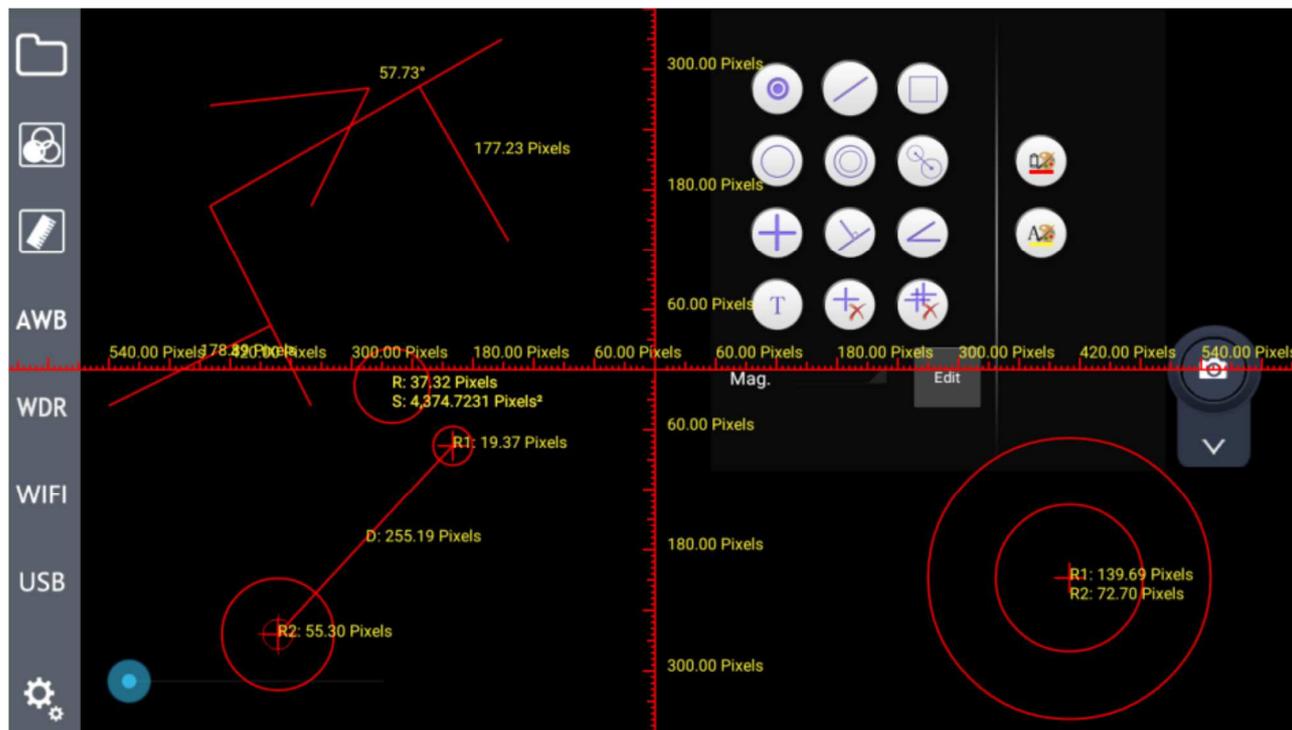

Once the system is calibrated, the user can select the unit for the displaying of the drawn lines (microns, millimeters, and so on)

#### 7.4 AWB (Automatic White Balance)

Click on the  button on the left menu to start AWB (Automatic White Balance).

When an image is taken under fluorescent bulb, it will appear with a greenish tinge; under halogen light it will be yellowish, while under sun light takes a blue tinge: the reason lies in the setting of "white balance".

The role of the white balance is to recover the images' normal colour in these scenes.


AWB is based on image algorithm, so when you start the automatic white balance with the microscope, you should choose white areas to establish the "white".

Put a white paper under the microscope: the system will automatically calculate the colour temperature and the results of the correction is that the image's average colour is white.


#### 7.5 WDR Function

WDR, acronym for **Wide Dynamic Range**, means that brightest and darkest parts of the image can be observed clearly at the same time. When WDR function is activate, system will improve the brightness of too dark area, especially under the stereomicroscope. Click on the button  on the left menu bar, the button will

change to **WDR**, and there is a prompt message for opening the WDR function in the center of the screen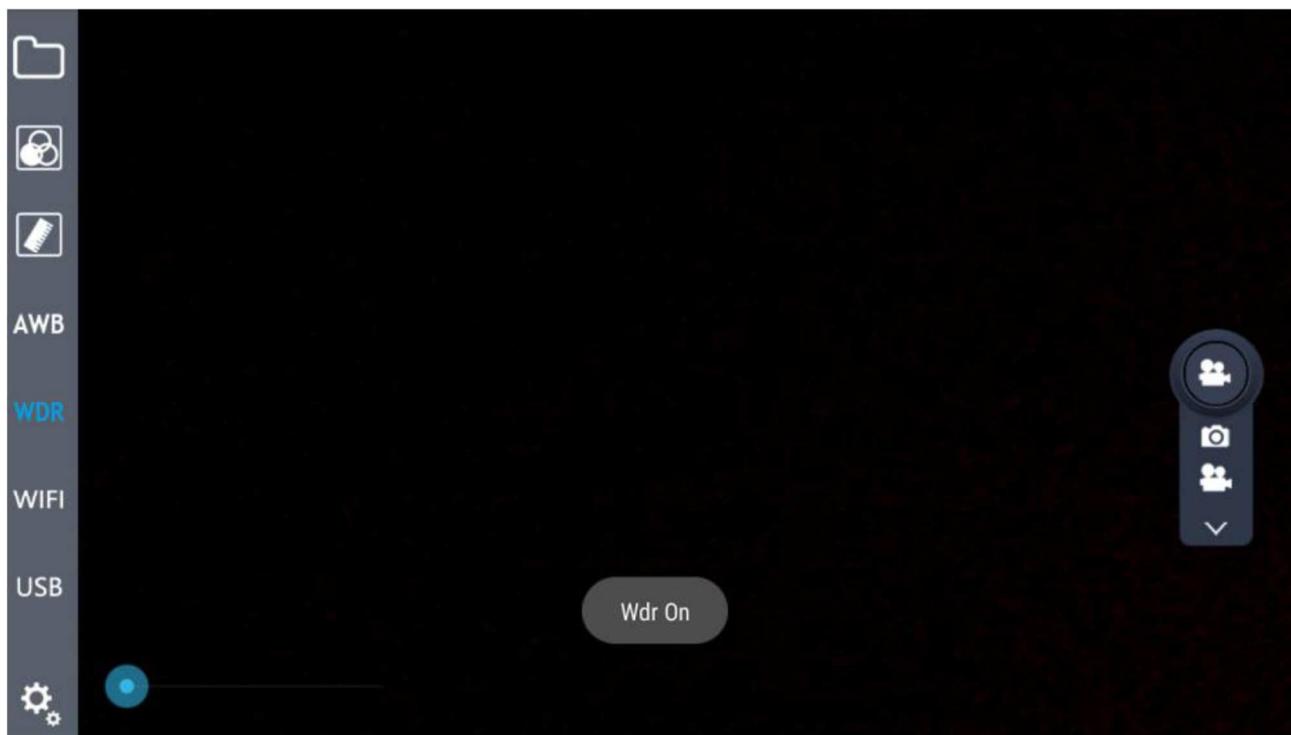


WDR function is active

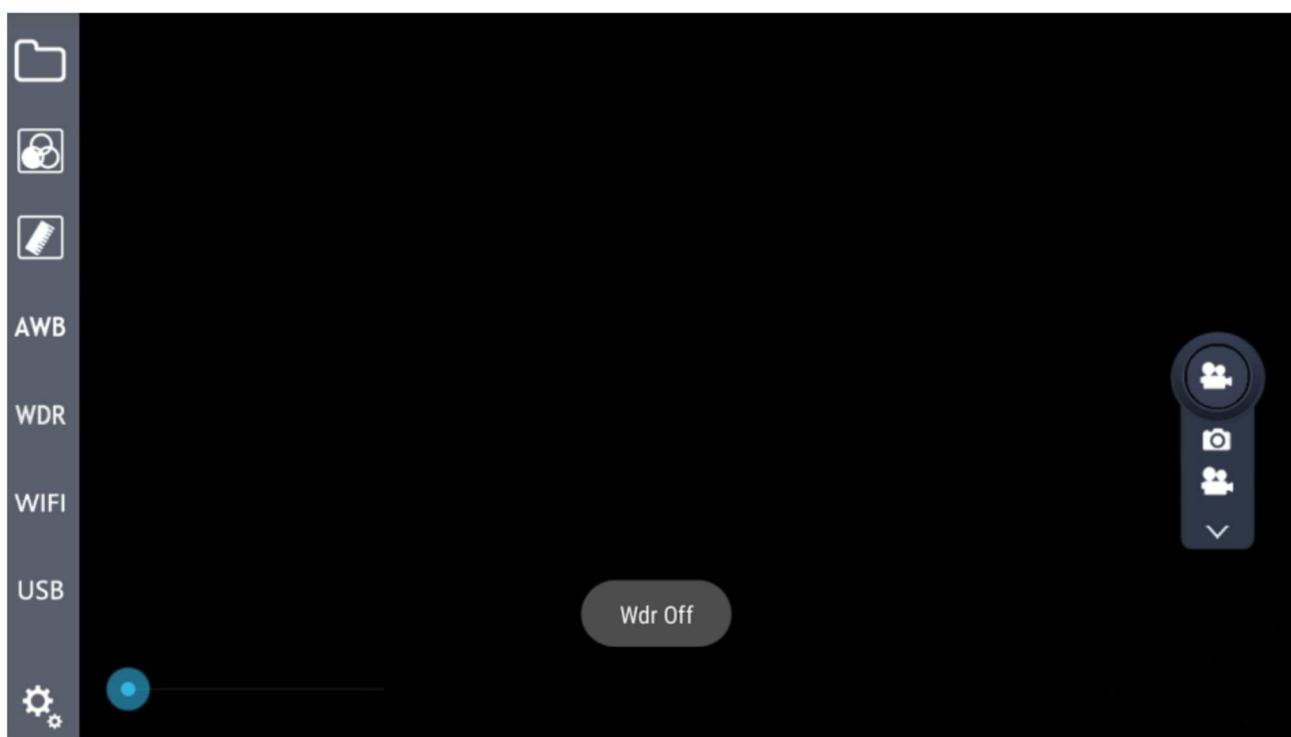WDR function is not active

## 7.7 Setting Button

To open the setting function, click on  button: a translucent window appears as shown: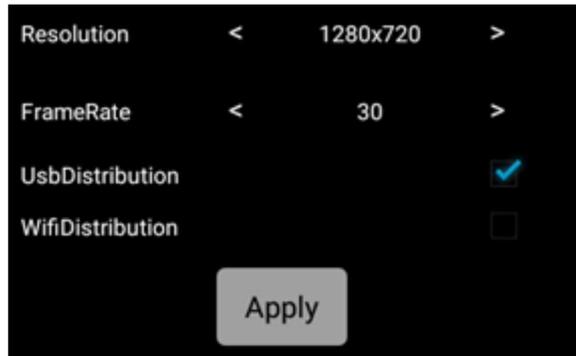

Setting module includes Resolution, Frame Rate and Transmission mode.

The Resolution **Resolution** has 4 options, respectively: 2688\*1520, 1920\*1080, 1280\*720, 672\*376.

The Frame Rate value **FrameRate** can be selected among 0FPS / 15FPS / 30FPS.

The Transmission mode has two options: WiFi **WifiDistribution** and USB transmission **UsbDistribution**.


## 7.8 Image real-time transmission

### 7.8.1 WiFi transmission



The software can not only capture and display the images in local, but also can do real-time transmission by external devices. WiFi hotspot transmission is not only one-to-one, also support multi-user to share. It only requires the shared equipment and system are connected to the same network address. For it, we can through spontaneous hotspot or through external WiFi hotspots.

1. In case of spontaneous hotspot, you need to open the network share hotspot service in system settings. For further details please refer to the Step 1:

Step 1: Spontaneous hotspot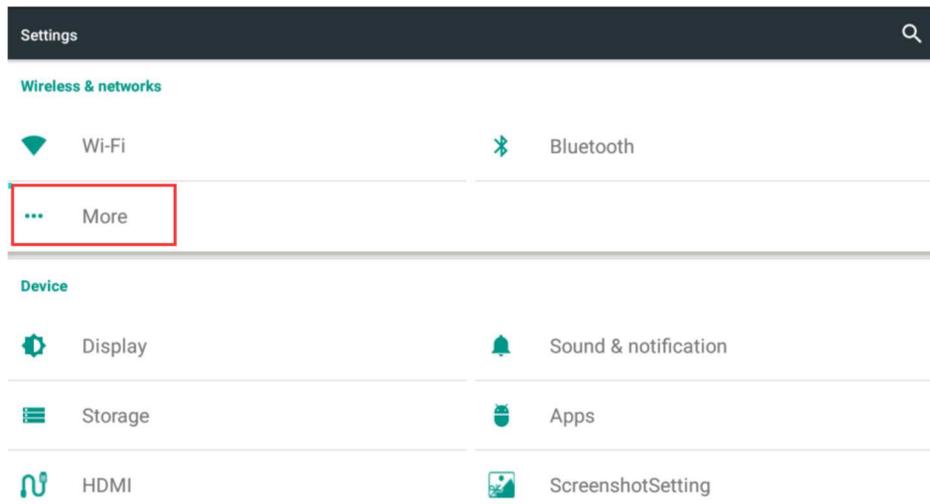


Open the “More” option in settings

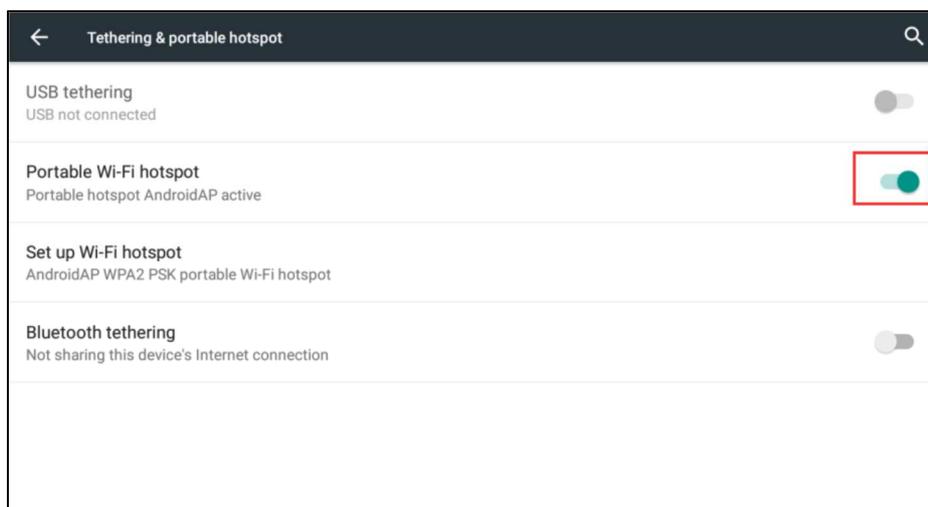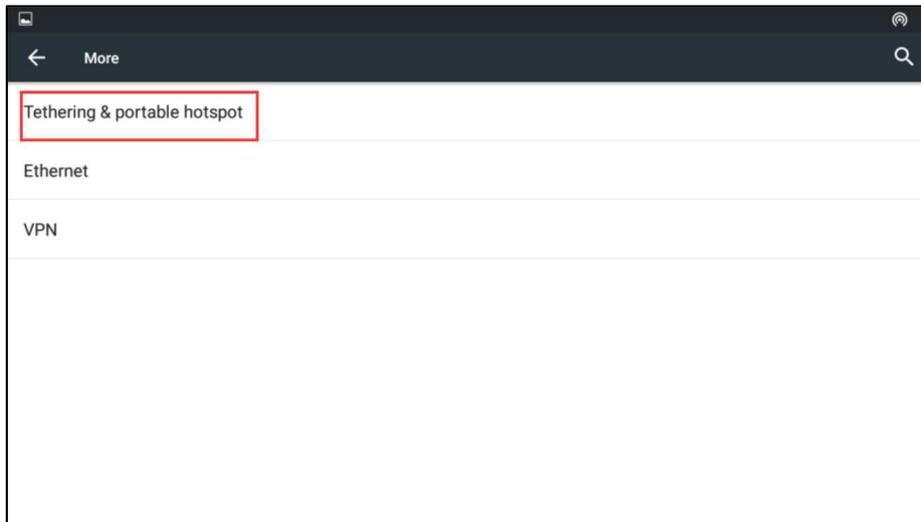
Open the portable WLAN hotspot option

2. In case of an external WiFi hotspots, you may need to open the system Settings of WLAN option, even on the need to connect the WiFi. For further details please refer to the Step 2:

#### Step 2: External WiFi hotspot

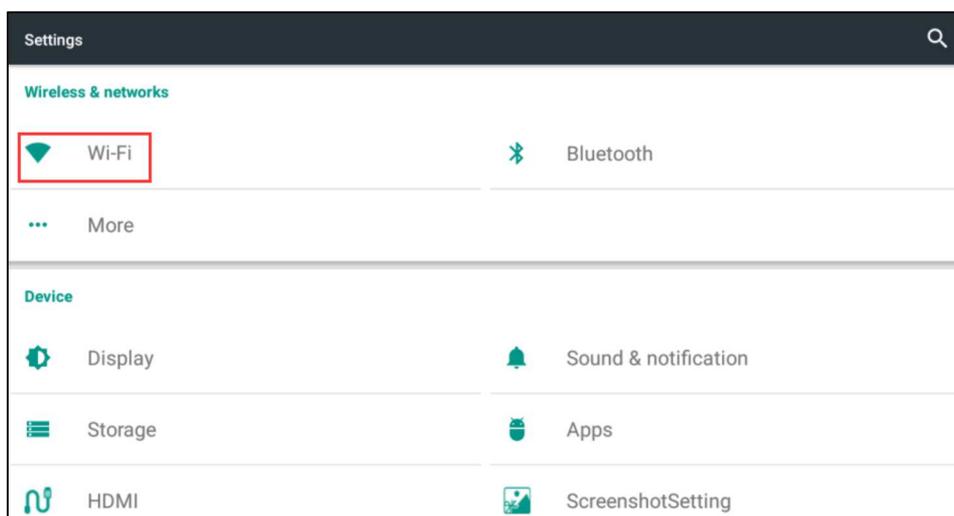Open the "WLAN" in settings

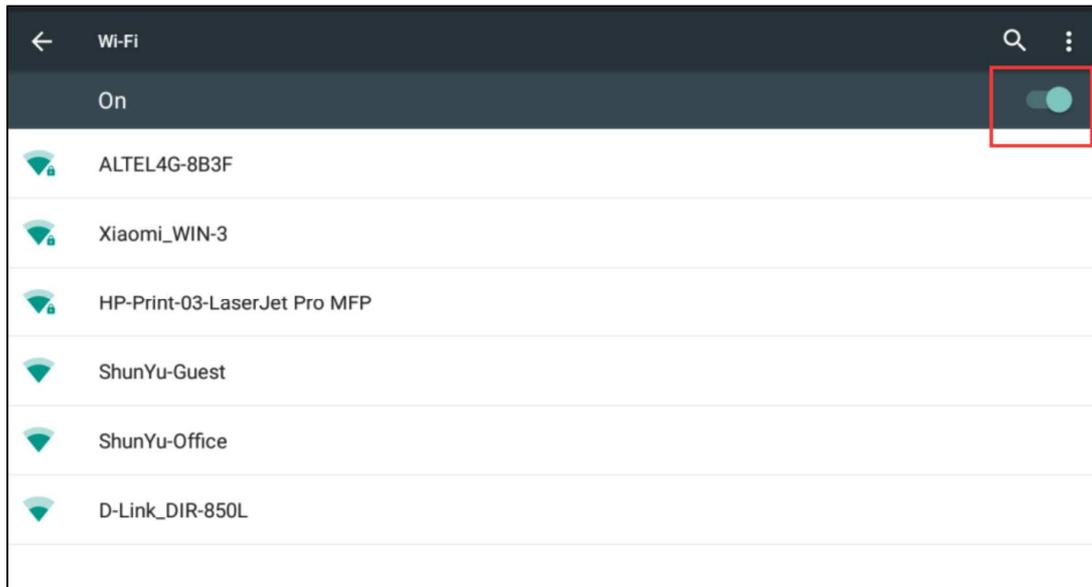
Choose the shared network after “WLAN”opened.

Now open **WifiDistribution** or WiFi shortcut key **WiFi** on the left menu bar, click on **WifiDistribution** in Settings: a string of IP address will appear **rtsp://192.168.43.1:8554/0/main**.

External devices through player software can directly enter the IP address in input field, then click play, the devices can get synchronized images as Inspection System.

### 7.8.2 USB Transmission

Real-time image transmission can be obtained through the USB to data transfer: the first operation is connect the external mini USB interface of the module to the USB interface of receiver device, then click on the USB output option **UsbDistribution** in the Settings,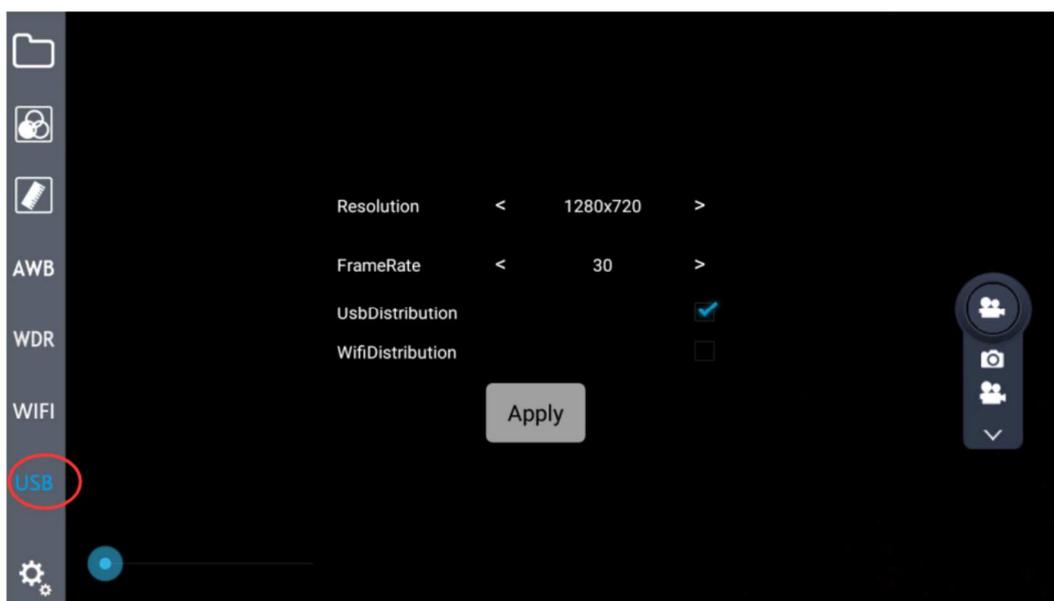

---

or click on the shortcut button  on the left menu bar, and then open the installed software in receiver device, choose the official version of the video capture program AMCap software, now you can use real-time image transmission function.


## 7.9 Image / Video Switch

### 7.9.1 Image function


Choose the image button in Image / Video selection box . Click on the Image icon . If the screen is showing the pictures' storage place, it means that the image acquisition is done. You can directly access to the gallery application to view the pictures, and also can click the local browse button  to view.

### 7.9.2 Video Function

Like the Image function, choose the Video button in Image / Video selection box .Click on the Video icon : now the icon becomes red , and the screen will show . This is the information that the video recording is active. Clicking once again the icon it will return to its original status  and the screen will show . This means that the video recording has been completed. You can go to the gallery or click local browse  to view.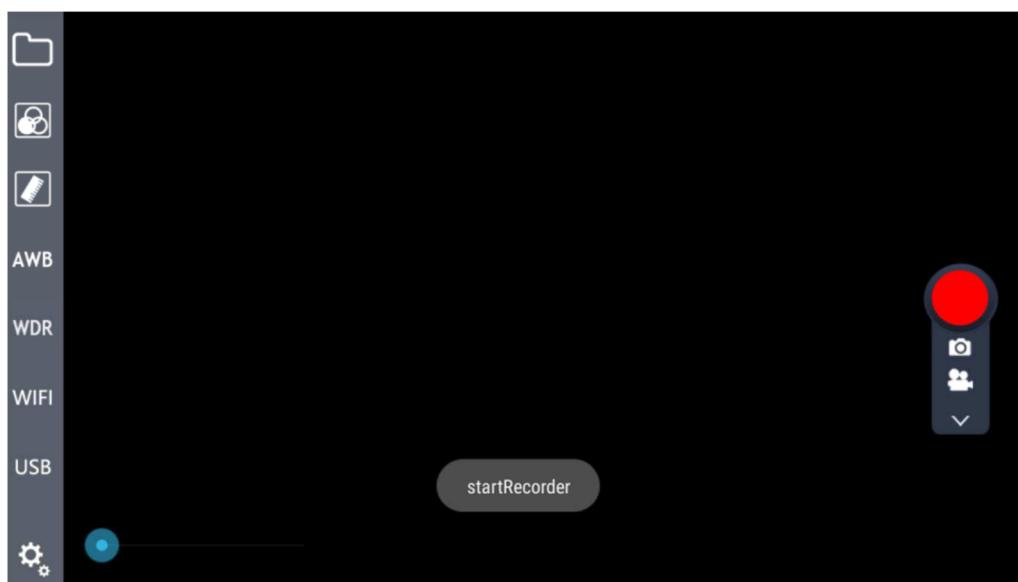

Start Recording

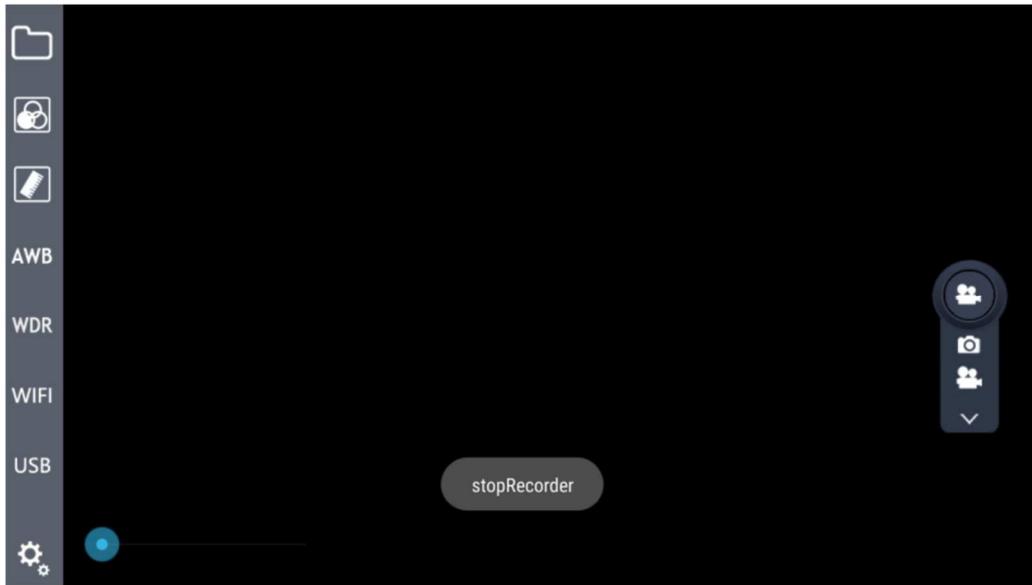Finish Recording

### 7.10 Zoom Picture

You can enlarge the picture when browsing. This function is done through the bar-type control , by moving from left to right the blue dot.

While performing measurement, zoom function has to be set at its minimum: zooming the image will result in errors during measurements.

---

## 8 Troubleshooting

As the performance of software cannot play fully due to unfamiliar operations, the table below can provide some solutions.

<b>Problem</b>	<b>Cause</b>	<b>Solution</b>
1) Clicking on WiFi button has no effect and does not appear the IP address.	Hotspot sharing not open	Open the hotspot sharing in settings (refer to step 1)
	WLAN is not open.	Open the WLAN in settings (refer to step 2)
	Resolution ratio (1520p) does not support WiFi transmission	Change to the feasible resolution in the program Settings
2) Camera program is not responding or forced termination	Too frequent and irregular operation cause system error response	Restart
3) Image distortion when opening AWB	Wrong settings	Start the automatic white balance in the correct area that close to white colour.

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**OPTIKA S.r.l.**