SZP Series

INSTRUCTION MANUAL

Model
ST-170

Ver. 1.0 2019

Contents

1. Warning	2
2. Symbols and conventions	2
3. Safety Information	2
4. Intended use	2
5. Overview	3
6. Unpacking	4
7. Assembling	4
7.1 Assembling procedure	5
8. Use of the attachment	6
8.1 Light path selection	6
Equipment disposal	7

1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

2. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

3. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

4. Intended use

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

5. Overview

This manual is only for trinocular port attachment.
For information on the microscope, consult the manual provided with the microscope.

PHOTO/TV
PORT

LIGHT PATH
SELECTOR

TRINOCULAR
PORT

6. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

7. Assembling

Once opened the box, the microscope parts are the following:

- ① Trinocular port attachment
- ② Photo port for projection lens
- ③ Photo adapter for "C"-mount adapters

7.1 Assembling procedure

1. Install the trinocular port attachment attachment on the microscope body, aligning the guide under the attachment with the pin on the microscope body.
2. Tight the locking screw ① using the provided Allen wrench. (Fig. 1)

3. Install the observation head aligning the guide under the head with the pin on the microscope body and tight the locking screw ② using the provided Allen wrench. (Fig. 2)


Installing camera with projection lens

4. Insert the bottom part of the projection lens ③ in the empty hole of the photo/TV port and screw the fixing knob ④. (Fig. 3)

Installing camera with "C"-mount adapter

5. Remove the photo/TV port unscrewing the fixing screws ⑤.
6. Insert the photo adapter for "C"-mount ⑥ and lock the fixing screws ⑤. (Fig. 4)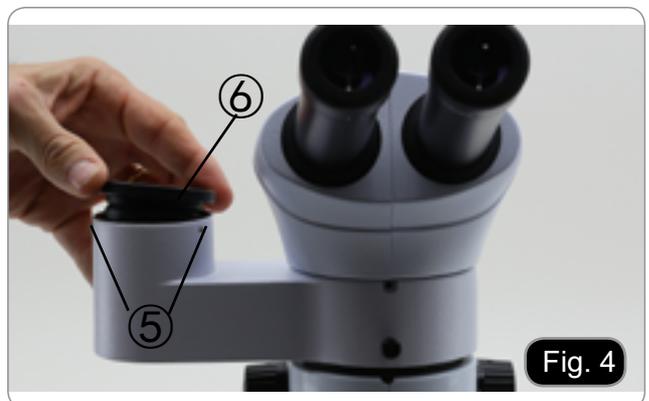

7. Install “C”-mount and camera in the empty hole (as explained in the camera instruction manual) and lock the fixing screw ⑦ . (Fig. 5)

8. Use of the attachment

8.1 Light path selection

1. Fully insert the light path selector ⑧ to split the light to binocular tube and to camera (50%/50%). (Fig. 6)
2. Pull out the selector to have 100% of the light to binocular tube, thus excluding light to trinocular port.

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.