Modular Series

INSTRUCTION MANUAL

Model
SZ-STLX

v 1.0 2017

SUMMARY

- 1. Warning**
 - 2. Intended use**
 - 3. Package Contents**
 - 4. Overview**
 - 5. Assembly**
 - 6. Using the base**
- Equipment disposal**

1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use.

We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users.

We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.


2. Intended use

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

3. Package Contents

DESCRIPTION	QUANTITY
Base	1
Vertical pillar	1
Drop prevention ring	1
Horizontal sliding arm	1
Head holder	1
Allen wrenches	2

4. Overview

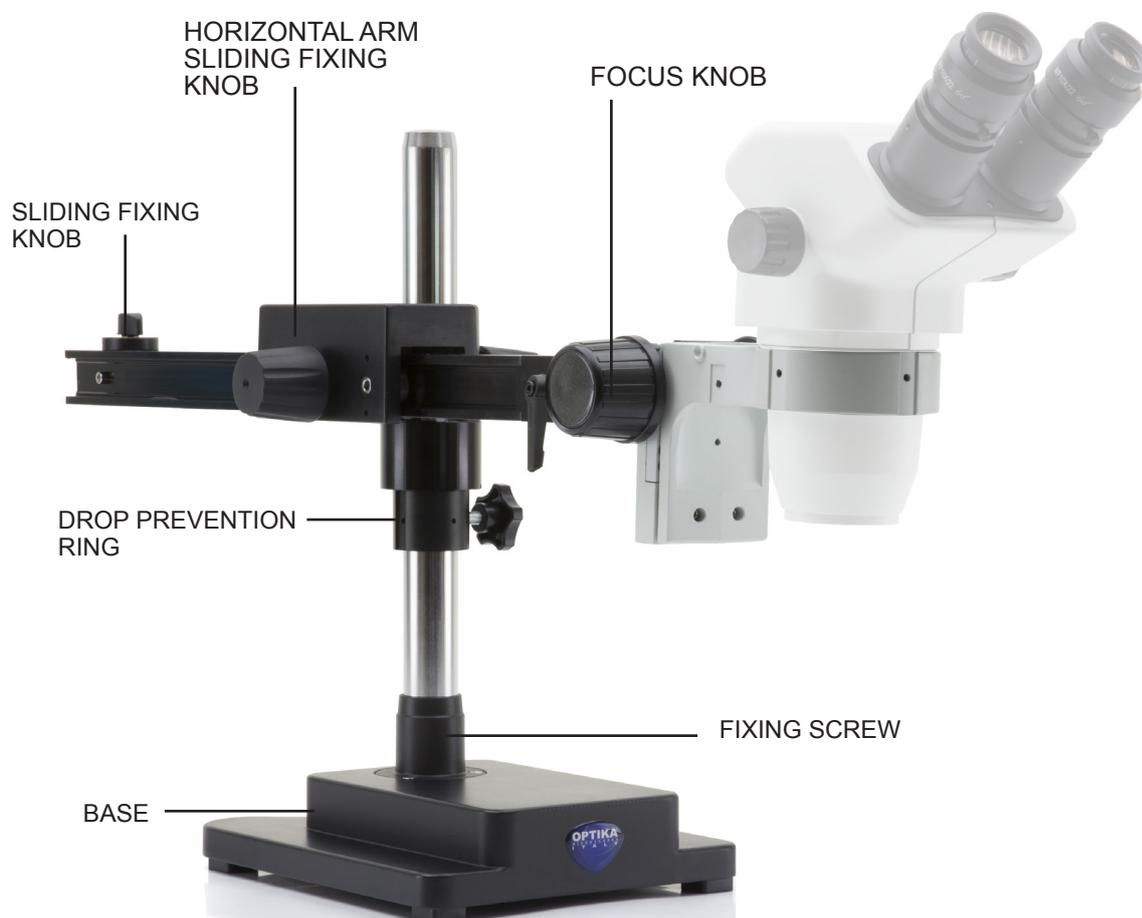5. Assembly

1. Screw the pillar on the base (Fig. 1)
2. Tighten the screw to lock the pillar (Fig. 2)
3. Insert the drop preventing ring. (Fig. 3)
4. Insert the horizontal arm and lock it with the locking knob (Fig. 4 & 5)

5. Install the head holder. Head holder can be installed in several ways according to specific need of the user.

Installation mode n° 1: (Fig. 6-7-8)

Insert the back part of the head holder (black round part) ① from above in the hole of the horizontal arm and screw the prevention knob ② from below.

- ▶ This mode allows the installation of heavy heads (or with heavy cameras installed on top of the head).

A system of stoppers (fig. 8) prevents the drop of the head.

Installation mode n° 2: (Fig. 9)

Insert the back part of the head holder (black round part) ① from below in the hole of the horizontal arm and screw the prevention knob ② from above.

- ▶ These two modes allow the tilting of the head (fig. 9).

Fig.9

Installation mode n° 3:

Insert the back part of the head holder (black round part) ① (Fig. 10) in the hole at the end of the horizontal arm and tighten the knob ③ (Fig. 11).

Fig.12

- ▶ This mode allows the swivel of the head (Fig. 12).

Install the head and lock the fixing knob (Fig. 13)

Fig.7

Fig.8

Fig.10

Fig.11

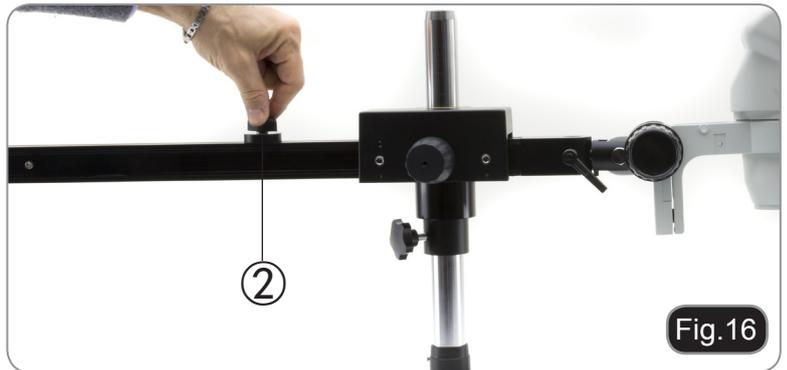
Fig.13

6. Using the base

Moving the horizontal arm (Fig. 14).
Unlock the knob on the left side of the horizontal arm ①.

The arm can be extended or shortened according to specific needs. (Fig. 15)

Setting the horizontal stopper (Fig. 16)
Unlock the fixing knob of the stopper ② and move it in a position suitable to user's needs.
Lock the fixing knob to set the movement limit.

Tilting of the head holder. (Fig. 17 - 18)
Loosen the allen wrench ③ on the top of the head holder.
The tilting black part can be moved. Once the desired angle is achieved, tighten the wrench again.

Swivel the head (Fig. 19 - 20)

Unlock the fixing knob ④ and rotate the head to the desired swivel angle (left or right), then fix the knob again.

Rotating the horizontal arm (Fig. 21)

Loosen the horizontal arm fixing knob ⑤ and rotate the microscope.


- ▶ **NOTE: 180° rotation of the microscope with respect to the base could cause a rollover of the entire system.**

Adjusting the tension of the focus knob (Fig. 22)

- ▶ **This adjustment allows to increase or decrease the tension of the knob by avoiding an involuntary descent of the microscope body under its own weight. Adjust the tension just above the point where the focus is stable.**

Grab the knobs ① with both hands and, holding right knob rotate the left one. Tension increases or decreases according the sense of rotation of the left focus knob.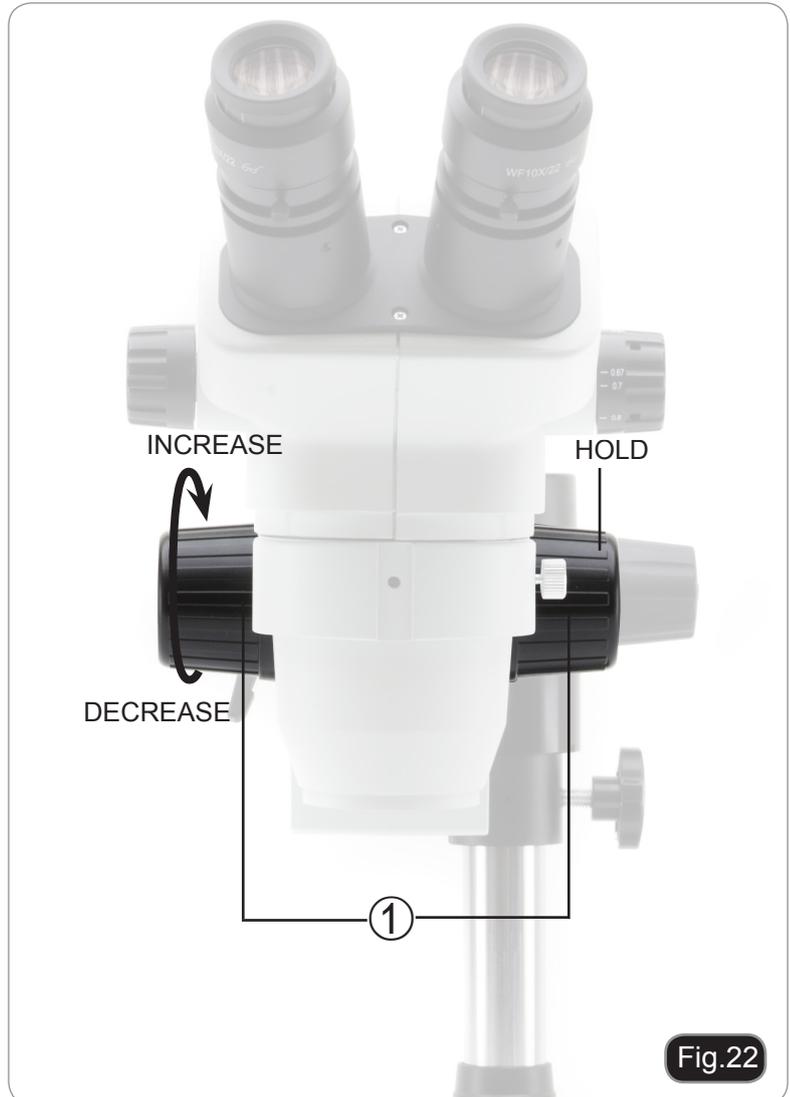

Equipment disposal

Art.13 Dlsg 25 july 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.