SZR Series

INSTRUCTION MANUAL

Model
SZR-180

Ver. 1.4 2026

Table of Contents

1.	Warning	3
2.	Safety Information	3
3.	Package content	4
4.	Unpacking	5
5.	Intended use	5
6.	Symbols and conventions	5
7.	Instrument description	6
8.	Assembling	8
9.	Using the microscope	10
9.1	Adjusting interpupillary distance	10
9.2	Diopter compensation	10
9.3	Magnification	11
9.4	Use of aperture diaphragm	11
9.5	Use of trinocular port	11
9.6	Use of transmitted light illumination	12
9.7	Use of conversion lens	13
10.	Microphotography	14
10.1	Installing the camera	14
10.2	Parfocality adjustment	14
11.	Maintenance	15
12.	Troubleshooting	16
	Equipment disposal	17

1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

2. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

3. Package content

- ① Base
- ② Microscope body
- ③ Objective
- ④ Observation head

- ⑤ Eyepieces
- ⑥ Power supply
- ⑦ Dust cover

4. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

5. Intended use

Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

6. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION


This symbol indicates a potential risk and alerts you to proceed with caution.

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

7. Instrument description

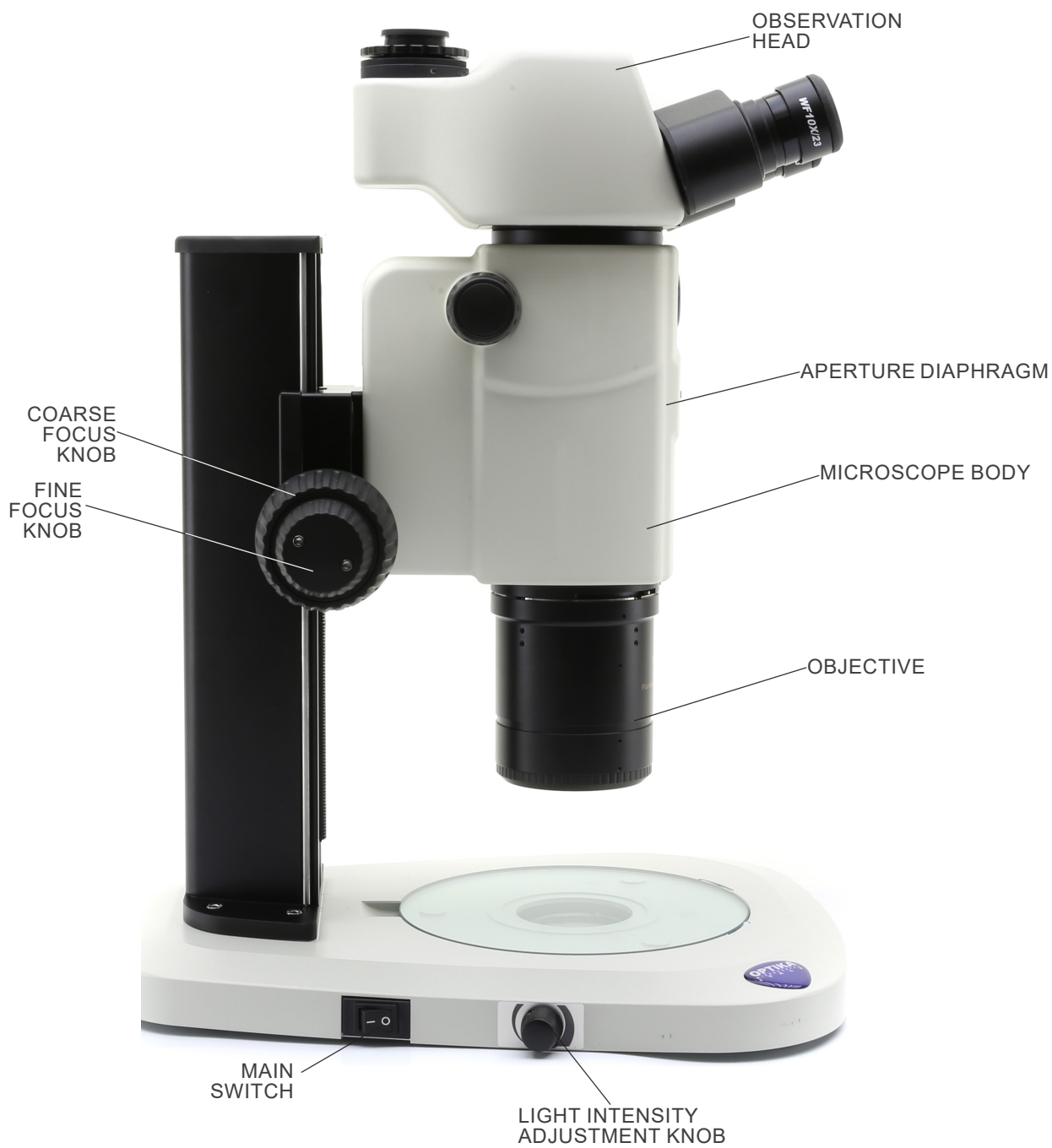
Opposite side


8. Assembling


1. Put the microscope base ① flat on the workbench. (Fig. 1)
2. Align the holes of the connecting plate ② of the stand column with the holes of the microscope base.
3. Use the provided fixing screws to connect the base and the column firmly.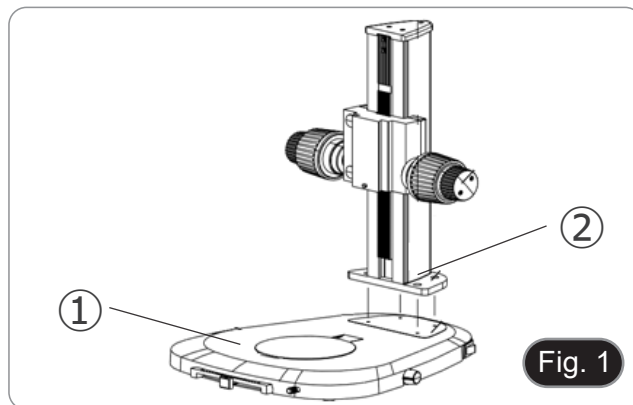


4. Insert the zoom body dovetail (3) into the dovetail of the focus system (4). (Fig. 2).
5. Lock the fixing screws (5).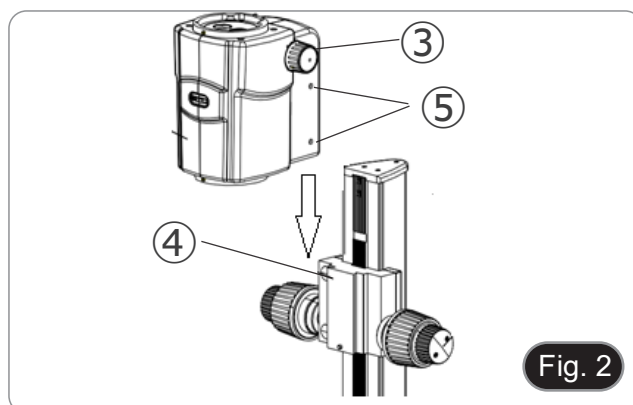


6. Install the head by aligning the round dovetail of the head with the hole of the microscope body. (Fig. 2)
7. Lock the fixing screw (6).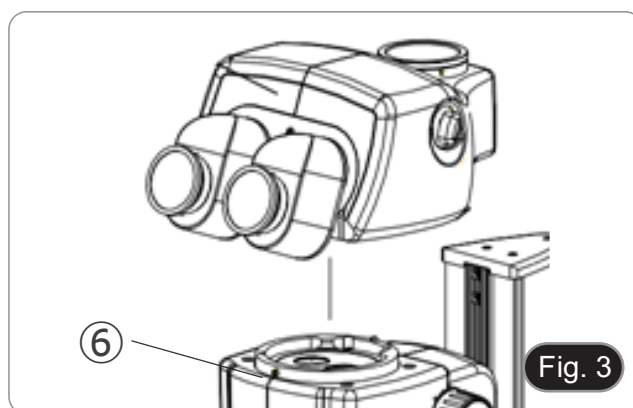

8. Insert eyepieces into empty eyepieces sleeves. (Fig. 4)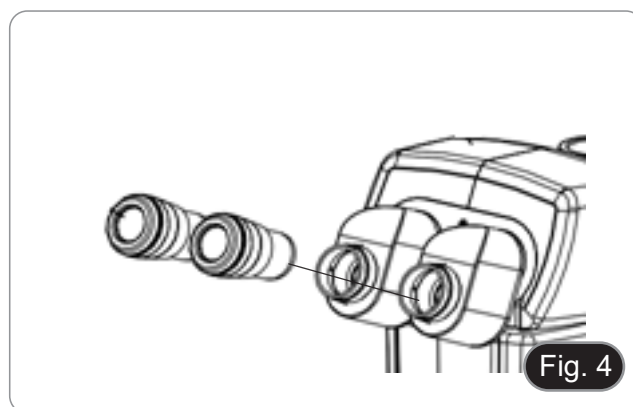

9. Screw the objective in the bottom part of the microscope body. (Fig. 5)
10. Lock the objective by screwing the fixing screw ⑦.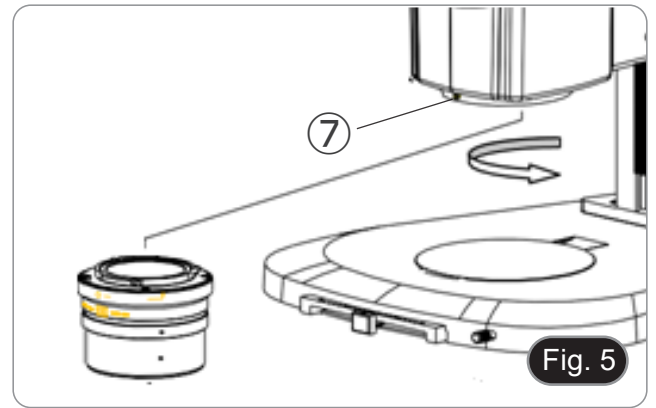

11. Plug the jack of the power supply in the socket of the microscope base.

9. Using the microscope


9.1 Adjusting interpupillary distance

1. Hold the right and left eyepiece tube with both hands and adjust the interpupillary distance by moving the two parts until one circle of light can be seen. (Fig. 6)
- If two circles appear, the interpupillary distance is too big.
- If two overlapped circles appear, the interpupillary distance is too small.

9.2 Diopter compensation

- **This compensation makes possible for people wearing glasses to adjust the microscope to their eyes and use the microscope without glasses.**
1. Put the zoom down to the lowest magnification ① and focus the specimen with the focusing knobs ②. (Fig. 7)
 2. Put the zoom to the maximum magnification and repeat the focusing.
 3. Return to the lowest magnification: the specimen will be out of focus.
 4. Adjust the diopter compensation ring of the right eyepiece ③ (Fig. 8) until the image of the right eyepiece is clear and sharp. Repeat the procedure for the left eyepiece ④.
 5. Check the focus of the image for the whole zoom range. It should be perfectly parfocal (focus is maintained during the change of magnification).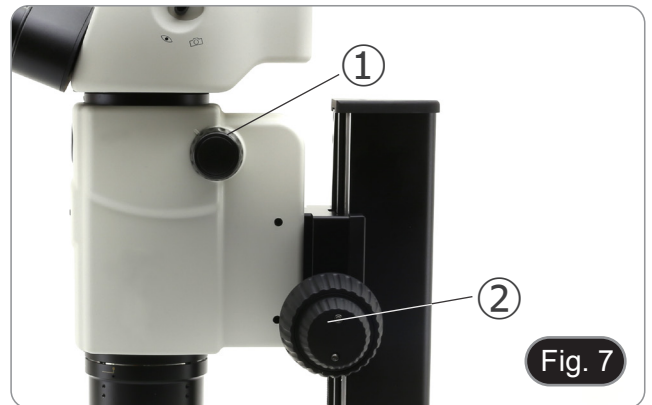

9.3 Magnification

1. Select the desired magnification by adjusting the zoom knob ①. (Fig. 9)
- Change the eyepieces if a different magnification range is needed.
 - The microscope body is equipped with a “click-stop” function that allows to obtain a precise setting of the desired magnification.
 - Total magnification used can be calculated as:
Eyepiece magnification * Zoom magnification * Objective lens magnification.

Eyepiece		10x	
Field number (mm)		24	
Objective	Total Mag	F.O.V (mm)	
1X	7.5X - 135X	30.67 - 1.7	

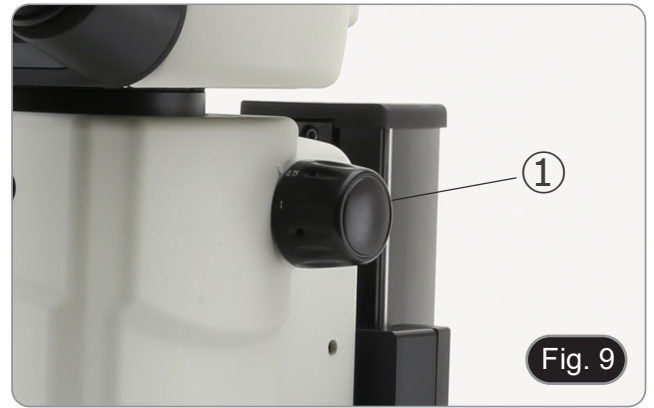
Fig. 9

9.4 Use of aperture diaphragm

By varying the diaphragm aperture, the user can adjust the image to achieve better contrast and greater depth of field. (Fig. 10)

1. Move the aperture diaphragm lever ② to the right or left. Rotation towards “○” opens the diaphragm while rotation towards “⊕” closes it.
 2. Adjust the diaphragm during observation to optimize the effect on contrast and depth of field.
- **Do not close the diaphragm too tightly so as not to have a poor image quality and not to have an excessively dark image.**

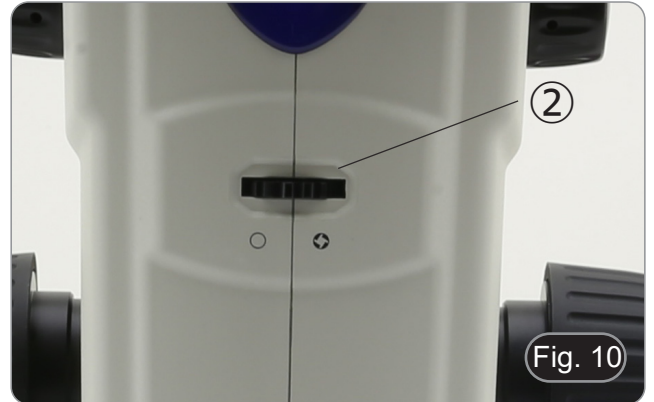
Fig. 10

9.5 Use of trinocular port

Observation head is equipped with a light split selector ③. (Fig. 11)

1. Moving the selector to the “eye” position ④ allows the binocular observation.
 2. Moving the selector to the “camera” position ⑤ allows the trinocular observation.
- **When the selector is in the “camera” position the right eyepiece will be completely dark. This is not a defect.**


Fig. 11

9.6 Use of transmitted light illumination


1. Set the main switch ① to "I" to turn on the base. (Fig. 12)
2. Turn the light intensity adjustment knob ② to adjust the brightness to the sample under observation.

3. Move the lever ③ located on the right side of the base. (Fig. 13)
 - This lever controls a black foil which slides horizontally along the light exit.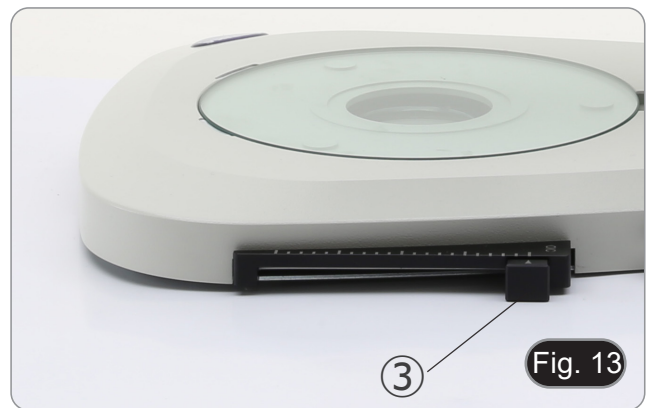

- Using the OIC (Oblique Illumination Contrast) system allows you to modulate the light to obtain three different types of observation.
4. When the foil is outside the optical path (Fig. 14), brightfield (BF) observation is obtained.

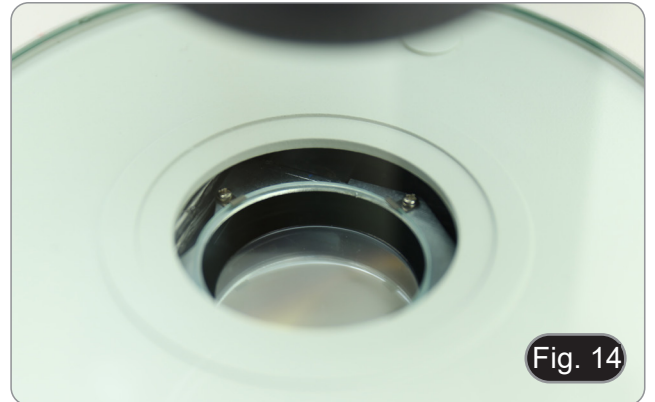
5. Moving the foil to an intermediate position (Fig. 15) results in the modulation of light called Oblique Illumination (OIC).

9.7 Use of conversion lens

- The conversion lens is used to optimize the light intensity according to the objective in use. (Fig. 16)
1. Fully insert the lever ① when using 1.6x or 2x objectives.
 2. Pull lever ① all the way out when using 0.5x or 1x objectives.
- **There are currently no additional objectives available other than the standard 1X objective.**

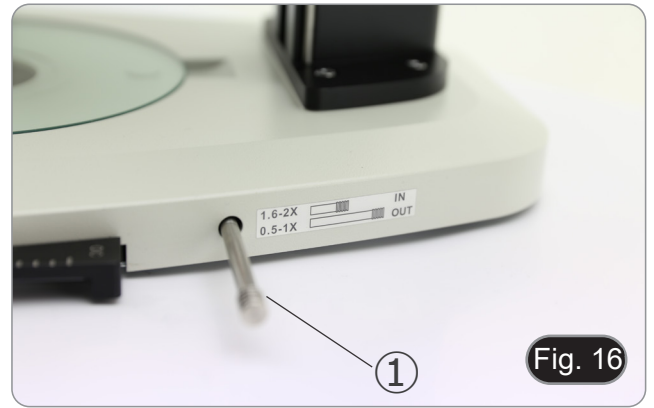10. Microphotography

10.1 Installing the camera

1. Screw the camera on the provided C-mount adapter.
2. To obtain a perfect alignment of the camera with respect to the eyepieces, loosen the fixing screw ① on the C-mount adapter, rotate the camera until correct alignment is achieved, then tighten the screw. (Fig. 17)

10.2 Parfocality adjustment

In order to have the same focus when observing the specimen through the eyepieces and on the screen/monitor, please check that the microscope is properly installed and follow the instructions below.

1. Use a low power magnification and focus the specimen.
2. Reach the highest magnification available using the zoom knob and then focus the specimen again.
3. Turn on the live-view on the camera, without changing the focus on the microscope.
4. While observing the image on the screen/monitor, adjust the focus by turning the knurled knob ② on the C-mount adapter. (Fig. 18)

11. Maintenance

Operative environment

This microscope is recommended to be used in a clean, dry and shock free environment with a temperature of 5°-40°C and a maximum relative humidity of 75 % (non condensing). Use a dehumidifier if needed.

To think about when and after using the microscope

- The microscope should always be kept vertically when moving it and be careful so that no moving parts, such as the eyepieces, fall out.
- Never mishandle or impose unnecessary force on the microscope.
- Never attempt to service the microscope yourself.
- After use, turn off the light immediately, cover the microscope with the provided dust-cover, and keep it in a dry and clean place.

Electrical safety precautions

- Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off-position.
- Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Cleaning the optics

- If the optical parts need to be cleaned try first to: use compressed air.
- If that is not sufficient: use a soft lint-free piece of cloth with water and a mild detergent.
- And as a final option: use the piece of cloth moistened with a 3:7 mixture of ethanol and ether.
- **Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.**
- Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics.
- Do not disassemble objectives or eyepieces in attempt to clean them.

For the best results, use the OPTIKA cleaning kit (see catalogue).

If you need to send the microscope to Optika for maintenance, please use the original packaging.


12. Troubleshooting

Review the information in the table below to troubleshoot operating problems.

PROBLEM	CAUSE	SOLUTION
I. Optical Section:		
The illumination is ON, but the field of view is dark	The plug is not connected to the illumination	Connect the cable
	The brightness is too low	Adjust to a proper setting
The edge of the field of view is vignetted or the brightness is asymmetric	The incident illuminator is not correctly oriented	Change the angle of the incident illuminator
Dust and stains can be seen in the field of view	There are stains and dust on the specimen	Clean the specimen
	There are stains and dust on the eyepiece	Clean the eyepiece
Poor image quality: <ul style="list-style-type: none"> The image is not sharp The contrast is not high The details are not clear Image glares 	The lenses (additional lens, objective, eyepieces) are dirt	Thoroughly clean all the optical system
One side of the image is out of focus.	The specimen is out of place (tilted)	Place the specimen flat on the stage
II. Mechanical Section:		
The focus knob is hard to turn	The tension adjustment collar is too tight	Loosen the tension adjustment collar
The focus is unstable	The tension adjustment collar is too loose	Tighten the tension adjustment collar
III. Electric Section:		
The LED doesn't turn on	No power supply	Check the power cord connection
The brightness is not enough	The brightness adjustment is low	Adjust the brightness
The light blinks	The power cord is poorly connected	Check the power cord
IV. Viewing tube assembly:		
The field of view of the two eyes is different	The interpupillary distance is not correct	Adjust the interpupillary distance
	The diopter correction is not right	Adjust the diopter correction
	The viewing technique is not correct, and the operator is straining the eyesight	When look into the eyepieces, do not stare at the specimen but look at the whole field of view. Periodically, move the eyes away to look at a distant object, then back into the eyepieces
V. Microphotography and video:		
The image is unfocused	Incorrect focusing	Adjusting the focus system as in the present manual
Bright patches appear on the image	Stray light is entering the microscope through the eyepieces and through the camera viewfinder	Cover the eyepieces and the viewfinder with a dark cloth

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

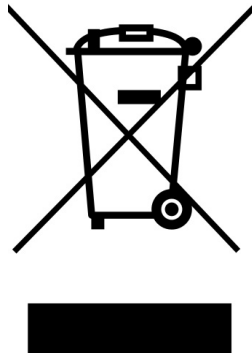The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.