“All-in-one” series

# INSTRUCTION MANUAL

Model
TB-6PRO4K

Ver. 1.1    2024

---

## Table of contents

1.	Warning	3
2.	Safety Information	3
3.	Package content	4
4.	Unpacking	5
5.	Intended use	5
6.	Symbols and conventions	5
7.	Instrument description	6
8.	Assembling	8
8.1	Camera assembly	8
8.2	Mini-PC assembly	8
8.3	Cable connection	8
8.4	Microscope assembly	9
8.5	Parfocality adjustment	10
9.	Use of camera	11
10.	Micrometric Slide M-005	11
	Equipment disposal	12

---

## 1. Warning

This camera is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

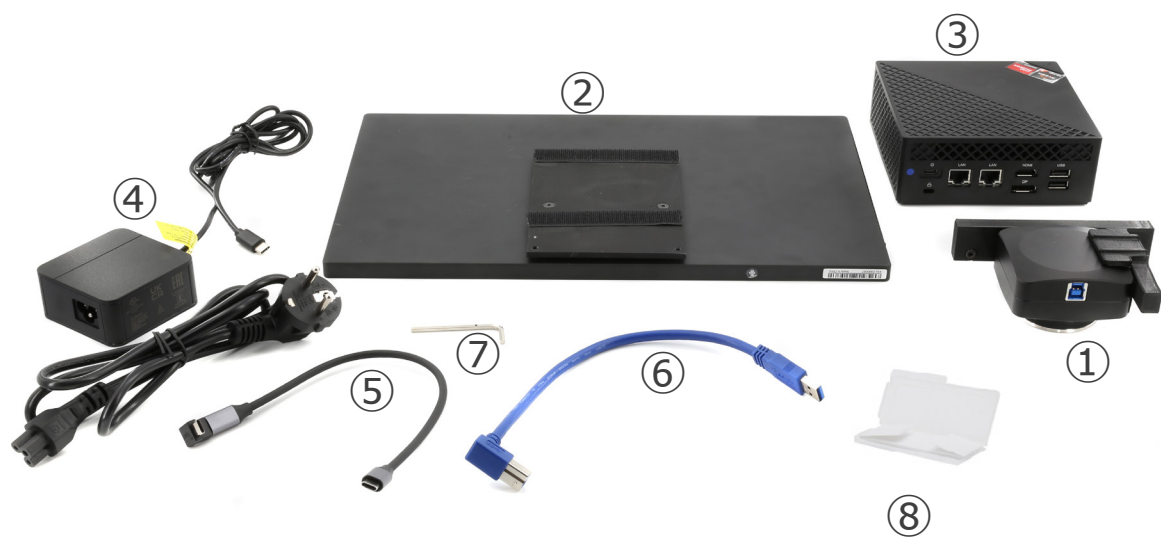
## 2. Safety Information

### Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

### 3. Package content

- ① Camera with mounting bracket
- ② Monitor
- ③ Mini-PC
- ④ Power supply + power cord for mini-PC

- ⑤ USB-C cable
- ⑥ USB3.0 cable
- ⑦ 3mm Allen wrench
- ⑧ Micrometric slide

---

## 4. Unpacking

The device is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items fall out and get damaged.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

## 5. Intended use

### Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

### IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

## 6. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

### CAUTION

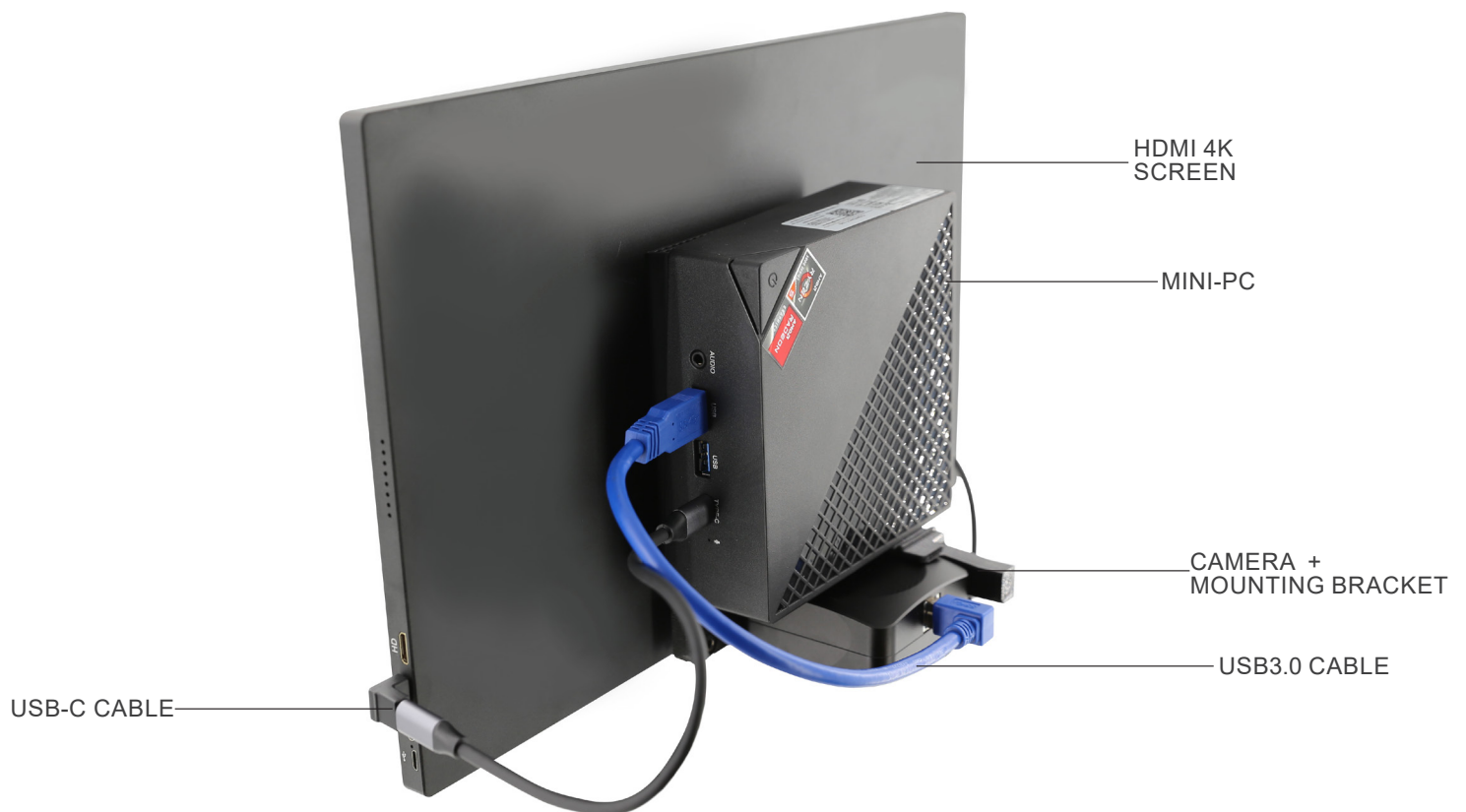
This symbol indicates a potential risk and alerts you to proceed with caution.

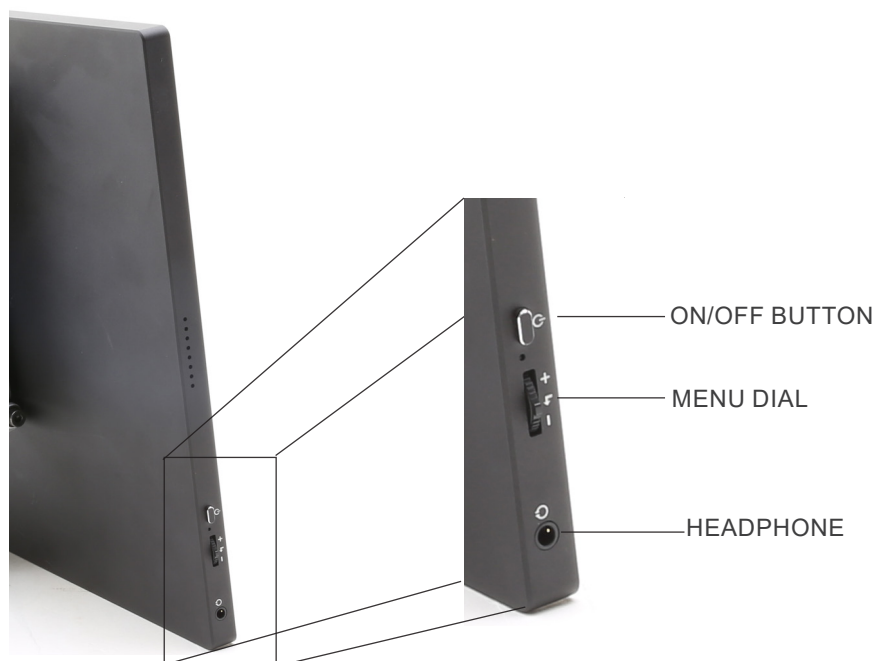
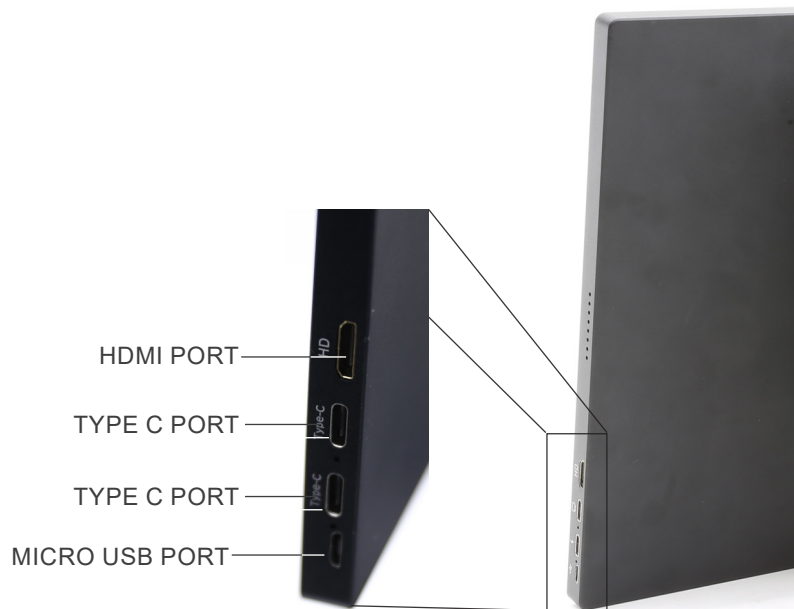
### ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

## 7. Instrument description

## 8. Assembling

### 8.1 Camera assembly

- **USB camera is delivered with the mounting bracket already assembled.**
1. Attach the mounting bracket to the monitor, using the supplied screws. (Fig. 1)

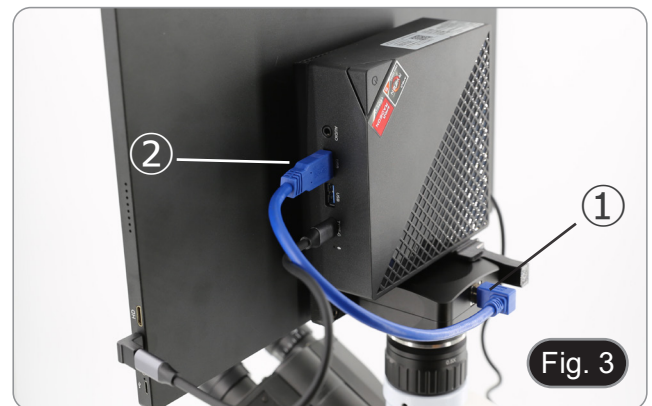
### 8.2 Mini-PC assembly

1. Using the supplied Velcro adhesive, attach the mini PC to the back of the monitor. (Fig. 2)
- For better system stability, it is recommended to rest the mini-PC on the mounting bracket.

### 8.3 Cable connection

1. Connect the USB3.0 cable from the camera ① to one of the USB3.0 ports on the mini PC ②. (Fig. 3)


2. Connect the USB-C cable ③ in the USB-C port of the mini-PC. (Fig. 4)
  3. Connect the other end of the cable to the upper USB-C port on the monitor ④ to power the monitor and to activate the monitor's "touch" mode. (Fig. 4)
- The monitor is equipped with "touch screen" functionality. Connecting the USB cable from the mini-PC to the monitor allows the operator to be able to work normally using all the functions of the PC by simply touching the icons on the monitor.
  - However, a keyboard and mouse (not included) can also be connected to the mini-PC.

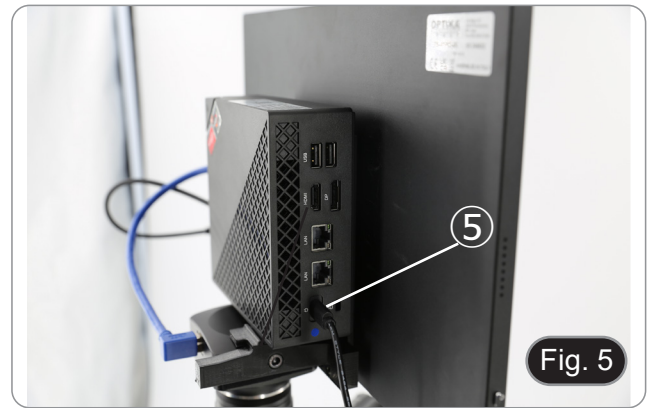4. Connect the USB-C plug of the mini-PC power supply in the USB-C port on mini-PC ⑤ to power it. (Fig. 5)
5. Connect the power cord to the power supply.
6. Plug the power cord into the wall socket.


#### 8.4 Microscope assembly

1. Remove the protective cap from the camera and screw on the C-mount. (Fig. 6)

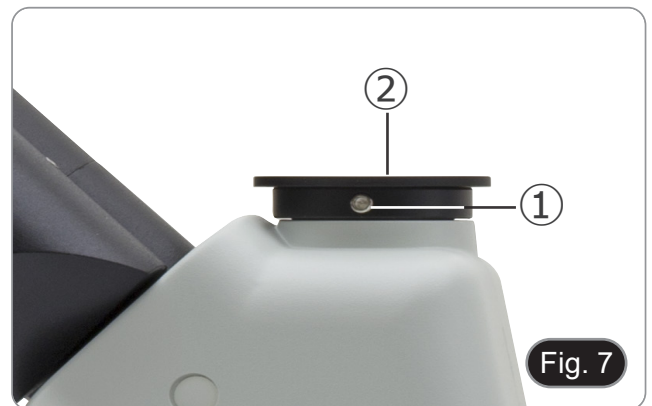2. Remove the protective cap ② from the third port of the observation head by loosening the fixing screw ① located on the side of the third port. (Fig. 7)

3. Insert the dovetail of the C-mount into the third port, properly align the whole system with respect to the microscope, and then tighten the fixing screw ①. (Fig. 8)

## 8.5 Parfocality adjustment

In order to have the same focus when observing the specimen through the eyepieces and on the monitor, please check that the microscope is properly set and follow the instructions below.

1. Use a low power objective and focus the specimen using the microscope focus knobs.
2. Switch to the highest dry objective available on the microscope (40x or 60x) and focus the specimen again.
3. Turn on the live-view on the camera, without changing the focus on the microscope.
4. While observing the image on the monitor, adjust the focus by turning the knurled knob on the C-mount adapter. (Fig. 9)

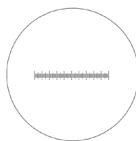
## 9. Use of camera

1. Tap the icon for the ProView software (or double-click the icon with the mouse). The software starts.
  2. In the "Camera List" panel, the "C-P6" item is shown.
  3. Tap on the C-P6 item (or click with the mouse): the live image will be displayed on the main window of the software.
  4. Adjust the camera parameters by acting on the exposure time ("Exposure and Gain" panel) and the white balance ("White Balance" panel).
  5. Once the first adjustments have been made, you can operate normally.
- The software's User Manual is available in PDF format within the application itself and can be opened using the "F1" function key. The manual contains all the operating instructions for using the camera and for the various functions of the software.
  - You must have Acrobat Reader installed to view the manual.


## 10. Micrometric Slide M-005

**Micrometric slide, 26x76mm, with 2 scales**  
**(1mm/100div. for biological microscopes / 10mm/100div. for stereomicroscopes)**

1 DIV=0.01mm

**For biological microscopes calibration**

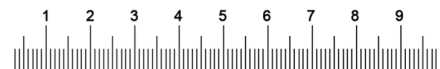
1 DIV=0.1mm

**For stereo microscopes calibration**

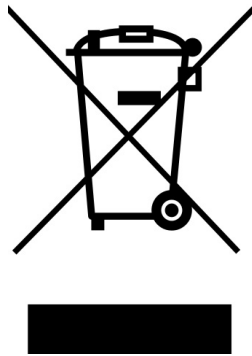---

## Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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**OPTIKA® S.r.l.**