OPTIStar

By **OPTIKA**

OPTISTAR By OPTIKA Series

INSTRUCTION MANUAL

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Ver. 1.0 2021

Table of contents

1.	Warning	3
2.	Safety Information	3
3.	Unpacking	3
4.	Intended use	3
5.	Symbols and conventions	3
6.	Instrument description	4
6.1	USB Eyepiece camera 2MP / 5MP	4
6.2	USB/HDMI C-mount camera 21MP	4
7.	Assembling	5
7.1	USB Eyepiece camera 2MP / 5M	5
7.1.1	Monocular or binocular model	5
7.1.2	Trinocular model	5
7.2	USB/HDMI C-mount camera 21MP	5
7.2.1	Use with PC	5
7.2.2	Use with HDMI monitor	6
7.3	Parfocality adjustment	6
7.3.1	OS-B Series	6
7.3.2	OS-C Series	6
7.3.3	OS-SZ Series	6
8.	Use of camera	7
8.1	USB Eyepiece camera 2MP / 5MP	7
8.1.1	Operating Instructions	7
8.1.2	Detailed Operating Instructions	10
a)	Menu Bar	10
b)	Tool Bar	26
8.1.3	Sidebar	27
8.2	USB/HDMI C-mount camera 21MP	30
8.2.1	Use of camera in HDMI mode	30
a)	Language selection	30
b)	Line setting	30
c)	Exposure setting	32
d)	White Balance setting	33
e)	Color setting	33
f)	Light frequency setting	33
g)	Formatting the SD card	34
h)	Date and Time setting	34
i)	Image size setting (size of image in recording)	34
j)	Mirror image setting	35
k)	Date Stamp setting	35
l)	System setting	35
m)	Electronic enlargement and reduction	36
n)	Switch camera / image mode	36
o)	Remote control and key explanation	36
8.2.2	Use of camera in USB mode	37
a)	User Interface	37
b)	Camera Control Panel	38
c)	Capture Images	41
d)	Measurement	42
e)	Calibration	43
	Equipment disposal	44

1. Warning

This camera is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

2. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

3. Unpacking

The device is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items fall out and get damaged.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

4. Intended use

Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

5. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

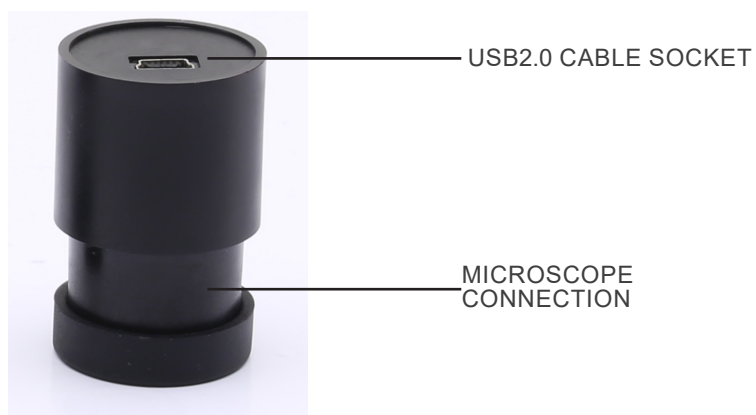
ELECTRICAL SHOCK


This symbol indicates a risk of electrical shock.

6. Instrument description

6.1 USB Eyepiece camera 2MP / 5MP

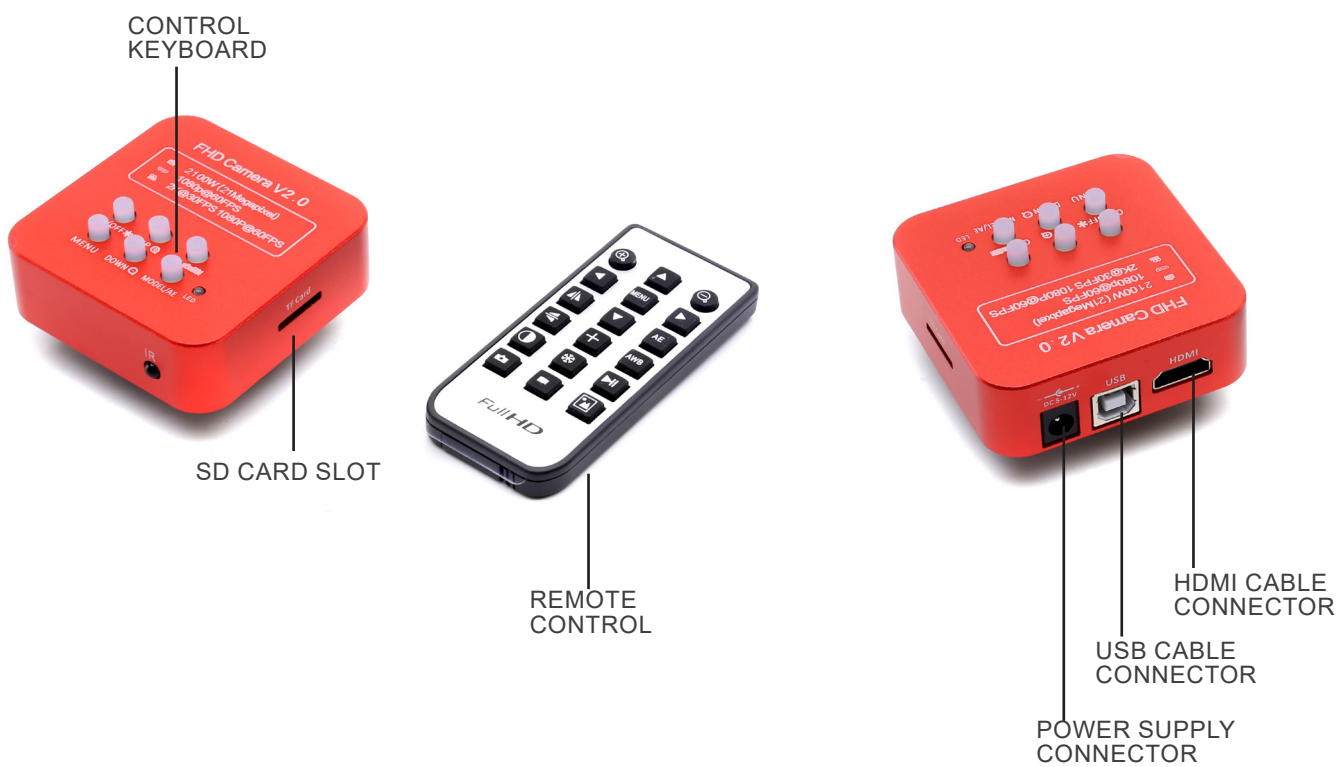6.2 USB/HDMI C-mount camera 21MP


7. Assembling

7.1 USB Eyepiece camera 2MP / 5MP

7.1.1 Monocular or binocular model

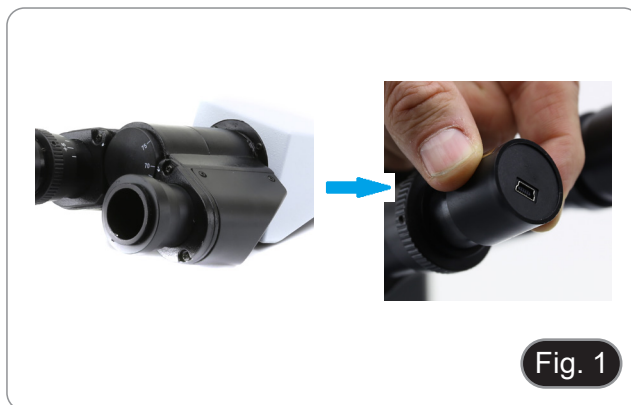1. Remove the eyepiece from the observation head and insert the camera into the empty hole of the observation head. (Fig. 1)
2. Connect the USB cable.
3. Install the software.


7.1.2 Trinocular model

1. Insert the camera into the empty hole of the photo tube. (Fig. 2)
2. Connect the USB cable.
3. Install the software.

7.2 USB/HDMI C-mount camera 21MP

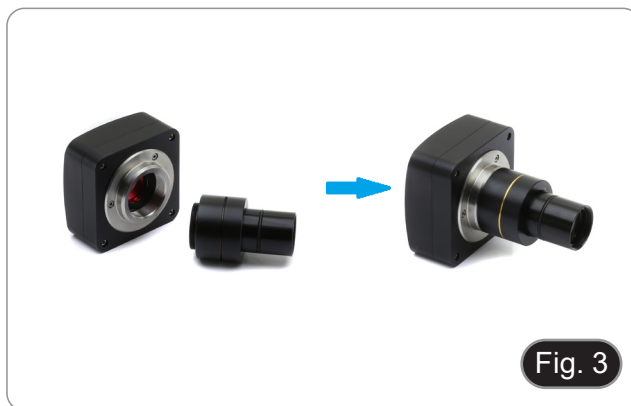1. Remove the dust cover from camera and from projection lens.
2. Screw the projection lens to the camera's thread. (Fig. 3)

3. Insert the final part of the projection lens into the empty hole of the photo tube. (Fig. 4)

7.2.1 Use with PC

1. Connect the USB cable.
2. Install the software.

7.2.2 Use with HDMI monitor

1. Insert the SD card into its slot. (Fig. 5)
2. Connect the HDMI cable.
3. Connect the power supply.

7.3 Parfocality adjustment

The proper parfocality adjustment is obtained when the same focus is reached when looking into the eyepieces and on the screen/monitor.

7.3.1 OS-B Series

1. Use a low power objective and focus the specimen.
2. Switch to the highest dry objective available on the microscope (40x or 60x) and focus the specimen again.
3. Turn on the live-view on the camera, without changing the focus on the microscope.
4. Rotate the knurled ring ① (Fig. 6) and while observing the image on the monitor, adjust the focus.

7.3.2 OS-C Series

1. Use a low power objective and focus the specimen.
2. Switch to the highest dry objective available on the microscope (40x or 60x) and focus the specimen again.
3. Turn on the live-view on the camera, without changing the focus on the microscope.
4. Unlock the knob ② (Fig. 7) and while observing the image on the monitor, adjust the focus by moving up or down the camera. (Fig. 8)
5. Lock the fixing knob ②.

7.3.3 OS-SZ Series

1. Use a low power magnification and focus the specimen.
2. Reach the highest magnification available using the zoom knob and then focus the specimen again.
3. Turn on the live-view on the camera, without changing the focus on the microscope.
4. Unlock the knob ② (Fig. 7) and while observing the image on the monitor, adjust the focus by moving up or down the camera. (Fig. 8)
5. Lock the knob ②.

8. Use of camera

8.1 USB Eyepiece camera 2MP / 5MP

These models are driven by IMAGE STAR (either for Windows and for IOS).

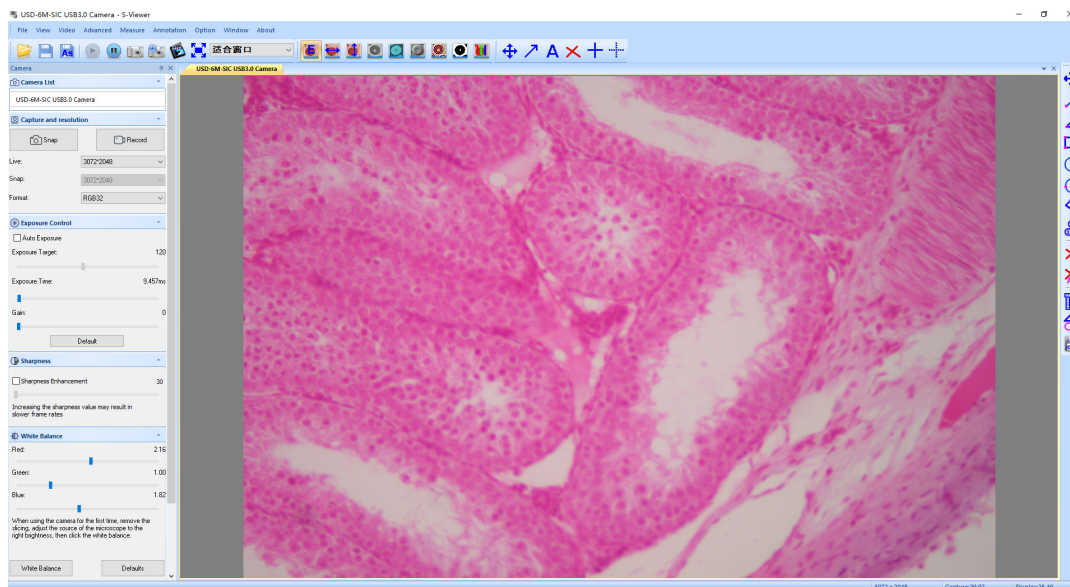
- The software can be downloaded from this link: http://www.optikamicroscopes.com/OPTIStar/IMAGE_STAR.zip
- **NOTE: no driver installation is required for the cameras. The software setup procedure automatically installs all the needed drivers for the correct operation of the camera.**

1. Run IMAGE STAR by clicking the icon  on the desktop.
2. Click on the camera name in the section *Camera List* on the upper left corner of the screen.
3. Live image starts automatically.


8.1.1 Operating Instructions

Basic operation on Live image


1. The first time the software is started after installation, an image capture device will be found for image preview.

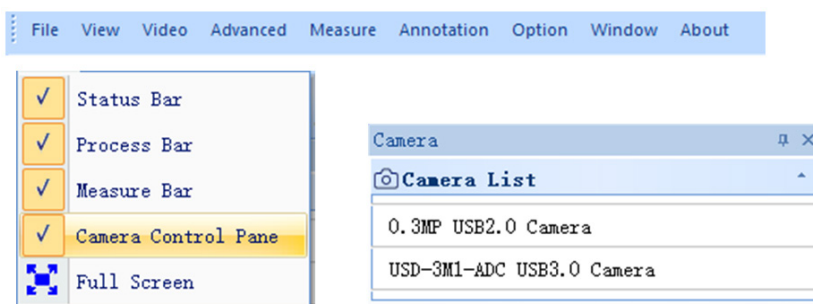
2. Select the desired device.
When more than one imaging device appears on computer, this software allows you to select the desired imaging device. Select *Menu --> View --> Camera Control Panel*. Pop up the left sidebar to select the target imaging device in the device name list.

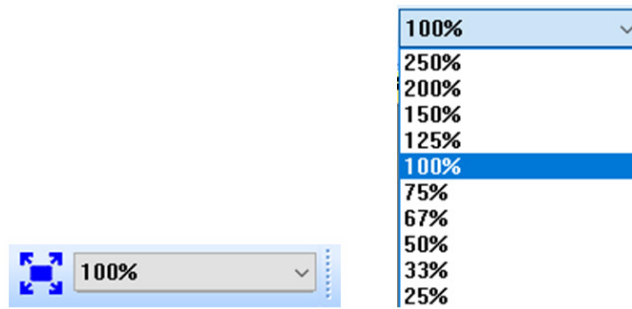3. Set image previewing size.
In the toolbar below the menu, change the image preview size by selecting full screen or by selecting the option in the output size list.

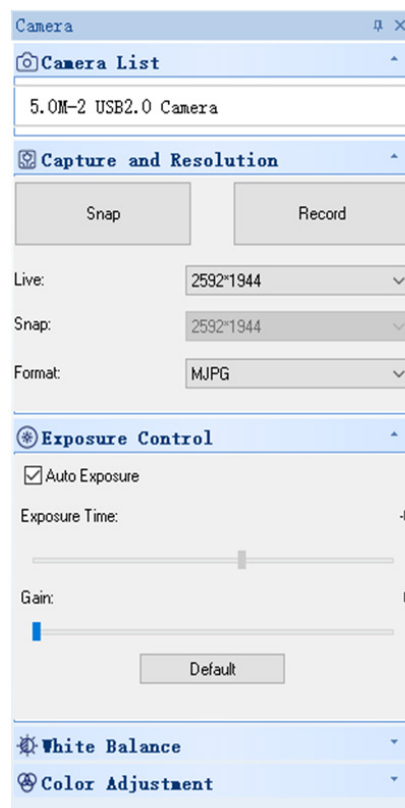4. Set image attribute.
In the left sidebar, adjust the basic parameters of the image. (See 8.3 Sidebar for details)

5. Capture the image.

Save: select *Menu --> File --> Save*, it will save automatically according to the information in “set to be saved”.

Save as: select *Menu-->File-->Save As*, then the dialog box of “Save As” pops up.

Save Option [X]

File Name

Prefix
Image

Start Suffix
69

Folder
C:\Users\tstx06\Documents\Swift-EasyView ...

Image Type

Bitmap Jpeg

Save With

Title Data/Time

Save With(Still Measurement)

Measure Ruler

Description

Title
[]

Author
[]

OK Cancel

8.1.2 Detailed Operating Instructions

a) Menu Bar

File

1.  **O**pen... **Ctrl+O**

Load images from local disk, and displays them in windows; operations like dynamic measurement can also be performed.

2. **C**lose

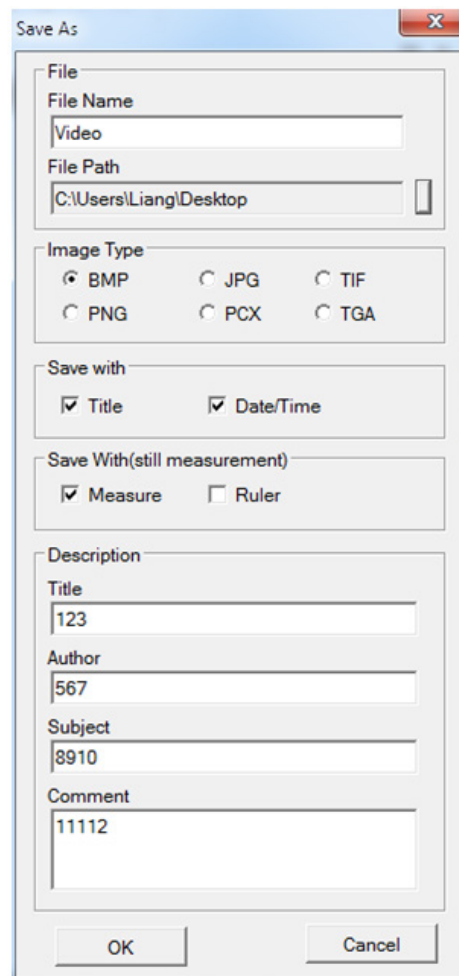
Closes active live or image windows.

3.  **S**ave **Ctrl+S**

Automatically saves the current live image or saves current images in the default path.

4. **As** Save **As...**

See the dialog box of “Save As” below.

5. **Print...** **Ctrl+P**

Print current image.

6. **Print Preview**

Preview the image of the printed image on A4 paper.

7. **Print Setup...**

Some settings for printing.

8. **Select Source...**

Unused with this kind of camera.

9. **Acquire...**

Unused with this kind of camera.

10. List of recently opened files.

11. **Exit**

Quit the application. (Close the software)

View

1. **S**tatus Bar

Show or hide the status bar.

2. **P**rocess Bar

Show or hide the process bar.

3. **M**easure Bar

Show or hide the measuring tool bar.

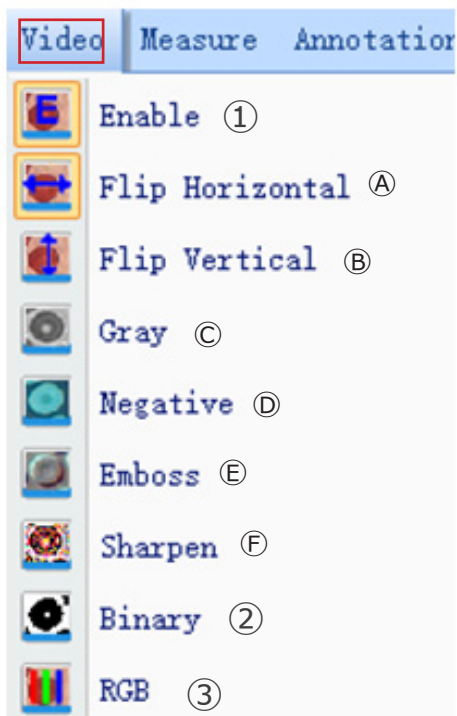
4. **C**amera Control Pane

Show or hide the camera control panel.

5. **F**ull Screen


Show image in full screen.

Video

1. Enable

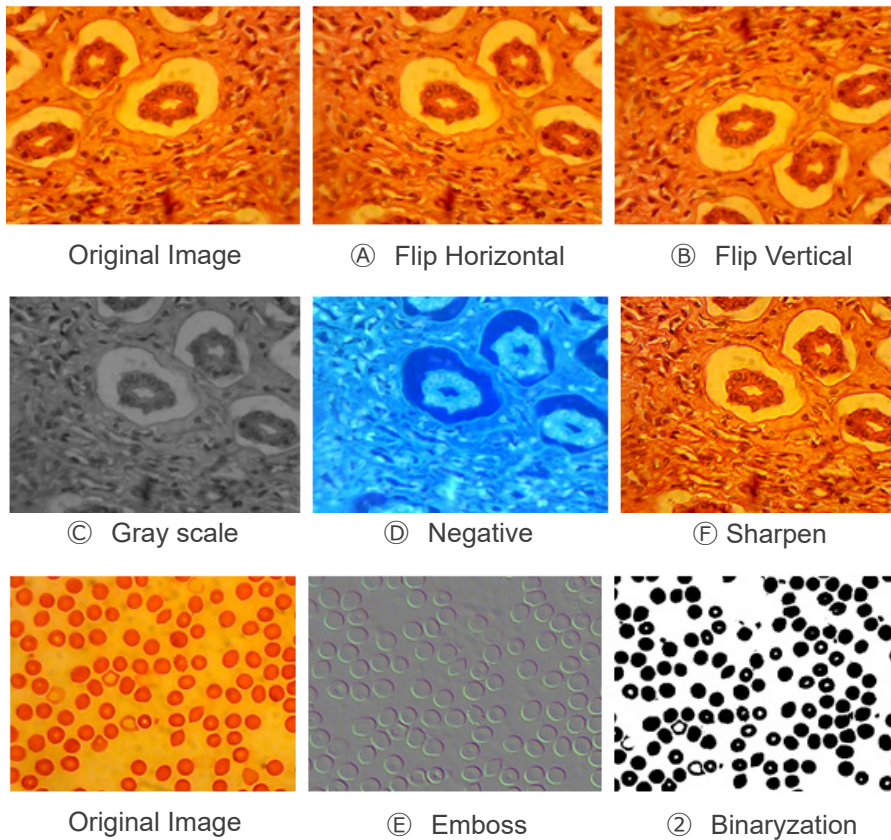Activate the opening or closing real-time processing or real-time measurement.
All of the effects of processing are as follows.

2.  Binary

If image binarization is needed, user can adjust the threshold value of binarization (1 to 255) to display dynamically. Start using check box to use or close the binarization real-time processing effect.

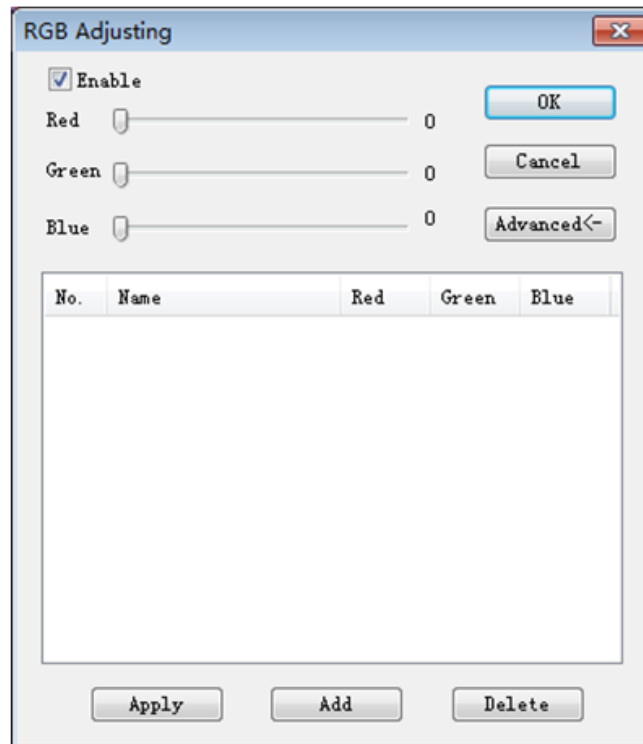
3.  Rgb

Start or close RGB real-time adjusting on image.

Enable : Start or close RGB adjustment

: Display advanced settings panel


Pitch on check box, adjust the red-green-blue sliders for image tricolor match.


Apply: Select options from list box and then apply them to the current image.

Add: Click "Add", input option names, add the adjusted red-green-blue value to the database, and it will display in the list.


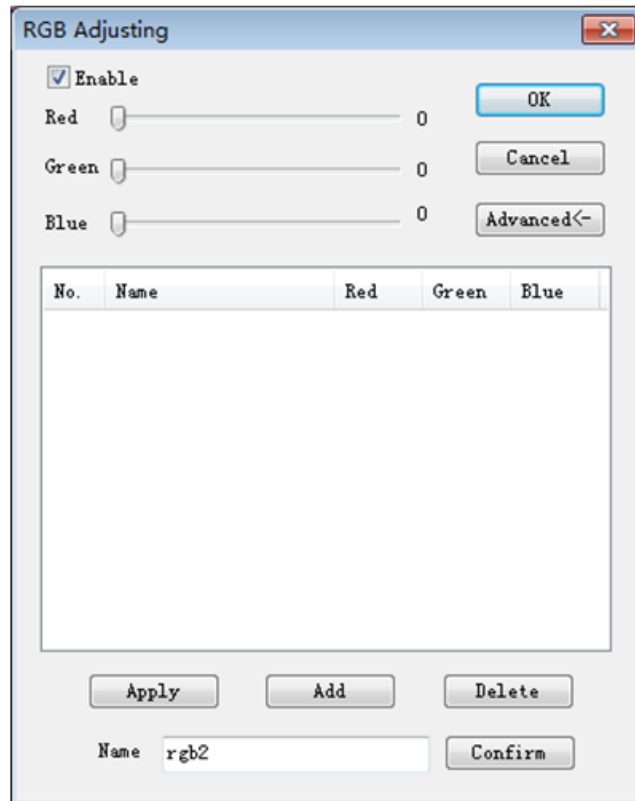
Delete: Select options from list box, then delete options from the list or database.

Advanced

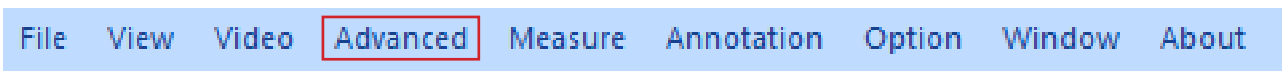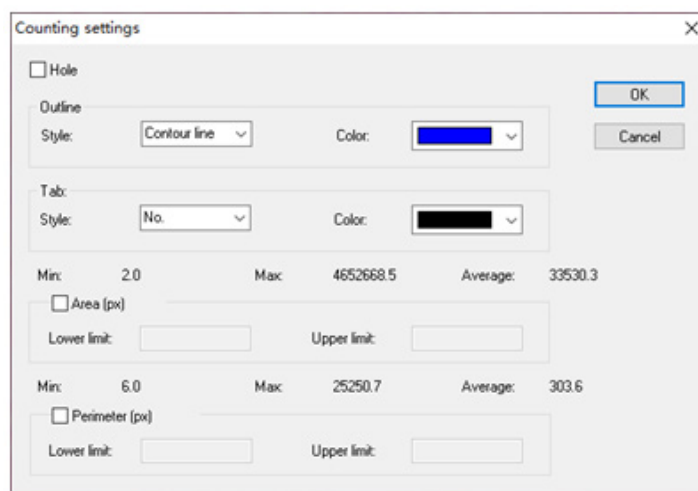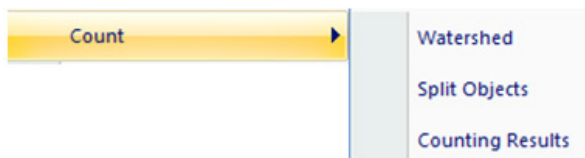1) Count

Watershed:

watershed algorithm is more suitable for the image with single background or large difference between the target and the background. The image with complex background is not suitable.

Holes:

when selecting holes, app will mark the outer contour of the segmentation object with contour color, and also mark the holes in the object with contour color and output its parameters (holes are defined as pixels whose gray or color is close to the background in the target object).

Outline:

use the outline to mark the segmentation object, and the color can be determined by the user.

Style: there are three options for the style.

Contour line: use the outline to mark the segmentation object, and the color can be determined by the user;

Fill: fill the divided objects with custom colors;

Nothing: do not mark the contour of the segmented object.

Color:

customize the color used for marking.

Tab:

an information will be appended to the measured object.

Style: there are three options for the style.

No: no informations will be appended to the measured object;

Area: information of area will be displayed;

Perimeter: information of perimeter will be displayed.

Color:

customize the color used for marking.

Area (px):

Lower limit: set the lower limit of the area of the object to be segmented, that is, objects smaller than the lower limit will not be divided or counted.

Upper limit: set the upper limit of the area of the object to be segmented, that is, objects larger than the upper limit will not be segmented and counted.

Perimeter (px):

Lower limit: set the lower limit of the perimeter of the object to be segmented, that is, objects smaller than the lower limit will not be segmented or counted.

Upper limit: set the upper limit of the perimeter of the object to be segmented, that is, objects larger than the upper limit will not be segmented and counted.

Split Object

Segmentation: when two single objects are counted into one object, these can be separated by segmentation and counted again.


Counting results

Statistical results: Displays statistics and exports them.

No.	core	Area	Perimeter
1	(1117.0,2046.0)	3.0	8.0
2	(260.0,2046.0)	2.0	6.0
3	(374.0,2043.0)	111.0	59.8
4	(274.0,2041.0)	42.0	41.3
5	(1930.0,2040.0)	192.5	59.2
6	(1524.0,2031.0)	38.5	24.2
7	(1548.0,2035.0)	364.0	82.3
8	(322.0,2037.0)	641.0	214.6
9	(426.0,2018.0)	61.5	29.9
10	(387.0,2010.0)	56.0	29.3
11	(369.0,2012.0)	177.0	61.1
12	(392.0,2000.0)	8.5	11.4
13	(918.0,1999.0)	82.5	43.6
14	(380.0,1984.0)	387.5	114.3
15	(925.0,1975.0)	145.5	54.0
16	(381.0,1958.0)	8.5	11.4
17	(919.0,1956.0)	143.5	58.4
18	(391.0,1959.0)	88.5	48.4
19	(916.0,1920.0)	24.0	19.7
20	(893.0,1899.0)	6.0	10.0


2) *Stitching*

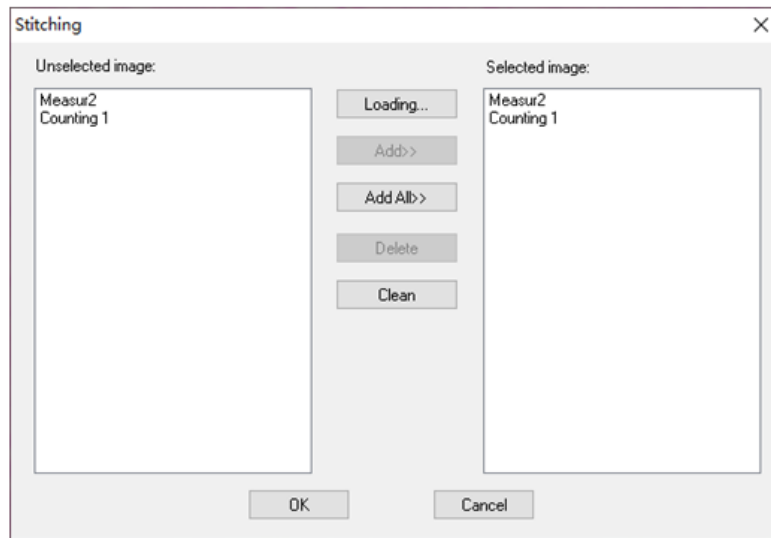Click this function to load the opened image to the list on the left.
Click *Loading...* to import the image to be stitched from outside. Select the image to be stitched, click *Add* to import the list on the right, and click *OK* to splice the selected image.

3) *EDF*

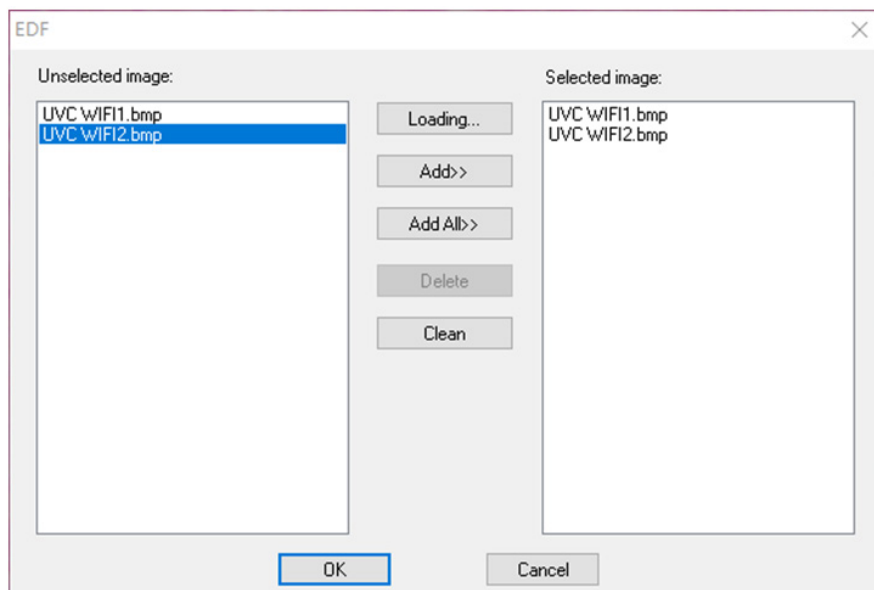
Click this function to load the opened image to the left list.
Click *Loading...* to import EDF image from outside. Select the image that needs EDF, click *Add* to import the right list, and click *OK* to get the result image.

4) HDR

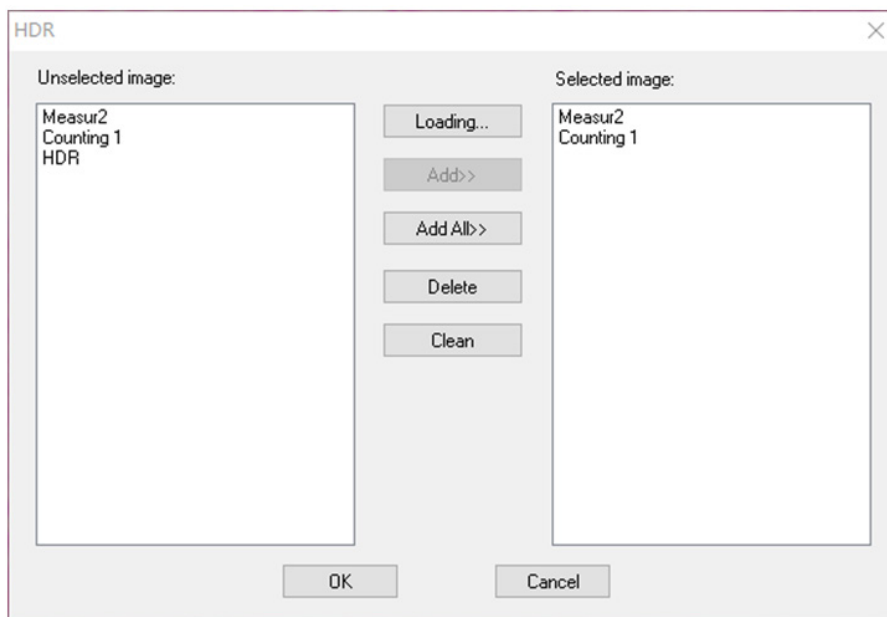
Turn off the camera auto exposure, dim the light source to the appropriate brightness, click *Start HDR*, and then slowly adjust the brightness of the light source.

Click this function to load the opened image to the left list.
Click *Loading...* to import HDR image from outside. Select the image that needs HDR, click *Add* to import the right list, and click *OK* to synthesize the selected image by HDR.


5) *MIX*

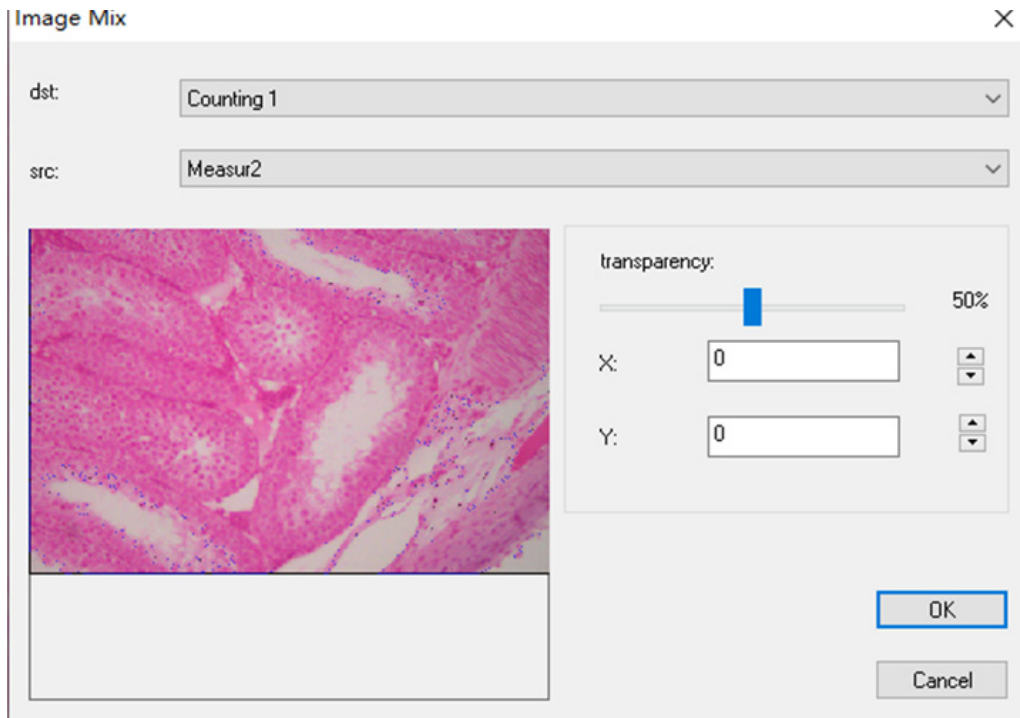You can merge two images according to different transparency ratios.


6) *Automatic Stitching* (stitching error may be caused by frame rate or roller shutter exposure):

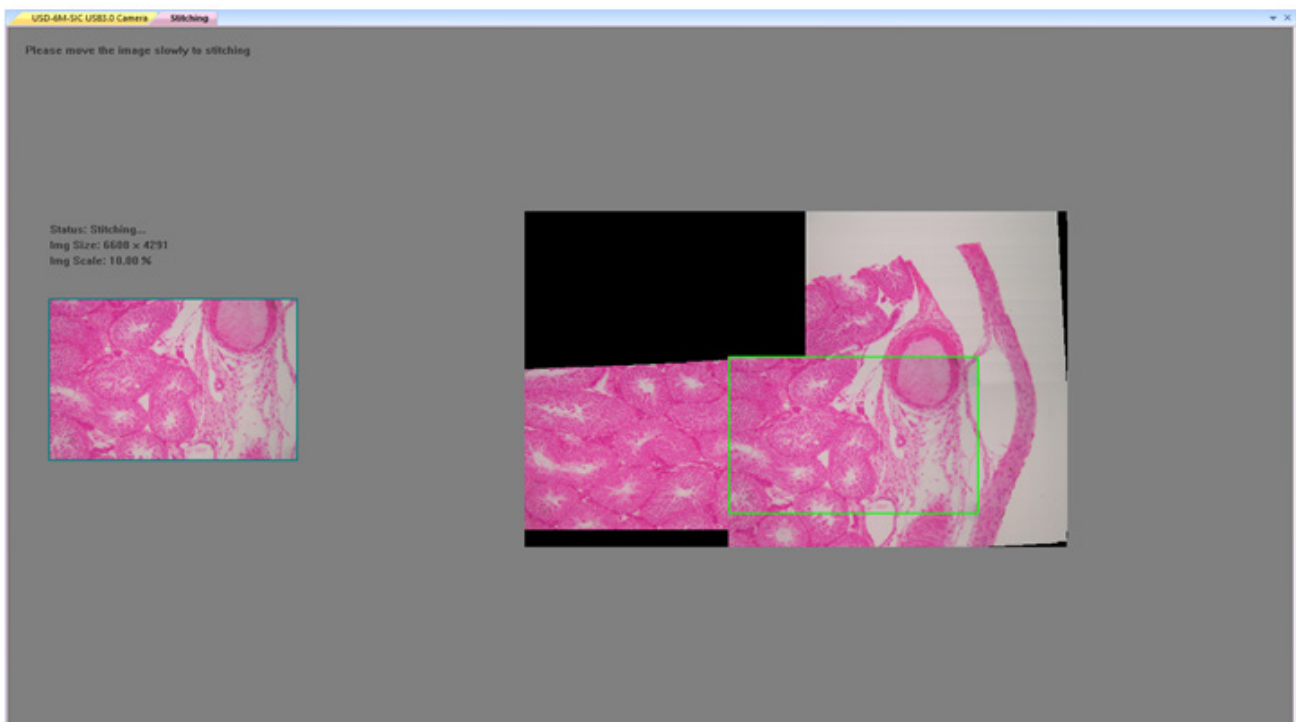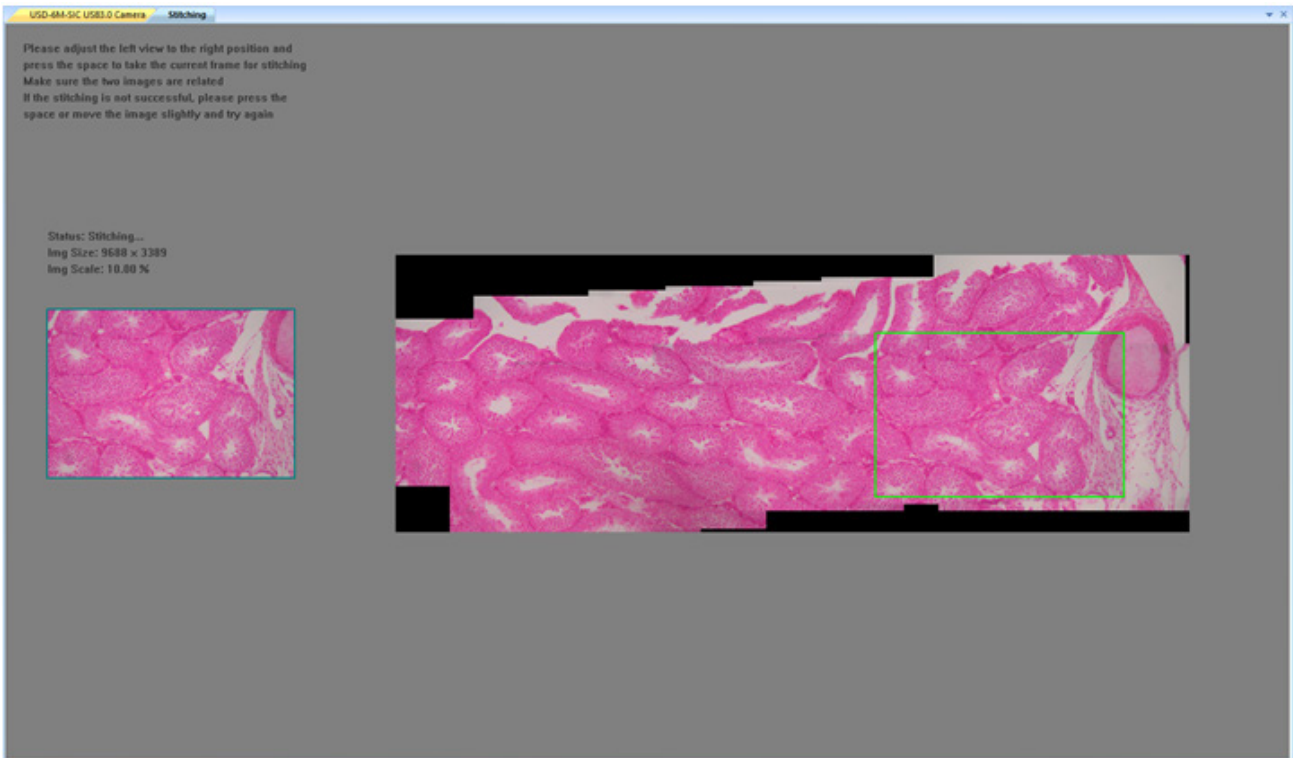It can be turned on in dynamic view, and on when auto exposure and sharpness are turned off. Image can be stitched automatically by slowly moving the stage when the two frames are related. The small image on the left is a real-time microscopic image, and the right is a mosaic image. If the stitching is not successful, please move the image back and try again.

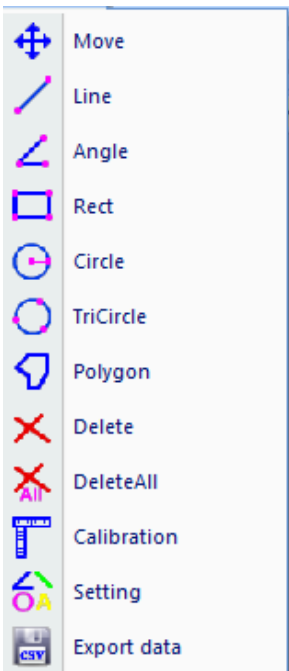
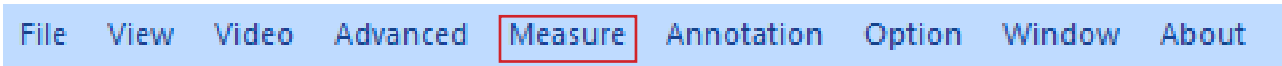
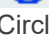
Semi automatic stitching (more accurate)

It can be opened in dynamic view, and it can be opened when auto exposure and sharpness are turned off. When the two frames are related, press the space bar to select the current frame for stitching. If the stitching is not successful, please move the image back or long press the space bar to try again.

Measure

Operation Type	Operation Instruction	Displaying Value
 Move Movement	First, select the image to be moved, now the image displays in dotted line, then move the object	
 Line Line-measurement	Select the starting point, click the left button of the mouse and keep it, then move it till the end; finally, release the mouse	Display the length of the line
 Angle Angle-measurement	Click three points in different positions with the mouse to form an angle	Display the angle
 Rect Rectangle-measurement	Select the top left corner of a rectangle, click the left button of the mouse and keep it, then move it till the lower right corner of the rectangle; finally, release the mouse	Display the area of the rectangle
 Circle Circle-measurement	Select the center of a circle, click the left button of the mouse and keep it, then move it till the point on the arc; finally, release the mouse	Display the radius
 TriCircle Three-point Circle-measurement	Click three points in different positions with the mouse to form a circle	Display the radius
 Polygon Polygon-measurement	Click any point in different positions with the mouse to form a polygon, then finish with the right mouse button	Display the area of the polygon
 Delete Delete	Click to select and delete a measurement	
 DeleteAll Eliminate	Click to delete all measurement	

 Calibration

The screenshot shows a software interface for calibration. On the left, a yellow background displays two vertical lines and a horizontal line. A blue circle with a crosshair is positioned at the intersection of the horizontal line and the left vertical line. A red circle with a crosshair is positioned at the intersection of the horizontal line and the right vertical line. A horizontal line segment connects these two red circles, with the text '100.00um' next to it. A vertical line segment extends from the bottom of the left vertical line to the horizontal line, with the text '100.00um' next to it.

On the right, a 'Ruler Setting' dialog box is open. It contains the following fields and buttons:

- Real Results:**
 - X-Ruler: um/pixel
 - Y-Ruler: um/pixel
- Physical Ruler:**
 - Ruler Unit:
 - X-Ruler Unit: um
 - Y-Ruler Unit: um
- Image Ruler(pixel):**
 - X-Start: Y-Start:
 - X-End: Y-End:
 - X-Lenth: Y-Lenth:

Buttons on the right side of the dialog box include: OK, Y = X, SaveTo, LoadFrom, and Cancel.

Unit: $\mu\text{m}/\text{mm}/\text{cm}/\text{inch}$;

X Direction: the length in the X direction of the ruler;

Y Direction: the length in the Y direction of the ruler;

Demarcation Adjustment: Move the movable points (in the blue circle) of demarcating ruler on the Image, the value of trim size (unit length per pixel) will change;

Y=X: Make X-Ruler is equal to Y-Ruler;

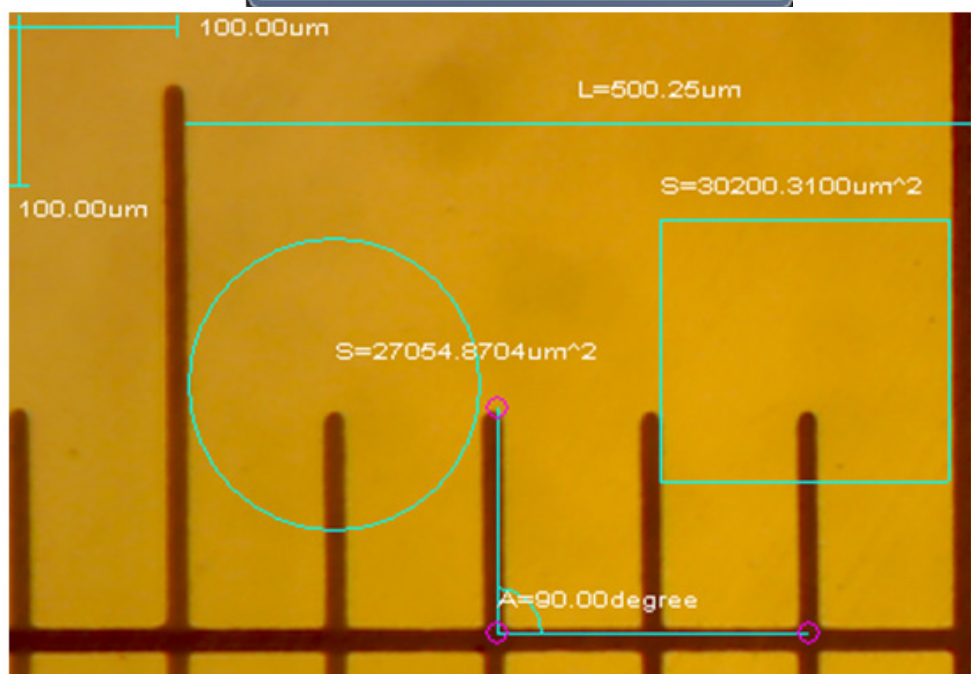
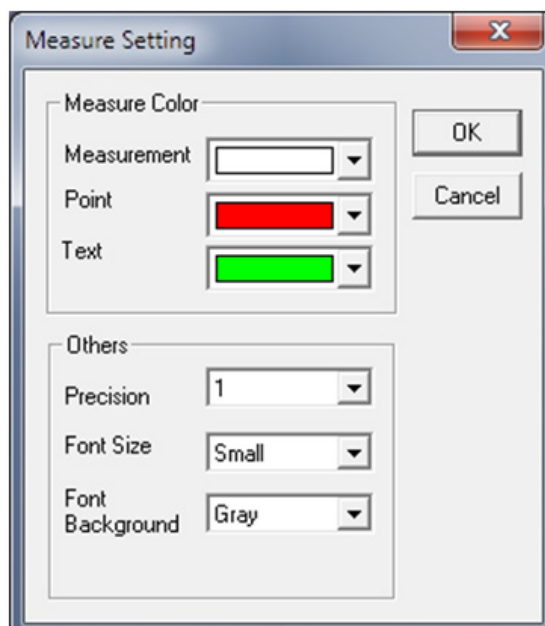
Save: Save the currently demarcated parameter;

Load: Load the demarcated-parameter file.

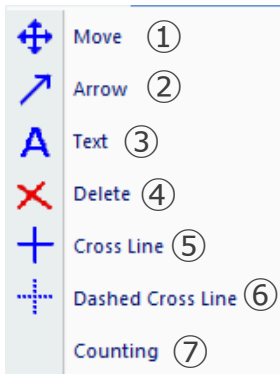
Setting

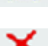
: Set the colors of measuring lines, movable points and texts.

Annotation

File View Video Advanced Measure **Annotation** Option Window About

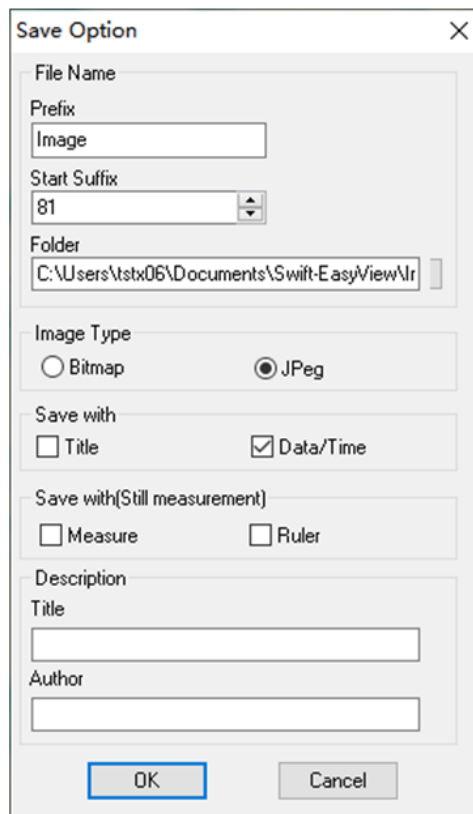
- 1)  **Move** : The added graphics and text can be moved;
- 2)  **Arrow** : Arrows can be added;
- 3)  **Text** : Text can be added;
- 4)  **Delete** : Delete added graphics and text;
- 5)  **Cross Line** : Solid cross line can be added. Click again to cancel;
- 6)  **Dashed Cross Line** : Dotted cross line can be added. Click again to cancel;
- 7) **Counting** : The serial number can be marked in the image when counting.

Option

File View Video Advanced Measure Annotation **Option** Window About

1) *Save Option*: It is used to set the parameter information when saving. See the dialog below for details.

File name: prefix + suffix + file format, such as *image1.bmp*, *image2.jpg*, etc;

File path: select location to save;

File format: BMP or JPEG file format;

Save with Title: writes the title in the description to image;

Save date / time: writes system date and time to image;

Save measurement: writes measurement data to image;

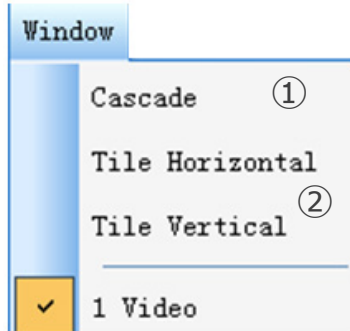
Save ruler: writes the calibration ruler to image;

Description: describes the title and author's content written to the file summary information.


2) *Language*: several languages can be selected.

Window

File View Video Advanced Measure Annotation Option **Window** About

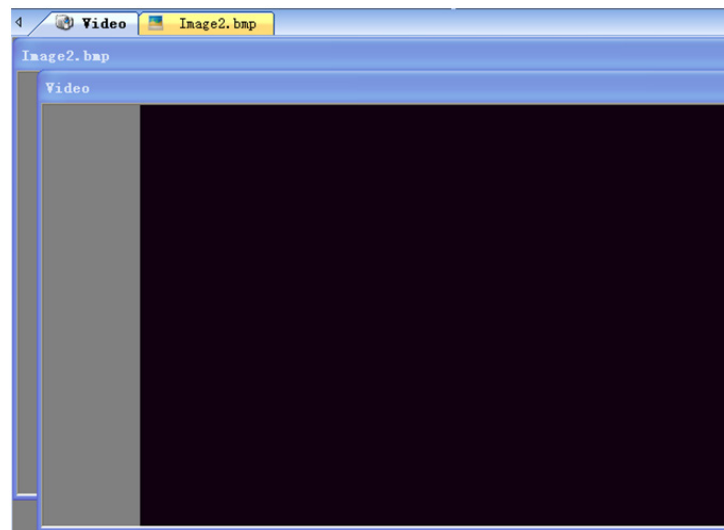
1) **Cascade** : Effects of Cascading


2) **Tile Horizontal** : Effects of Tiling

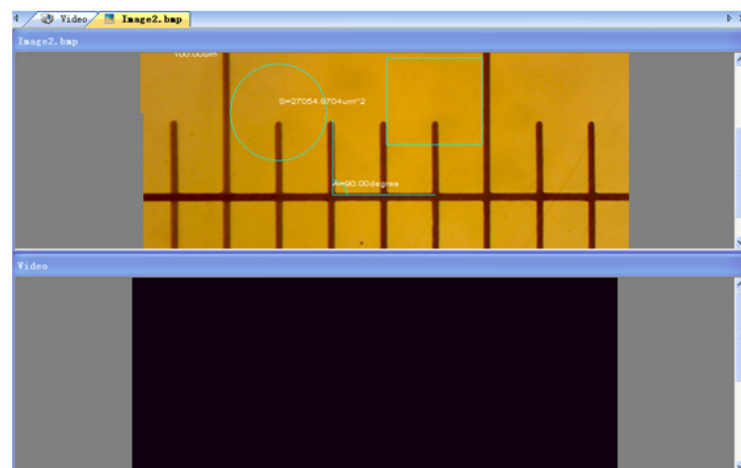
Tile Vertical

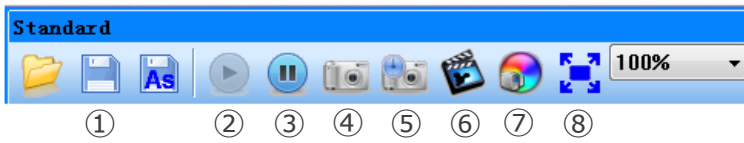b) Tool Bar

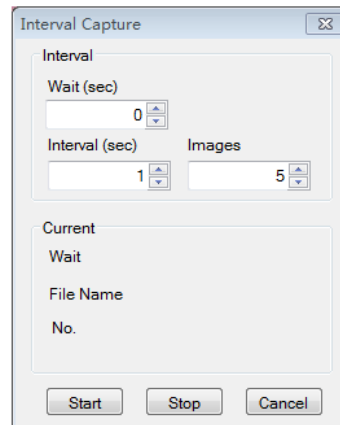- 1) : See the *File* sub menu;
- 2) : Start the live image;
- 3) : Pause the live image;
- 4) : Capture image and display it in a new window;
- 5) : Interval Capture

Wait (sec): Set after how many seconds the system starts acquisition;
Interval: Set the acquisition interval between images;
Images: The number of images to be taken.

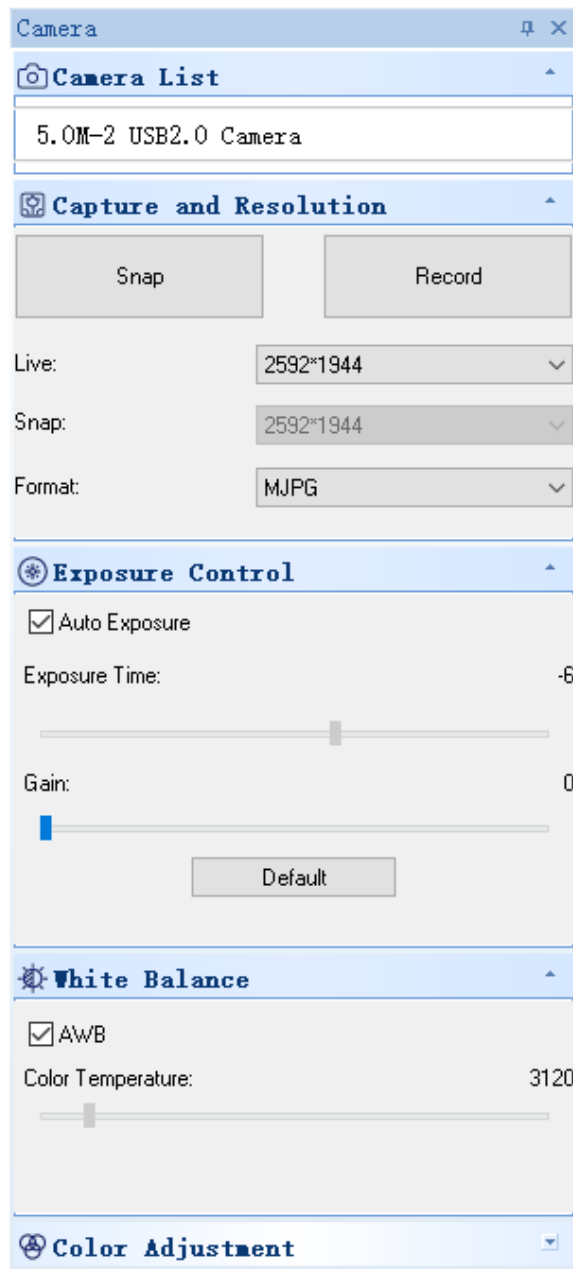
- 6) : Record

Path: Select the address for saving image files and file names;
Start: Start recording;
Pause: Pause the video;
Stop: Stop recording and the video will be automatically saved in the selected path.


- 7) : Image attribute adjustment, see "Image adjustment" in Image sub menu.
- 8) : Display live image in full screen mode. (Or double click live display window to enter into or quit full screen mode).

8.1.3 Sidebar


1. Camera List.

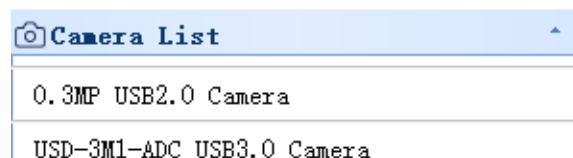When more than one camera is connected to the computer, the list of available sources appears. Select the one you are interested in.

2. Capture and Resolution.

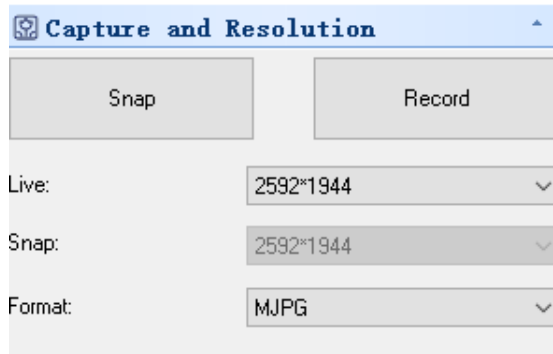
You can take photos and record images. Here you can set the resolution of live image, the resolution of snapshot and the image format.

3. Exposure Control.

The default is *Auto Exposure*.

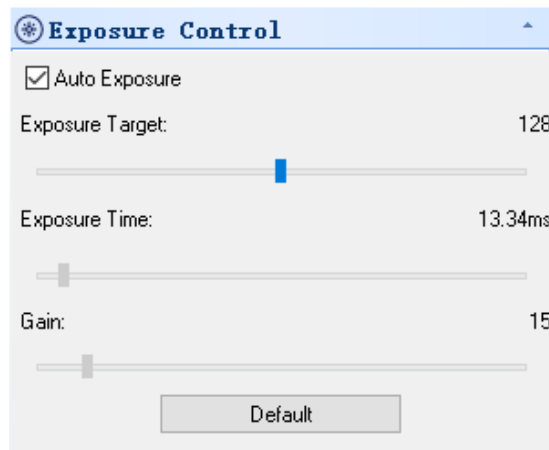
If you deselect the *Auto Exposure* check box you can manually set the exposure time and gain by yourself. Click *Default* to restore the default *Auto Exposure*.

4. White Balance.


RGB color temperature can be adjusted.

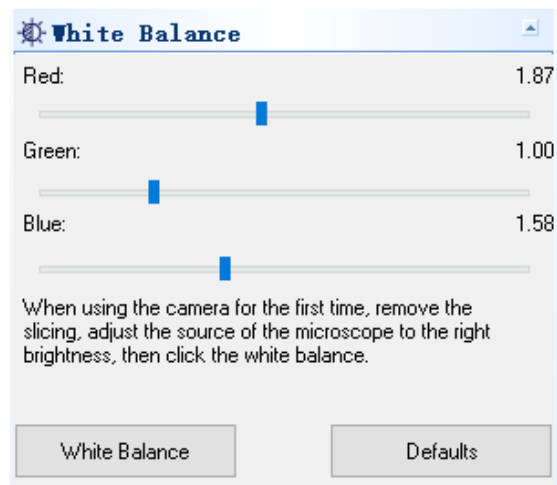Click on the *White Balance* button to set the desired white balance. Click on *Defaults* to restore the default value.

5. Color Adjustment.

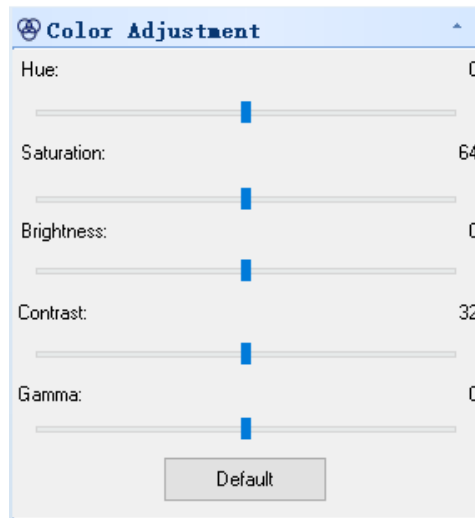Adjust the hue, saturation, brightness, contrast and gamma value of the image. Click the *Default* button to restore the default value.

8.2 USB/HDMI C-mount camera 21MP

The camera is driven by S-EYE (either for Windows and for IOS) when used in USB mode.

- The software can be downloaded from this link: http://www.optikamicroscopes.com/OPTIStar/S_EYE.zip

When used in HDMI mode the camera has a built-in software.

- **NOTE: no driver installation is required for the cameras. The software setup procedure automatically installs all the needed drivers for the correct operation of the camera.**

8.2.1 Use of camera in HDMI mode

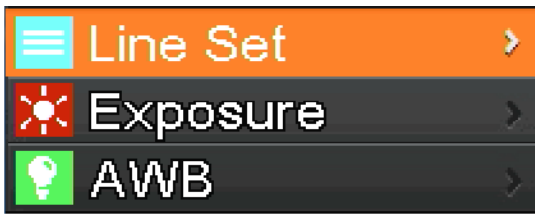
a) Language selection

1. Press *MENU* on the camera to enter the main menu, and then press *DOWN*; the *Language* option is highlighted.
2. Press *OK* to enter the language settings, use *DOWN* to select *English*, press *OK* to return to the main menu.

b) Line setting

1. Press *MENU* on the camera to enter the main menu, and then press *DOWN*; the *Line Set* option is highlighted.
2. Press *OK* to enter the line setting menu.

3. You can use *DOWN* and *UP* to select options that you want to set.

4. When the entry *Cross Hair* is highlighted, press *OK* to turn on and off the central cross line.

5. When the *Line* entry is highlighted, press *OK* or *MODEL/AE* to select number of lines (1-8).

Cross Hair	ON
Line	2
Direction	OFF
Position	100
Color	Red
Thickness	1

6. When the *Line Direction* entry is highlighted, press *OK* or *MODEL/AE*; you can choose line direction (horizontal, vertical or none).

Cross Hair	ON
Line	2
Direction	V
Position	100
Color	Red
Thickness	1

7. When the *Line Position* entry is highlighted, press *OK* or *MODEL/AE* to move the line.
- To quickly move the line keep pressed the key.

Cross Hair	ON
Line	2
Direction	V
Position	100
Color	Red
Thickness	1

8. When the *Line Color* entry is highlighted, press *OK* or *MODEL/AE*; the color of the line can be changed. Five colors are available: Red, White, Green, Black and Blue.

Cross Hair	ON
Line	2
Direction	V
Position	100
Color	White
Thickness	1

- When the *Line Thickness* entry is highlighted, press *OK* or *MODEL/AE*; the line thickness can be changed. Four types of thickness can be selected.

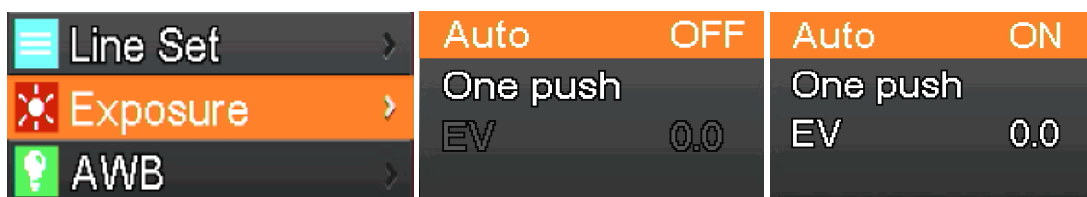
c) Exposure setting

- To quickly turn off automatic exposure, long press *MODEL/AE*.

- Press *MENU* on the camera to enter the main menu, press *DOWN* to move to the *Exposure* option, then press *OK* to enter the exposure settings. You can set the automatic exposure on/off and exposure values through the menu.
- Enter the *Exposure* menu as shown below. Use *UP* and *DOWN* to select the *Auto* option highlighting. Use *OK* and *MODEL/AE* to select automatic exposure ON or OFF.
- Select *One push*, and then press *OK*; the camera will perform a key exposure. After the key exposure, automatic exposure will be closed.

- Enter the *Exposure* menu as shown below. Select the exposure value (EV) by using *UP* and *DOWN*. Use *OK* and *MODEL/AE* to adjust the corresponding value.
- NOTE: the exposure value is allowed to be adjusted only when the automatic exposure is OFF.**

d) White Balance setting

1. Press *MENU* on the camera to open the main menu, press *DOWN* to move to the *AWB* (Auto White Balance) entry, then press *OK* to enter the *White Balance* menu.
2. Use *OK* and *MODEL/AE* to select the *AWB* (Automatic White Balance) ON or OFF.

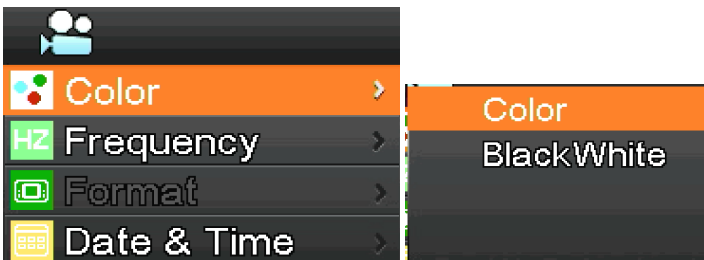
3. Enter the *White Balance* settings menu as shown below.
4. After disabling the Automatic White Balance, you can adjust Red, Green and Blue values.
5. Select items that need to be adjusted by using *UP* and *DOWN*.
6. Use *OK* to increase the value, use *MODEL/AE* to decrease it.
7. Select *One push*, and then press *OK*. The camera will operate a “one shot” white balance, and the key exposure after the completion of automatic white balance will be closed.

e) Color setting

1. Press *MENU* on the camera to open the main menu, press *DOWN* to move to *Color* option, then press *OK* to enter the *Color* menu.
2. Use *UP* and *DOWN* to select the color mode or black-and-white mode.
3. Press *OK* to confirm and exit or press *MENU* on the camera to abort operation and return to the main menu.

f) Light frequency setting

1. Press *MENU* on the camera to open the main menu, press *DOWN* to select the *Frequency* entry, then press *OK* to enter the *Light Frequency* menu.
2. Use *UP* and *DOWN* to select 50Hz or 60Hz.
3. Press *OK* to confirm and exit or press *MENU* on the camera to abort operation and return to the main menu.

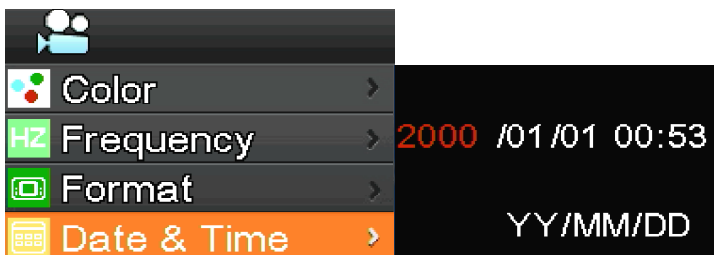
g) Formatting the SD card

1. Press *MENU* on the camera to open the main menu, press *DOWN* to move to the *Format* option, then press *OK* to enter the *Format* menu.
 2. Choose *No* or *Yes* by using *UP* or *DOWN*.
 3. Select *No* and press *OK* to abort the operation and return to main menu.
 4. Select *YES* and press *OK*: the camera will format SD card.
- **NOTE: the original file content in the SD card will be removed after the formatting operation. Please ensure the content of SD card has been backed up before this operation!**


h) Date and Time setting

1. Press *MENU* on the camera to open the main menu, press *DOWN* to select the *Date & Time* option, then press *OK* to enter the *Date & Time* menu.
2. Press *OK* to switch and select the items that need to be adjusted (such as the date).
3. When the item to be adjusted is red, press *UP* and *DOWN*: the red one indicates that you can adjust the value of selected item.
4. After adjustment, press *MENU* on the camera to save changes and return to the main menu.


i) Image size setting (size of image in recording)

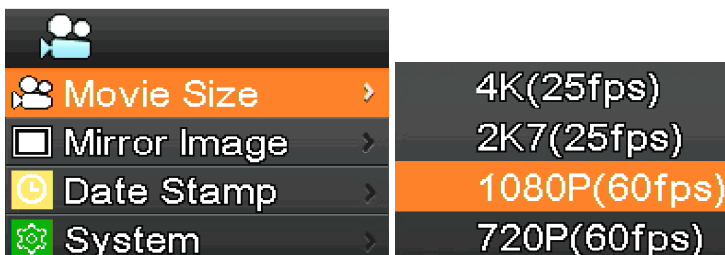1. Press *MENU* on the camera to open the main menu, press *DOWN* to select the *Movie Size* option, then press *OK* to enter the *Movie Size* menu.
2. Select required size by using *UP* and *DOWN*, press *OK* to save and return to the main menu.


j) Mirror image setting

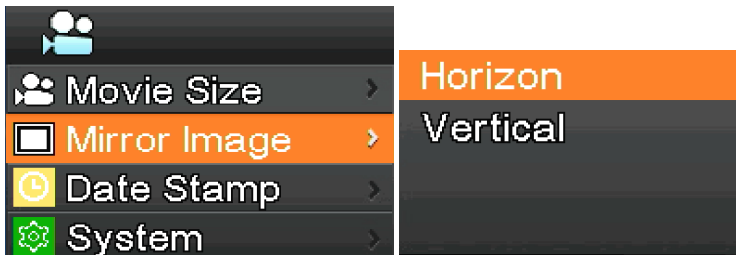1. Press *MENU* on the camera to open the main menu, press *DOWN* to select the *Mirror Image* option, then press *OK* to enter the *Mirror Image* menu.
2. Use *UP* and *DOWN* to select the *Horizontal* or *Vertical* mirror settings. Press *OK* to select *ON* or *OFF*, press *MENU* on the camera to save and exit.


k) Date Stamp setting

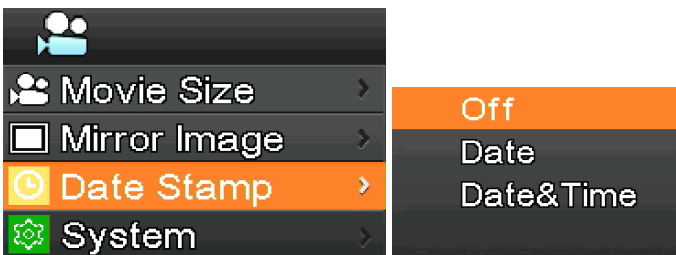1. Press *MENU* on the camera to open the main menu, press *DOWN* to select the *Date Stamp* option, then press *OK* to enter the *Date Stamp* menu.
2. Use *UP* and *DOWN* to select *Off*, *Date* or *Date & Time*, press *OK* to confirm and exit.
 - If you select *Off*, the image will not show the time at which it was captured.
 - If you select *Date*, the image will show in overlay the date on which it was acquired.
 - If you select *Date & Time*, the image will show in overlay the date and the time on which it was acquired.

l) System setting

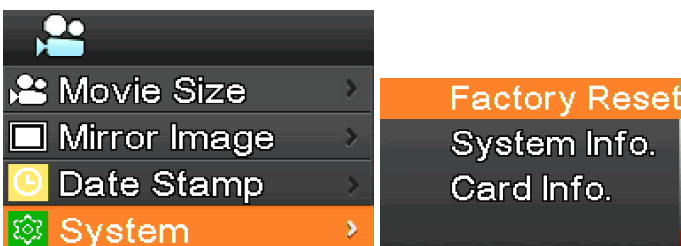1. Press *MENU* on the camera to open the main menu, press *DOWN* to select the *System* option, then press *OK* to enter the *System* menu.
2. Use *UP* and *DOWN* to select *Factory Reset*, *System Info.* and *Card Info.*
3. When selecting *Factory Reset*, and pressing *OK*, a dialog will remind the user whether to restore the factory settings. Use *UP* and *DOWN* to select *Yes* or *No*, press *OK* to confirm.
4. When you select *System Info.* and pressing *OK* again, the current version of the software is displayed. Press *OK* to accept, Press *MENU* on the camera to return to the main menu.
5. When selecting *Card Info.* and pressing *OK* again, the current SD card information are displayed. Press *OK* to accept, Press *MENU* on the camera to return to the main menu.

m) Electronic enlargement and reduction

1. Press *UP*, the image will be zoomed in electronically.
- On the left of the image, the zoom factor will be displayed. A maximum of 5X of electronic zoom is allowed.
2. After the electronic zoom, press *DOWN*. The image will be zoomed out.
- On the left of the image, the zoom factor will be displayed until the original size of the image is restored.

n) Switch camera / video mode

1. The system starts as defaults to camera mode. In this mode press *OK* to take pictures.
2. Press *MODEL/AE* to switch to video mode. In this mode press *OK* to start recording video, and then press *OK* again to stop recording.

o) Remote control and keys


	Zoom in		Up		Zoom out
	Left		MENU NENU		Right
	Mirror		Down		AE ON/OFF
	Flip		Center cross		AWB ON/OFF
	Mono		Frozen		Play/pause
	Photograph		Record		PAIayback

8.2.2 Use of camera in USB mode


1. Run S-EYE by clicking the icon  on the desktop.
2. Click on the camera name in the section *Camera List* on the upper right corner of the screen.
3. Live image starts automatically.

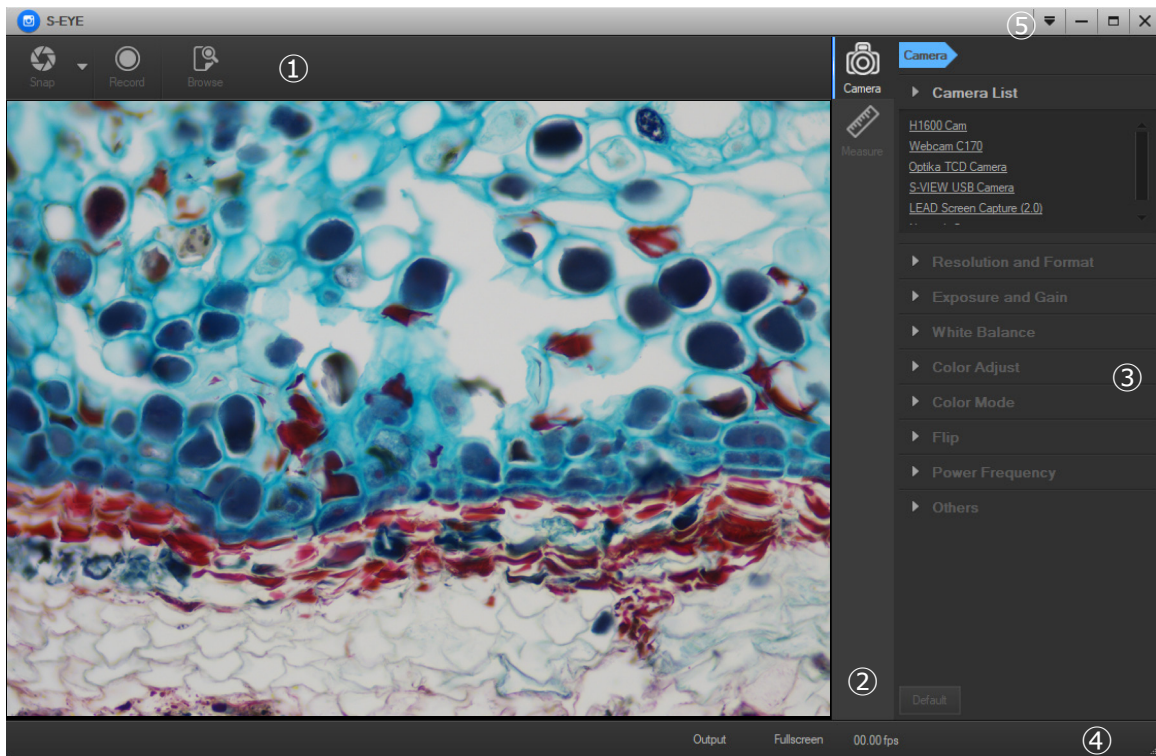a) User Interface


1. *Toolbar*: Take pictures/Recording image/Browse files.
2. *Sidebar*: Switching the rightmost control panels, currently provides camera control and measurement functions.
3. *Control panel*: Specific camera control and measurement operations.
4. *Status bar*: Display camera information, change user interface view, toggle measure result panel.
5. *Main menu*: Open setting dialog, show software information, etc.

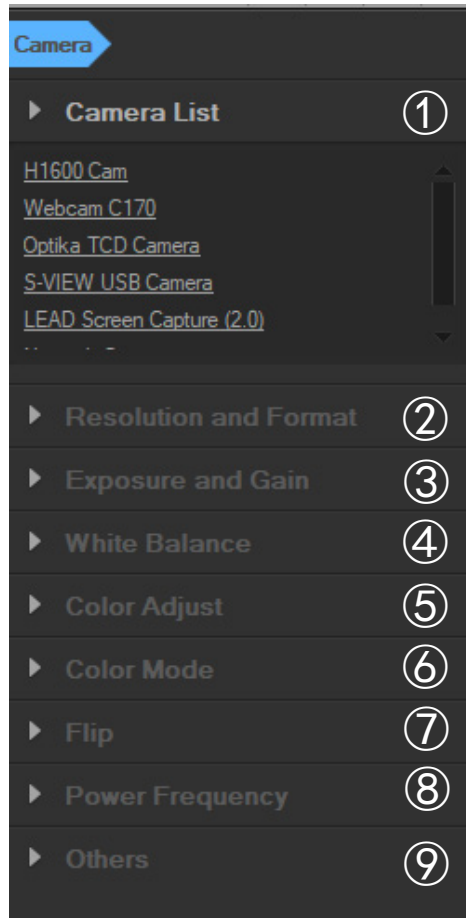b) Camera Control Panel


1. Camera List:

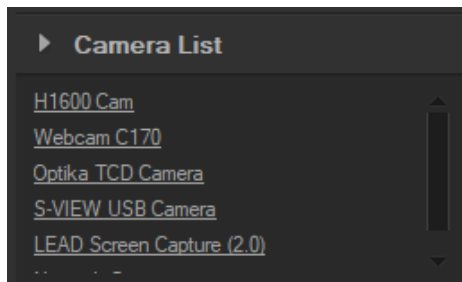When more than one camera is connected to the computer, the list of available sources appears. Select the one you are interested in.


2. Resolution and Format:

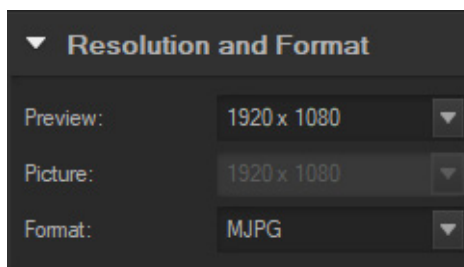User can take images and record video. Here you can set the resolution of live image, the resolution of snapshot and the image format.

3. Exposure and Gain:

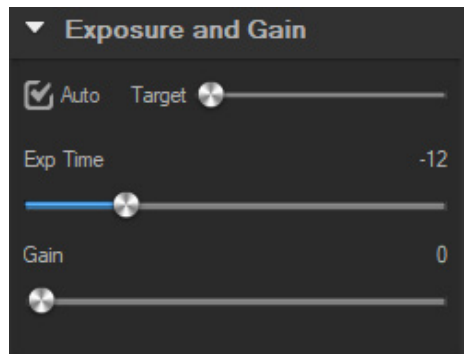
When *Auto* check box is selected, the camera will detect the best exposure time. When *Auto* check box is deselected, the user can manually set the exposure time.

4. White balance:


When *Auto* check box is selected, the software will detect the automatic white balance.

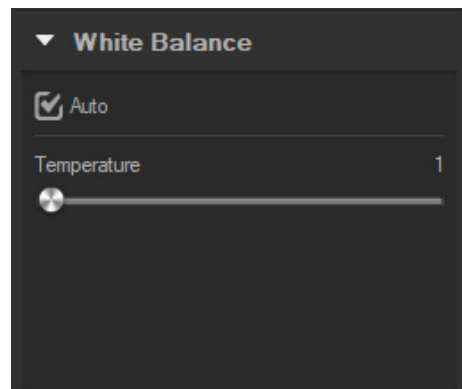When *Auto* check box is deselected, the user can change the color temperature operating on the cursor *Temperature*.



5. Color Adjust:

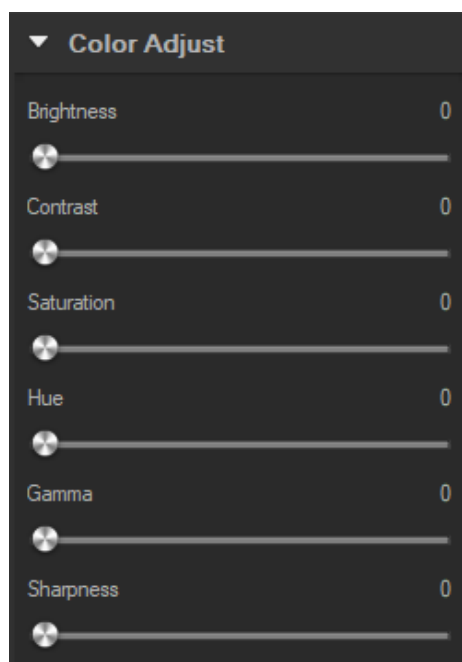Adjust the hue, saturation, brightness, contrast, gamma and sharpness value of the image.


6. *Color mode:*

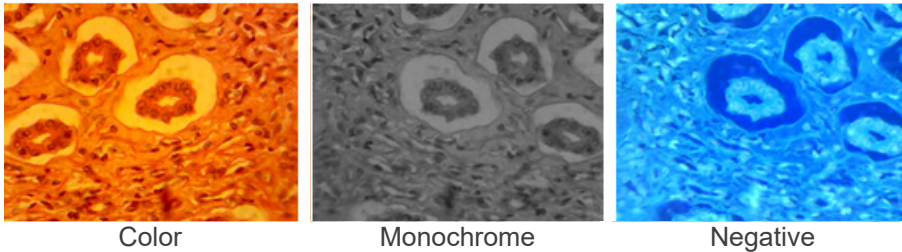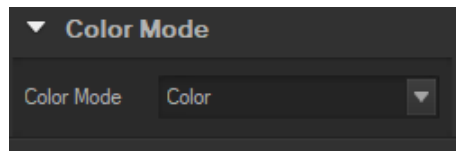Available options are *Color*, *Monochrome* or *Negative*.


7. *Flip:*

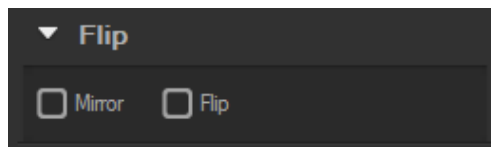Flip image vertically and horizontally.

8. *Power frequency:*

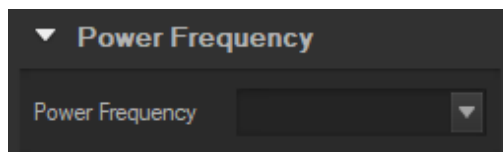To avoid image flickering, user may need change this parameter.


9. *Default:*

Reset camera parameters.


- **NOTE:** If the selected camera does not support some features, the corresponding interface controls will not be operating.

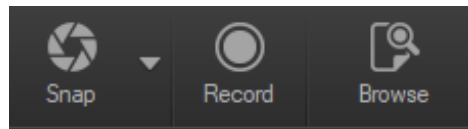c) Capture Images


User can capture images and videos.

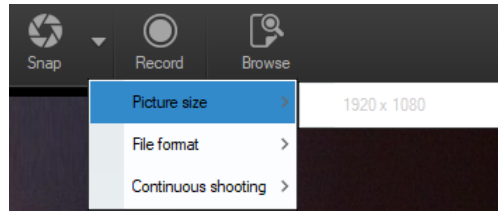1. Take an image: click *Snap* button to take an image.
Click the right arrow beside the *Snap* button to set *Picture size*, *File format* and *Continuous shooting*.

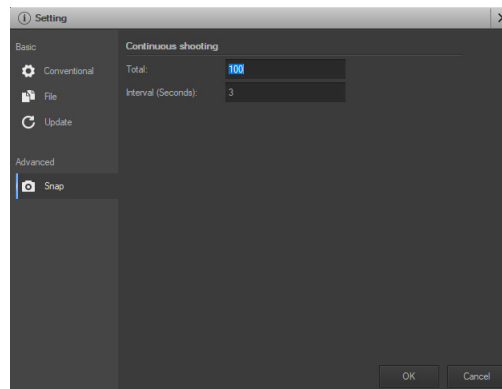When *Continuous shooting* is selected, user can start the Time Lapse acquisition or set the acquisition parameters. Selecting *Setting* the following dialog box appears:


Total: The number of images to be taken.

Interval (seconds): after the acquisition of one image, several seconds later, another image is available to be taken.


2. Recording: click *Record* button to start image recording, click again to stop recording. When recording, an animation indicator appears over the image.
3. Browse: click *Browse* button to open the folder where images and image are stored.

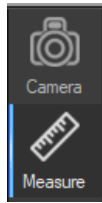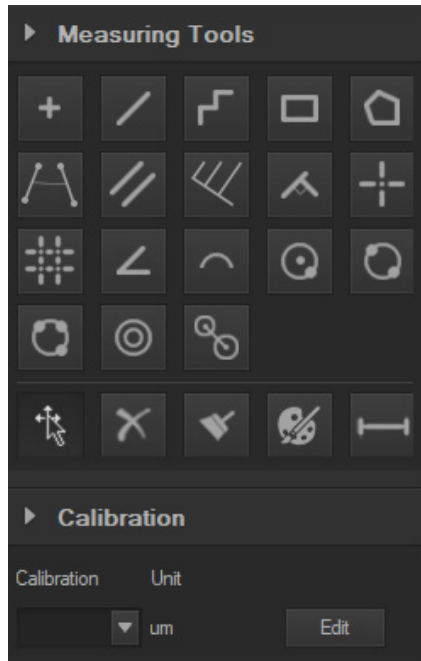d) Measurement

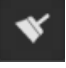
Click on *Measure* icon in the Sidebar.

The rightmost sidebar changes accordingly.


Icon	Function	Description
	Move	Move or adjust the measurement tool
	Point	Draw a marker over the image
	Line	Measure length of a line
	Break line	Measure length of break line, right click to finish drawing
	Rectangle	Measure perimeter and area of a rectangle
	Polygon	Measure perimeter and area of a polygon
	Circle	Measure a circle, the circle is determined by center point and radius
	Two Points Circle	Measure a circle, the circle is determined by two points



	Three Points Circle	Measure a circle, the circle is determined by three points
	Arc	Measure length of arc
	Angle	Measure angle
	Perpendicular	Measure the distance from a point to a line
	Delete	Delete one measurement
	Delete all	Delete all measurements
	Export	Export measurement result to file

e) Calibration

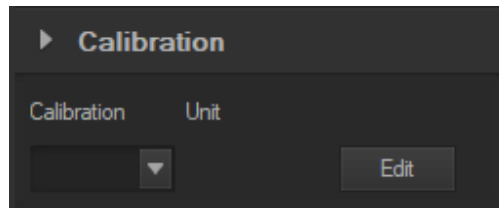1. Click on *Calibration* in the Sidebar.

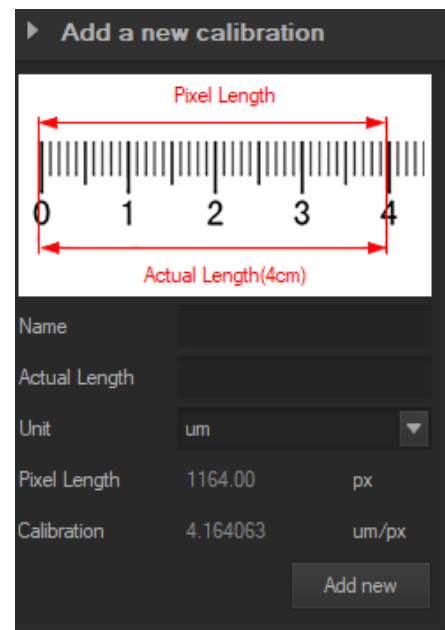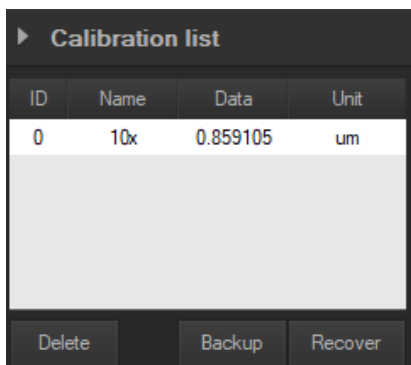2. Click on *Edit*. The live image will be added of a red ruler and the Sidebar changes accordingly.
 - The upper part is the *Calibration List*, where the existing calibration points are available.
 - The lower part is the *Add a new calibration* dialog where a new calibration can be performed.

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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The image shows a 'Save Option' dialog box with the following fields and options:

- File Name:**
 - Prefix: Image
 - Start Suffix: 81
 - Folder: C:\Users\tstx06\Documents\Swift-EasyView\Mr
- Image Type:**
 - Bitmap (unselected)
 - JPeg (selected)
- Save with:**
 - Title (unchecked)
 - Data/Time (checked)
- Save with(Still measurement):**
 - Measure (unchecked)
 - Ruler (unchecked)
- Description:**
 - Title: [empty text box]
 - Author: [empty text box]

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